
making a difference

BIOSCIENCE CATALOGUE

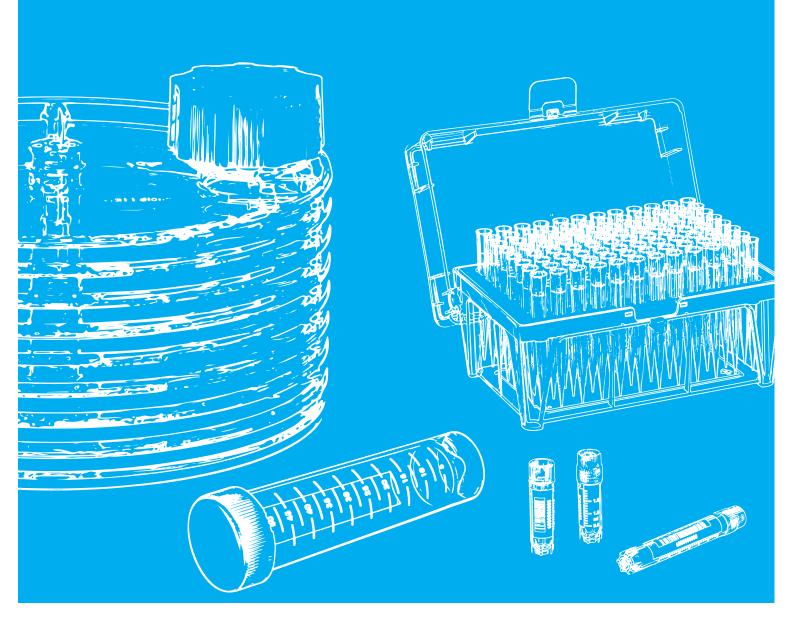

EDITION 03/2022



making a difference

To us, health is the highest personal priority.
This inspires our daily mission of serving life sciences and healthcare professionals with advanced and sustainable products and solutions.

It is the passion and dedication of our people that make the difference. With a straightforward approach our global team continuously strives for excellence and closely collaborates with our customers.

We empower people to make a difference in health.

トスツと

6 8 10 12









BIOSCIENCE

Making a difference in research and science

PREANALYTICS

Preanalytical specimen collection systems and safety solutions

OEM

Tailor made solutions

SERVICE

E-Commerce Solutions from Greiner Bio-One

14 30 240 272



VALUE PROPOSITIONS

About our value propositions



PRODUCTS

Product overview




TECHNICAL APPENDIX




NUMERIAL INDEX

出 い Z M O S O M

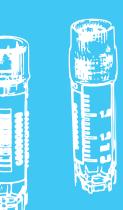



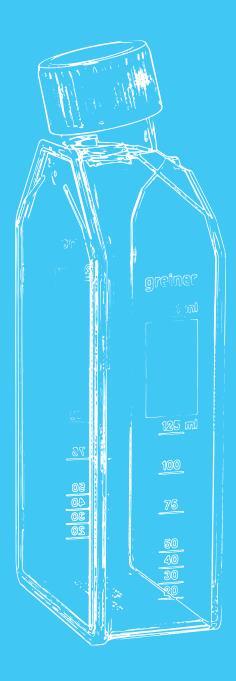

MAKING A DIFFERENCE IN RESEARCH AND SCIENCE


Greiner Bio-One is one of the leading suppliers of special products for the cultivation and analysis of cell and tissue cultures. Under the brand name CELLSTAR®, we offer cell culture flasks, dishes and plates in a wide variety of formats and surface modifications, so that you will always find exactly the right product for every application and all cell types.

In addition, we have been developing and producing microplates for high-throughput screening for more than 50 years, ena-

bling industry and research to carry out fast and most efficient drug testing. Among other things, Greiner Bio-One was the first manufacturer to introduce a microplate in 1536 well format to the market, setting new standards in terms of automation, performance and cost reduction.

Based on decades of experience in the cryogenic storage of samples, we also offer comprehensive solutions for automated storage systems in biobanks.

- / High-quality products for your lab
- / Special products for cell and tissue culture
- / Microplates for industry and research
- / Cryogenic storage systems

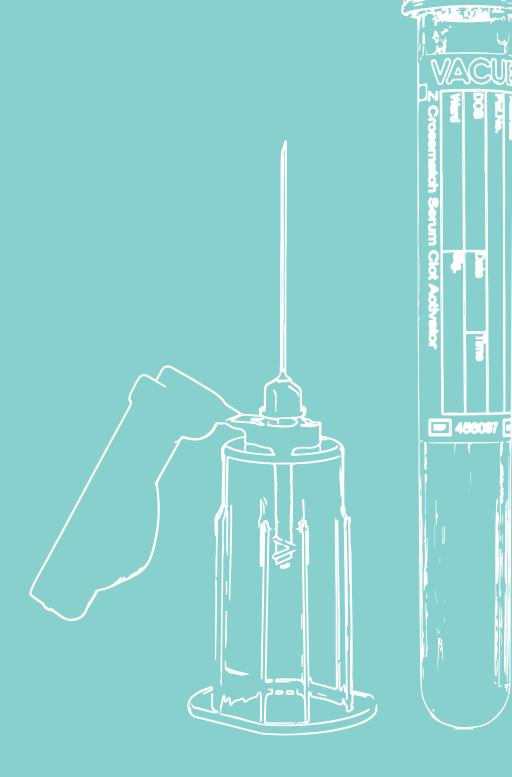

PREANALYTICAL SPECIMEN COLLECTION SYSTEMS AND SAFETY SOLUTIONS


Greiner Bio-One was the first company in the world to successfully launch a blood collection system made out of virtually unbreakable PET plastic including a safety cap. Many components of the VACUETTE® System have often been copied, but never reached our high quality standards.

As a trendsetter in the field

of sample collection, we offer solutions that meet the highest demands of our customers. Because your work is an important contribution to health protection.

Continuously new and improved product solutions, perfect functionality, maximum safety and high product quality are the result of decades of experience.

- / Provide safety
- / Simplify handling
- / Provide support
- / Increase efficiency
- / Save costs

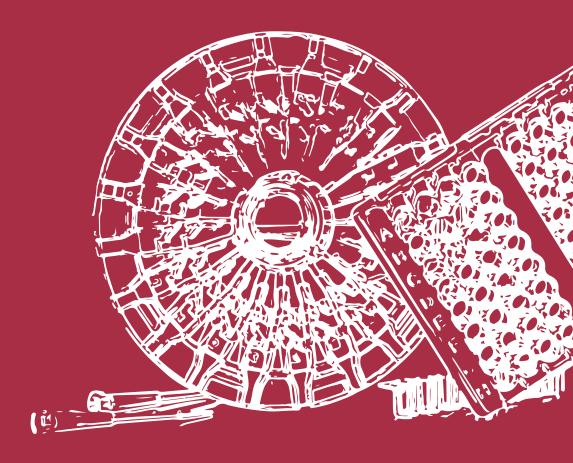


TAILOR MADE SOLUTIONS

Greiner Bio-One as an Original Equipment Manufacturer (OEM) is a long-term partner of the pharmaceutical industry, biotechnology, diagnostic and medical technology industries. The company manufactures numerous products through injection moulding and offers the entire product development and production process. From design,

via prototyping, through fully automated manufacturing, Greiner Bio-One delivers the complete solution with support of experienced scientists, engineers and specialists.

With a global sales network specifically established for the field of OEM, Greiner Bio-One offers custom-made solutions with personal service.

- / Research & development with diverse application and research laboratories
- / Construction of moulds, planning of plants and assembly lines
- / Modern, fully automated production
- / Post-production according to customer requirements
- / Comprehensive quality assurance and management system
- / Worldwide networked warehousing and customer service

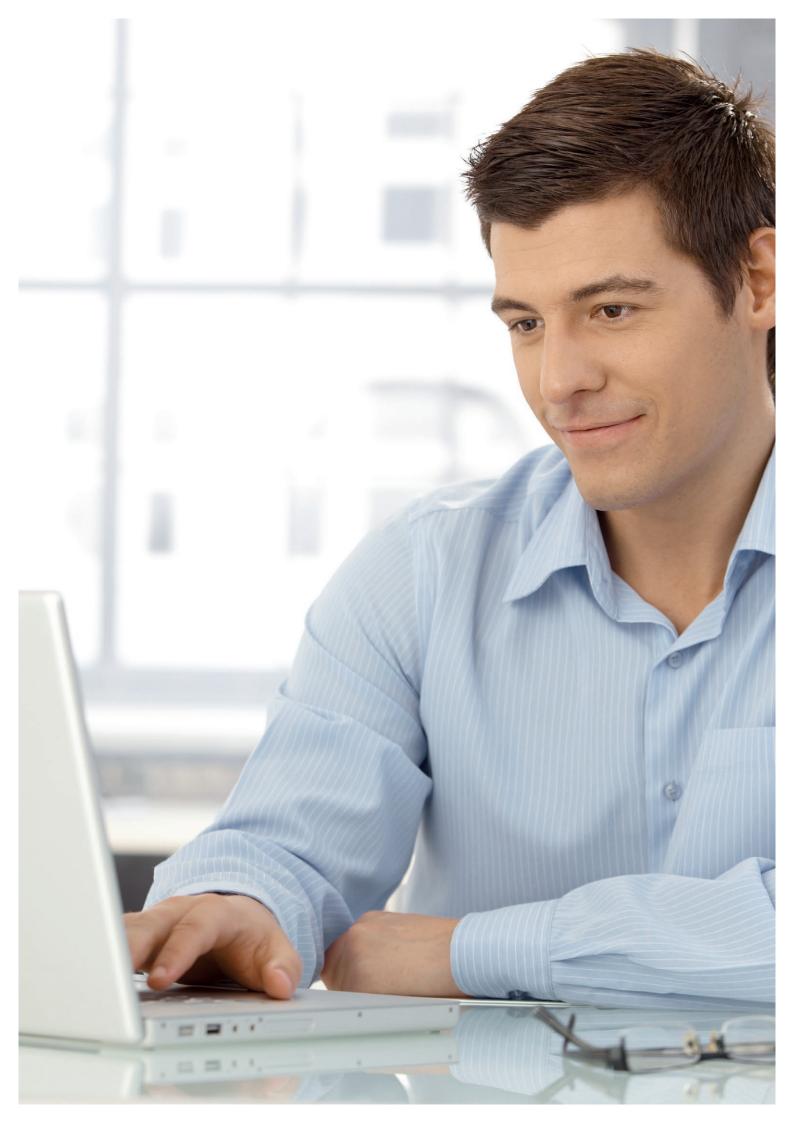

FAST. SECURE. EFFICIENT. E-COMMERCE SOLUTIONS FROM GREINER BIO-ONE

As your strong partner in the fields of biotechnology, diagnostics, medical devices and research, Greiner Bio-One already provides you with solutions on a daily basis. In addition, we can offer you a wide selection of e-business solutions to ensure the optimal integration of our products and services into your purchasing process. That in turn will significantly reduce your process-related costs and warehouse costs.

Take advantage of e-business connectivity and make your processes faster, more secure and more efficient.

Greiner Bio-One offers you numerous solutions and the necessary expertise to support your processes, regardless of which ERP system you use. Would you like to purchase via our online shop or are you interested in one of our eProcurement solutions such as EDI or OCI? Contact our experts: edi@de.gbo.com. In order to stay informed, you can subscribe to our Greiner Bio-One Newsletter and never miss any news about our products and events.



- / Online-Shop
- / EDI (Electronic Data Interchange)
- / OCI (Open Catalog Interface)
- / Electronic catalogs



We can find the right solution for you in all standard formats.

Please contact our specialists at: edi@de.gbo.com.



TRUE TRUST

Take advantage of our experience and passion for your projects and tasks

WE KEEP OUR PROMISES

For us, real trust means keeping promises. This is one of the core values that we are committed to as a company, both towards you as our customers and towards our employees.

As a family-owned company, we believe that trusting, active partnership is the key to sustainable and reliable processes. We

understand your requirements, speak your language and are happy to support you in advancing your scientific progress. We are at your side at every stage our competent team of experts is always on hand, from the first contact and flawless delivery of our products, to all forms of after-sales service.



PUT YOUR TRUST IN US:

- / Partner for sustainable and reliable processes
- / 50 years of experience in science and healthcare
- / Innovation with tradition
- Support during every work step



DISCOVER MORE:





QUALITY

Reproducible results you can rely on

CONSISTENTLY HIGH QUALITY FOR RELIABLE RESULTS

Quality has been our top priority for 60 years. This applies all over the world, as all of our production facilities are committed to the same high standard.

Our products are made from strictly controlled, certified raw materials. With state-of-theart production technology and

continuous in-process checks by specially trained employees, we ensure consistent quality of the products. Quality promise not only relates to the manufacturing process and therefore to product quality, but to every process and every interaction with you. We are ISO 9001 and ISO 13485 certified.




CONSISTENT HIGH QUALITY:

- / Locationindependent
- / State-of-the-art manufacturing technologies
- / Controlled and certified raw materials
- / ISO 9001 and ISO 13485 certified



DISCOVER MORE:

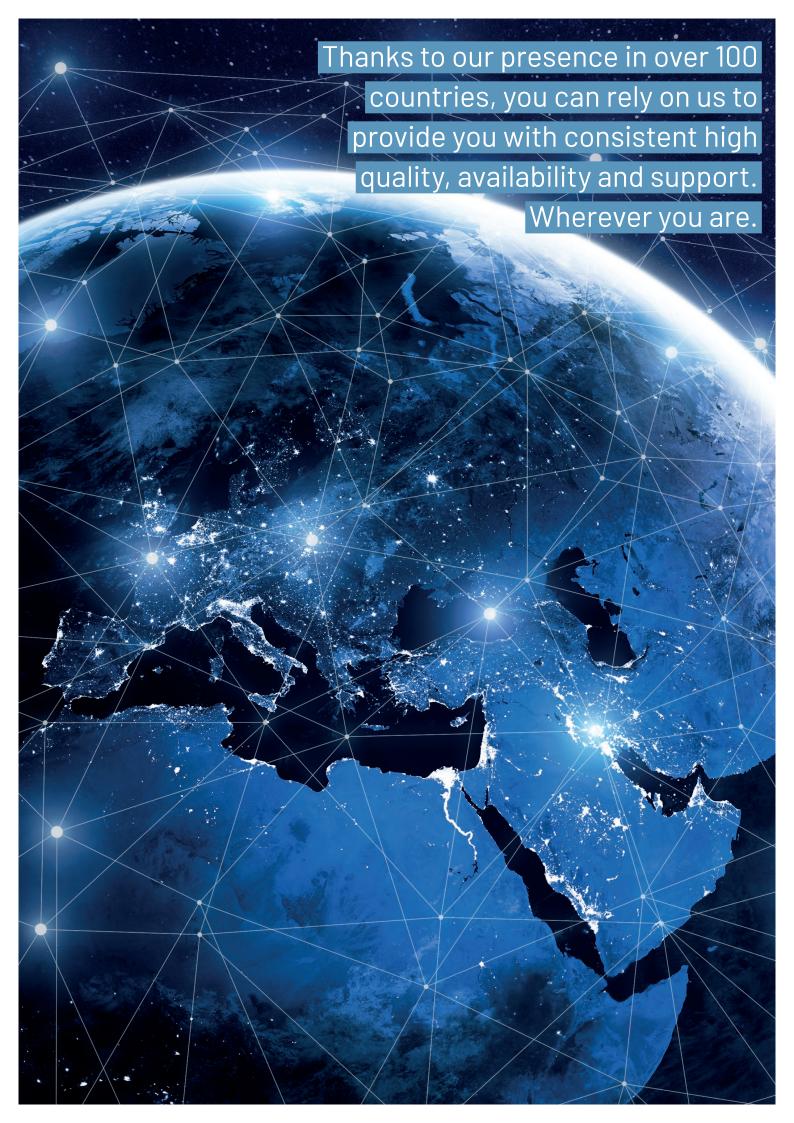

GLOBAL PRESENCE

There to support you - wherever you are

ON-SITE FOR YOU WORLDWIDE

By global presence we mean being there for you whenever and wherever you need us.

Our seven production sites in Europe, Asia and North and South America meet the same high quality standards and ensure low-risk delivery routes and a secure product supply. The support teams of our 28 offices and distributors in more than 100 countries are helpful and competent, and are always on hand when you need help or support.



WE PLACE IMPORTANCE ON:

- / Competent local support
- / Optimum product availability
- / Consistent high quality worldwide









IDEAS BECOME SOLUTIONS

For more than 50 years, our passion for plastic technology and manufacturing processes has driven us to constantly innovate to support you optimally in your daily work and ensure your success.

We have been instrumental in the development of consumables in the diagnostics and life sciences industries, as well as in the medical market, since 1963. We are proud of our rich pioneering history – from the production of the first plastic petri dishes, to Europe's first plastic vacuum sampling system or the invention of the world's first 1536-well microplates. We will continue to strive for new and innovative product solutions that are precisely tailored to your requirements and optimally support you in your daily work.




PROGRESS NEEDS PIONEERS:

- Product solutions for current and emerging trends
- / Innovative production technologies





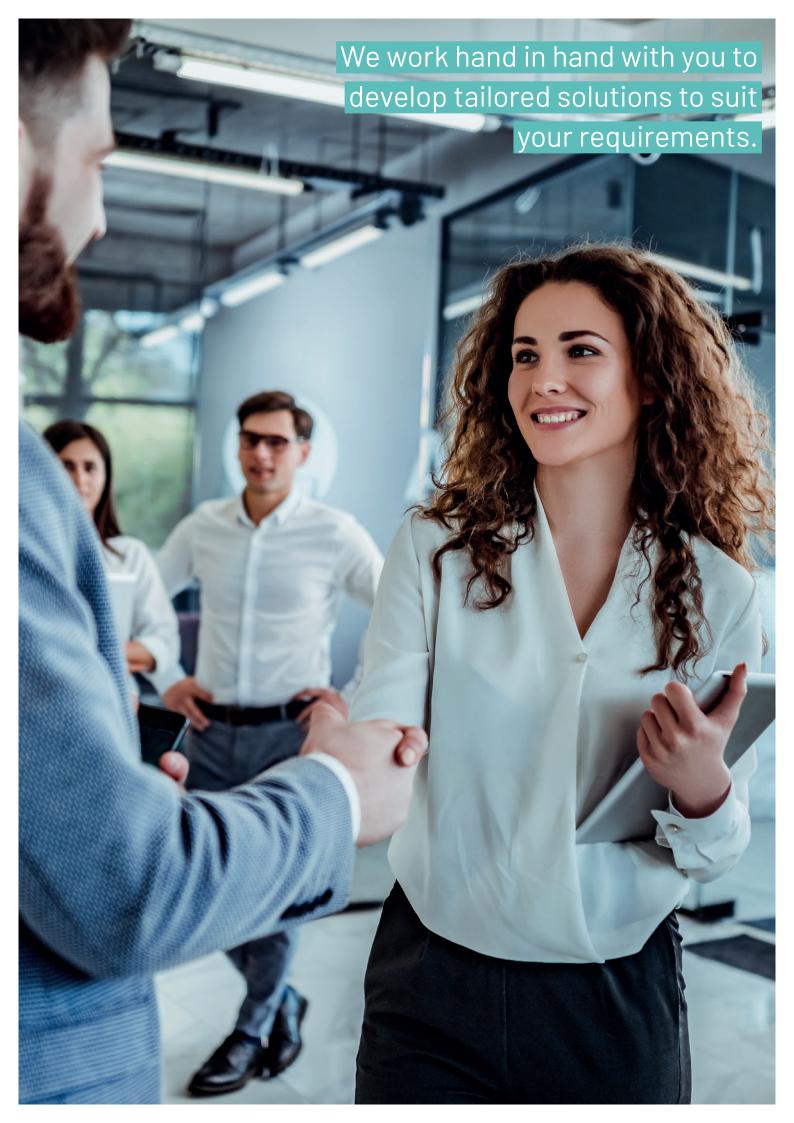

CUSTOMER FOCUS

We focus on you and your goals

WE SUPPORT YOU SO WE CAN MAKE BIG THINGS HAPPEN TOGETHER

Everything we do and the way we do it is designed to make your work easier. By understanding your wishes and needs, as well as your requirements and workflows, we are able to offer you real added value and thus trust in longterm and sustainable customer relationships.

We are here to provide you with the best products and solutions to suit your needs. Our dedicated team ensures that processes run smoothly and is always on hand when needed. Our multidisciplinary teams have many years of practical experience in science and laboratory work and understand the challenges of your daily work first hand. We are happy to use this expert knowledge to create training materials tailored to your specific needs.



WE OFFER:

- / Long-term customer relationship
- / Solutions to suit your requirements
- / Training and training materials







EXPERTISE

In all relevant market trends

WE SPEAK YOUR LANGUAGE

We understand your tasks and their workflows with all their requirements. Our experts are at your side with more than 50 years of experience – from the development of products and solutions to support with regulatory issues or individual training.

More than 80% of our sales and application specialists have a background in natural sciences

and practical experience in the laboratory. We would be happy to advise you on site – at the place where your requirements arise. You can rely on our expertise to quickly integrate our products into your processes, because we know and understand your workflow. We are happy to support you in every step of your individual workflow, for example with product training, support documentation and service.



WHAT WE OFFER:

- / Comprehensive portfolio for your workflow needs
- / Experienced team of industry experts
- / Best practices for optimal efficiency



DISCOVER MORE:





EFFICIENCY

We are your partner for increasing efficiency – inside and outside the laboratory

WE INCREASE YOUR EFFICIENCY

In the laboratory and day-to-day work, increasing efficiency is no longer just an option, but an essential requirement. As a strategic cooperation partner for day-to-day routine activities in and outside the laboratory, we support you in overcoming your current and future challenges.

We offer significant added value by supporting the efficiency of your workflows. For example, our digital solutions create a connection between pre-barcoded test tubes and patient data from the laboratory information system. This reliably rules out mix-ups and minimizes sampling errors. Your workflows can thus be measurably optimized, better results can be achieved with minimum time, resources, and work.



YOUR STRATEGIC PARTNER:

- / Increased efficiency
- / Digital solutions
- / Best practice recommendations







SUSTAINABILITY

the principle behind our processes for many years

BECAUSE IT'S REALLY IMPORTANT TO US

Sustainable action is the basic prerequisite for a future worth living in. As a plastics processing company, we strive to develop sustainable solutions that provide answers to the major challenges of our time.

Greiner AG's overarching "Blue-Plan" strategy focuses on the three pillars of climate, circular economy and people. Greiner Bio-One is also working intensively on these sub-areas.

As a producer of plastic medical products, it is our declared long-term goal to combine ecological, economic and social aspects. We attach particular importance not only to your excellent supply of high-quality products, but also to energy-efficient and sustainable operation as a plastics producer.



OUR GOALS UNTIL 2030:

- / Climate-neutral company
- / Comprehensive circular approach
- / Preparation of employees for the challenges of the future



DISCOVER MORE:




35	CELL CULTURE	
	Cell Culture Flasks	40
	CELLSTAR® Cell Culture Dishes	47
	CELLSTAR® Cell Culture Multiwell Plates	52
	ThinCert® Cell Culture Inserts for 6, 12 and 24 Well	
	Multiwell Plates	56
	CELLSTAR® Cell Culture Microplates	60
	CELLSTAR® Cell Culture Tubes	68
	EASYstrainer Cell Strainers	70
73	MASS CELL CULTURE	
	CELLMASTER Cell Culture Roller Bottles	74
	CELLdisc Cell Culture Device	77
25	3D CELL CULTURE	
03		
	CELLSTAR® Cell Culture Vessels Cell-Repellent Surface	86
	Magnetic 3D Cell Culture	89
	ThinCert® Cell Culture Inserts for 6, 12 and 24 Well	
	Multiwell Plates	94
99	PRODUCTS FOR MICROSCOPY	
	CELLview Dish Cell Culture Dish with Glass Bottom	104
	CELLview Slide Cell Culture Slide with Glass Bottom	106
	CELLview Plate Cell Culture Plate with Glass Bottom	108
	SCREENSTAR Microplates	110
	SensoPlate Glass Bottom Microplates	112
115	TUBES / MULTIPURPOSE BEAKERS	
	Tubes without closures	116
	Tubes with closures	119



	Separation Tubes Multipurpose Containers / Beakers	124 126
131	CRYOTECHNICS	
	Cryo.s Cryo Tubes Cryo.s with Barcode Cryo.s Biobanking Tubes	132 137 141
149	LIQUID HANDLING	
	Sapphire Pipettes Sapphire Pipette Tips CELLSTAR® Serological Pipettes	150 153 162
167	MICROBIOLOGY	
167	MICROBIOLOGY Dishes / Plates / Others	168
167173		168
	Dishes / Plates / Others HTS-MICROPLATES 96 Well Microplates	178
	Dishes / Plates / Others HTS-MICROPLATES 96 Well Microplates 384 Well Microplates	178 182
	Dishes / Plates / Others HTS-MICROPLATES 96 Well Microplates 384 Well Microplates 1536 Well Microplates	178 182 185
	Dishes / Plates / Others HTS-MICROPLATES 96 Well Microplates 384 Well Microplates	178 182
	Dishes / Plates / Others HTS-MICROPLATES 96 Well Microplates 384 Well Microplates 1536 Well Microplates Polypropylene Storage Plates	178 182 185 187
	HTS-MICROPLATES 96 Well Microplates 384 Well Microplates 1536 Well Microplates Polypropylene Storage Plates Microplates for Compound Storage Non-binding Microplates Streptavidin-coated Microplates	178 182 185 187 190 192 195
	HTS-MICROPLATES 96 Well Microplates 384 Well Microplates 1536 Well Microplates Polypropylene Storage Plates Microplates for Compound Storage Non-binding Microplates Streptavidin-coated Microplates UV-Star® Microplates	178 182 185 187 190 192 195
	HTS-MICROPLATES 96 Well Microplates 384 Well Microplates 1536 Well Microplates Polypropylene Storage Plates Microplates for Compound Storage Non-binding Microplates Streptavidin-coated Microplates	178 182 185 187 190 192 195

207	REACTION TUBES / ANALYSER CUPS	
	Reaction Tubes Analyser Cups	208
	Semi-micro / Macro Cuvette	212
215	IMMUNOLOGY / HLA	
	ELISA Microplates	216
	ELISA Strip Plates	218
	Immuno Tubes	221
	Terasaki Plates	223
227	MOLECULAR BIOLOGY	
	Sapphire PCR Tubes	228
	Sapphire PCR Microplates	232
235	LAB EQUIPMENT	
	Lab Equipment	236
2/1	TECHNICAL ADDENDIN	



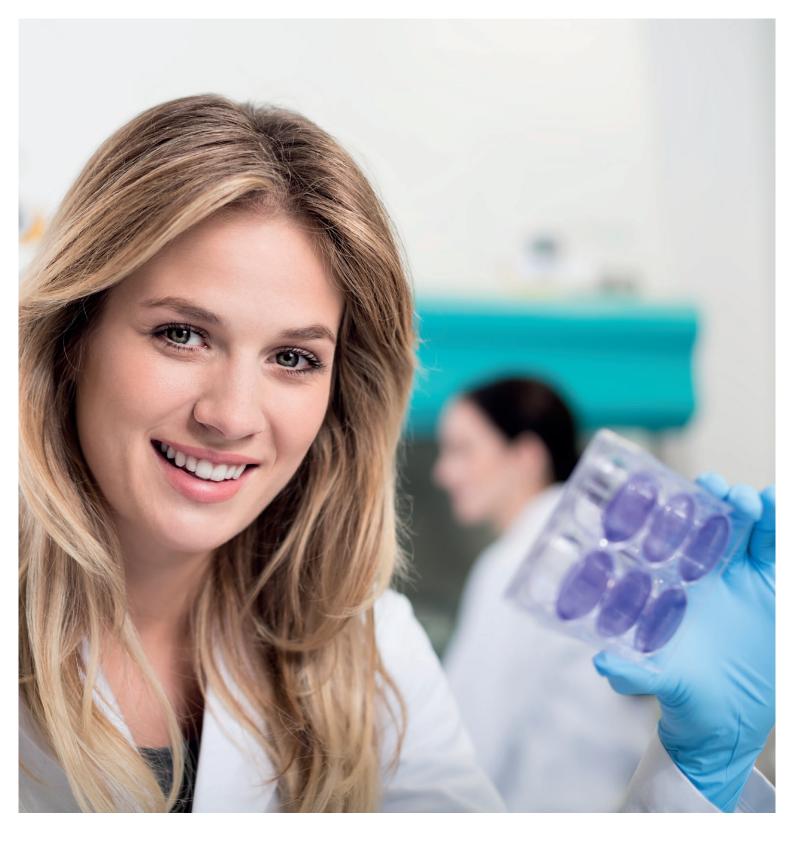

The secret of the continuing success of the CELLSTAR $^{\circ}$ product line from Greiner Bio-One lies in its high quality and reliability. For more than 50 years, Greiner Bio-One has been setting standards for future-oriented research and new technologies in the field of cell culture.

CELL CULTURE


/	Cell Culture Flasks40
	CELLSTAR® Standard Cell Culture Flasks 41
	Cell Culture Flasks - Advanced TC42
	Cell Culture Flasks - CELLCOAT®42
	Suspension Culture Flasks43
	Cell Culture Flasks Cell-Repellent Surface44
	AutoFlask - Cell Culture Flask For Automated
	Systems44
	Cell Scraper45
	Media Bottles45
	Disposal Bags46
/	CELLSTAR® Cell Culture Dishes47
	Cell Culture Dishes48
	Cell Culture Dishes - Advanced TC48
	Cell Culture Dishes - CELLCOAT®49
	Cell Culture Dishes Cell-Repellent Surface50
	CELLview Dish50
	OneWell Plate FourWell Plate51
/	CELLSTAR® Cell Culture
	Multiwell Plates52
	Multiwell Plates 6 / 12 / 24 / 48 Well Format53
	Multiwell Plates Advanced TC54
	Multiwell Plates CELLCOAT®54
	Multiwell Plates Cell-Repellent Surface55
,	ThinCert® Cell Culture Inserts for 6, 12
•	and 24 Well Multiwell Plates
	ThinCert® Cell Culture Inserts 6 Well
	ThinCert® Cell Culture Inserts 12 Well
	ThinCert® Cell Culture Inserts 24 Well58
	ThinCert® Plate 6 / 12 Well
	5016 1 1465 57 12 17611


/	CELLSTAR® Cell Culture Microplates	60
	Cell Culture Microplates 96 Well	6
	Cell Culture Microplates 96 Wel	
	-µClear®	62
	Cell Culture Microplates 96 Well	
	- Half Area	62
	Cell Culture Microplates 96 Wel	
	- CELLCOAT®	63
	Cell Culture Microplates 96 Well	
	- Advanced TC	63
	Suspension Culture Microplates 96 Well	64
	Cell Culture Microplates 384 Well	64
	Cell Culture Microplates 384 Well	
	- Small Volume	65
	Cell Culture Microplates 1536 Well	
	Lids	66
/	CELLSTAR® Cell Culture Tubes	
	CELLreactor	
	CELLSTAR® Cell Culture Tubes	69
/	EASYstrainer Cell Strainers	70
	EASYstrainer	


CELLSTAR®

Quality and reliability for over 50 years

- / Automated production processes
- / Consistent quality control system
- / High-quality raw material

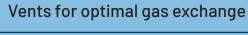

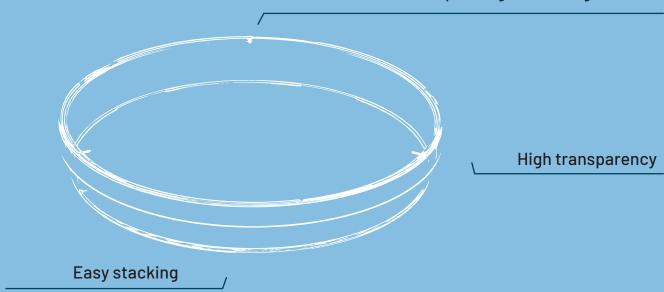

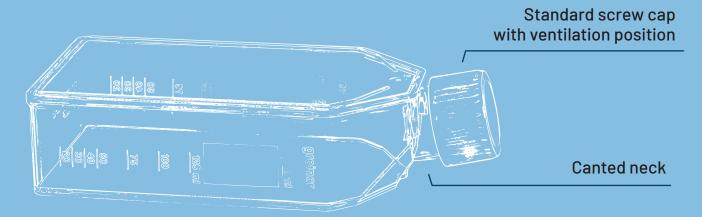

Graduation and writing area on both sides




OPTIMISED CELL CULTURE SURFACES FOR BETTER RESULTS

Cells cultured on plastic surfaces require interaction with molecules in the microenvironment to stimulate normal functioning and proliferation. Easily culturable cell lines are commonly grown on tissue culture treated surfaces like CELLSTAR® tissue culture or in suspension like CELLSTAR® suspension culture. The CELLSTAR® product line offers solutions for most applications in cell biology including the propagation of cells, the performance of cell-based assays, and imaging procedures.

CELLSTAR® cell culture vessels with a cell-repellent surface reliably prevent cell attachment in suspension cultures of semi-adherent and adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient. All cell culture products are labelled with expiry date and lot number in order to ensure transparency of product processes and retraceability of our products throughout the production process.

LITERATURE:

- / Application Note "siRNA-dependent gene silencing on various cell culture surfaces" (F071105)
- / Application Note "Improved cultivation / differentiation of embryonic stem cells" (F073117)
- / Application Note "Cultivation and differentiation of hADSCs with CELLSTAR® and CELLCOAT® products" (F073113)

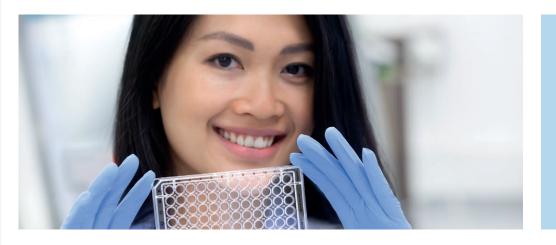

QUALITY CONTROL
Based on automated
production processes
with quality controls,
we provide immaculate,
high-quality products
for all areas of cell
culture.

THE BEST SCREW CAP FOR EVERY APPLICATION

The secured click-in ventilation position is reached, once the cap snaps in audibly. In addition, the correct position is indicated by a vertical tip of a triangle and the readable writing VENT. This allows visual verification of the aeration position, even when the flasks are stacked in the incubator. If the cap is turned clockwise to the end stop, the flask is closed gas-tight.

Filter screw caps for cell culture / suspension culture flasks, roller bottles and the CELLreactor have a patented hydrophobic capillary pore membrane. The defined and constant pore size of 0.2 µm is achieved with minimal variation by means of a specially developed, high-technology method.

The filter insert provides both optimal protection against contamination and efficient gas exchange. By using PET / PTFE which are responsible for the mechanical strength and hydrophobic properties of the membrane, these advantages are retained even if the inside of the cap is briefly wetted with medium.






HIGH-QUALITY RAW MATERIAL

Exclusively high-grade polystyrene and polyethylene terephthalate are used as raw materials for manufacturing our cell culture products.


Polystyrene (PS) is characterised by its high clarity, which greatly simplifies the optical control of cell growth in polystyrene flasks, tubes and roller bottles.

Polyethylene terephthalate (PET) is used for manufacturing roller bottles, media bottles and membranes, due to its beneficial chemical, optical and mechanical properties.

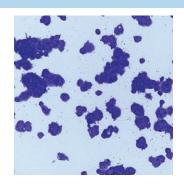

CELLSTAR®


CELLSTAR® TC is the standard surface for cultivation of adherent cells. The special physical surface treatment leads to the incorporation of polar groups such a carboxyl and hydroxyl residues, which functionalised the hydrophobic polystyrene surface resulting in improved and consistent cell attachment.

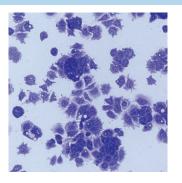

Advanced TC

The Advanced TC surface is based on polymer modification to result in a highly hydrophilic surface which positively influences cellular features and functions.

Advanced TC facilitates consistent and homogenous cell attachment, increasing the overall cell yield and reducing cell loss.

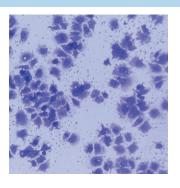

CELLCOAT®

The CELLCOAT® product line comprises cell culture vessels which are coated with biological or synthetic proteins of the extracellular matrix (Collagen Type I, Fibronectin, Laminin) or synthetic proteins (Poly-D-Lysine, Poly-L-Lysine).





/ This video shows how much fun you can have with our cell culture flasks in your lab:



CELL CULTURE FLASKS

Greiner Bio-One offers standard and filter cap cell culture flasks with different surfaces. All Greiner Bio-One cell culture flasks are made of high-grade polystyrene and are free of detectable DNases, RNases, human DNA, endotoxins and are non-cytotoxic.

The specific design facilitates an optimal access of the flask with a cell scraper or pipette. The stacking rim on the top of the flasks ensures firm standing and easy stackability in the incubator. Both sides have a printed graduation for easier

filling. In addition to the filter screw cap Greiner Bio-One offers a standard screw cap with secured click-in ventilation position which guarantees the gas exchange within the flasks.

For the maintenance of adherent cells Greiner Bio-One offers a specific physical surface treatment which ensures adhesion and proliferation of these cells whereas the novel Advanced TC surface provides optimal conditions for the cultivation of sensitive and fastidious cells or the usage of restricted growth conditions.

- / Standard or filter screw cap
- / Canted neck
- / Sterile and user-friendly packaging
- / Growth area for adherent cells: 25, 75 and 175 cm²
- / Filling volume for suspension cells: 50, 250, 550, 650 ml



CELLSTAR® Standard Cell Culture Flasks

For the maintenance of adherent cells Greiner Bio-One offers a specific physical surface treatment which ensures adhesion and proliferation of these cells.

/ Item No. 690170 with measuring grid

STERILE

FREE OF letectable DNase FREE OF detectable numan DNA

FREE OF detectable RNase





Raw material: PS, Surface treatment: TC

Item no.	Flask design	Growth area	Cap colour	Working volume	Total volume	Cap design	Triple -packed	Sterile	Oty. inner / outer
690160		25 cm ²	● red	5 ml - 10 ml	50 ml	screw cap		+	10 / 200
690170		25 cm ²	● red	5 ml - 10 ml	50 ml	screw cap		+	10 / 200
690160-TRI		25 cm ²	● red	5 ml - 10 ml	50 ml	screw cap	yes	+	10 / 130
690175		25 cm ²	● red	5 ml - 10 ml	50 ml	filter screw cap		+	10 / 200
690175-TRI		25 cm ²	● red	5 ml - 10 ml	50 ml	filter screw cap	yes	+	10 / 130
658170		75 cm ²	● red	15 ml - 38 ml	250 ml	screw cap		+	5 / 120
658170-TRI		75 cm ²	● red	15 ml - 38 ml	250 ml	screw cap	yes	+	5/90
658175		75 cm ²	● red	15 ml - 38 ml	250 ml	filter screw cap		+	5 / 120
658175-TRI		75 cm ²	● red	15 ml - 38 ml	250 ml	filter screw cap	yes	+	5/90
660160	flat	175 cm ²	● red	20 ml - 45 ml	550 ml	screw cap		+	5/50
660160-TRI	flat	175 cm²	● red	20 ml - 45 ml	550 ml	screw cap	yes	+	5 / 40
660175	flat	175 cm²	● red	20 ml - 45 ml	550 ml	filter screw cap		+	5/50
660175-TRI	flat	175 cm²	● red	20 ml - 45 ml	550 ml	screw cap	yes	+	5 / 40
661160	high	175 cm²	● red	20 ml - 85 ml	650 ml	screw cap		+	4/40
661175	high	175 cm²	● red	20 ml - 85 ml	650 ml	filter screw cap		+	4/40

[/] The item no. of triple-packed products is composed of the standard item no. plus -TRI (e.g. 604160-TRI). In general, all Greiner Bio-One cell culture products can be produced triple-packed. Please contact your Greiner Bio-One sales representative regarding customised triple-packed products.



Cell Culture Flasks - Advanced TC

The Advanced TC surface provides optimal conditions for the cultivation of sensitive and fastidious cells or the usage of restricted growth conditions.

- / Consistent and even cell attachment
- / Homogeneous and optimised cell growth














Raw material: PS, Surface treatment: Advanced TC

Item no.	Flask design	Growth area	Cap colour	Working volume	Total volume	Cap design	Sterile	Qty. inner / outer
690960		25 cm ²	blue	5 ml - 10 ml	50 ml	screw cap	+	10 / 200
690975		25 cm²	blue	5 ml - 10 ml	50 ml	filter screw cap	+	10 / 200
658970		75 cm ²	blue	15 ml - 38 ml	250 ml	screw cap	+	5 / 120
658975		75 cm²	blue	15 ml - 38 ml	250 ml	filter screw cap	+	5 / 120
660960	flat	175 cm ²	blue	20 ml - 45 ml	550 ml	screw cap	+	5/50
660975	flat	175 cm²	blue	20 ml - 45 ml	550 ml	filter screw cap	+	5/50

- / Forum No. 12: Advanced TC: An innovative surface improving cellular assays (F071104)
- / Application Report "Advanced TC for improving the cultivation / differentiation of embryonic stem cells" (F076036)

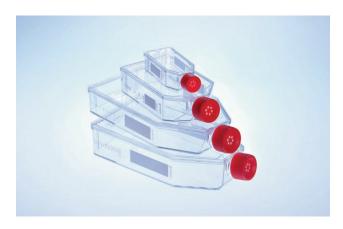

Cell Culture Flasks - CELLCOAT®

- / Canted neck for optimal access
- / Graduation on both sides
- / Improved adhesion and cell proliferation
- / Reduced-serum or serum-free cultivation
- / Improved growth of primary cells

Raw material: PS, Surface treatment: CELLCOAT®, Cap design: filter screw cap

Item no.	Flask design	Growth area	Protein coating	Cap colour	Working volume	Total volume	Qty. inner / outer
690950		25 cm²	Collagen Type I	● red	5 ml - 10 ml	50 ml	10 / 50
658950		75 cm ²	Collagen Type I	● red	15 ml - 38 ml	250 ml	5/50
661950	high	175 cm²	Collagen Type I	● red	20 ml - 85 ml	650 ml	5 / 40

Item no.	Flask design	Growth area	Protein coating	Cap colour	Working volume	Total volume	Qty. inner / outer
690940		25 cm ²	Poly-D-Lysine	● red	5 ml - 10 ml	50 ml	10 / 50
658940		75 cm ²	Poly-D-Lysine	● red	15 ml - 38 ml	250 ml	5/50
661940	high	175 cm ²	Poly-D-Lysine	● red	20 ml - 85 ml	650 ml	5/40
690920		25 cm ²	Fibronectin	● red	5 ml - 10 ml	50 ml	10 / 10
658920		75 cm ²	Fibronectin	● red	15 ml - 38 ml	250 ml	10 / 10
661920	high	175 cm ²	Fibronectin	● red	20 ml - 85 ml	650 ml	5/5
690910		25 cm ²	Laminin	● red	5 ml - 10 ml	50 ml	10 / 10
658910		75 cm ²	Laminin	● red	15 ml - 38 ml	250 ml	10 / 10
661910	high	175 cm ²	Laminin	● red	20 ml - 85 ml	650 ml	5/5



Suspension Culture Flasks

- / Hydrophobic surface, ideal for suspension cultures, hybridoma and embryonic stem cells
- / Graduation on both sides
- / Sterile and user-friendly packaging
- / Canted neck for optimal access

nonpyrogenic

Raw material: PS, Surface treatment: suspension

Item no.	Flask design	Cap colour	Total volume	Cap design	Triple-packed	Sterile	Qty. inner / outer
690190		○ white	50 ml	screw cap		+	10 / 200
690190-TRI		○ white	50 ml	screw cap	yes	+	10 / 130
690195		○white	50 ml	filter screw cap		+	10 / 200
690195-TRI		○ white	50 ml	filter screw cap	yes	+	10 / 130
658190		○white	250 ml	screw cap		+	5 / 120
658190-TRI		○ white	250 ml	screw cap	yes	+	5/90
658195		○white	250 ml	filter screw cap		+	5 / 120
658195-TRI		○ white	250 ml	filter screw cap	yes	+	5/90
660190	flat	○white	550 ml	screw cap		+	5/50
660190-TRI	flat	○ white	550 ml	screw cap	yes	+	5/40
661190	high	○white	650 ml	screw cap		+	4 / 40
661195	high	○ white	650 ml	filter screw cap		+	4 / 40
661195-TRI	high	○white	650 ml	filter screw cap	yes	+	4 / 28



Cell Culture Flasks Cell-Repellent Surface

A cell-repellent surface reliably prevent cell attachment in suspension cultures of semiadherent and adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.




STERILE



Raw material: PS, Surface treatment: cell-repellent

Item no.	Flask design	Cap colour	Total volume	Cap design	Sterile	Qty. inner / outer
690980	flat	○ white	50 ml	screw cap	+	10 / 20
690985	flat	○ white	50 ml	filter screw cap	+	10 / 20
658980	flat	○white	250 ml	screw cap	+	5 / 15
658985	flat	○ white	250 ml	filter screw cap	+	5 / 15
660980	flat	○white	550 ml	screw cap	+	5/5
660985	flat	○ white	550 ml	filter screw cap	+	5/5
661980	high	○white	650 ml	screw cap	+	4/4
661985	high	○white	650 ml	filter screw cap	+	4/4

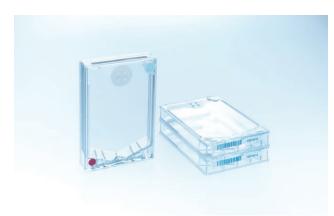

AutoFlask - Cell Culture Flask For Automated Systems


- Standard microplate footprint
- Compatible with a wide range of cell culture and liquid handling systems
- Hydrophobic filter membrane
- Different surface treatments
- Customised barcoding on request



Flask design: AutoFlask, Growth area: 83.6 cm², Raw material: PS, Total volume: 110 ml

Item no.	Barcode	Surface treatment	Product colour	Colour code	Working volume	Sterile	Qty. inner / outer
779160	yes	TC	○ clear	●red	20 ml - 40 ml	+	10 / 100
779190	yes	suspension	○ clear	○white	60 ml - 80 ml	+	10 / 100

[/] Further information on the AutoFlask: "Comparative cell growth study using the AutoFlask" (F072094)

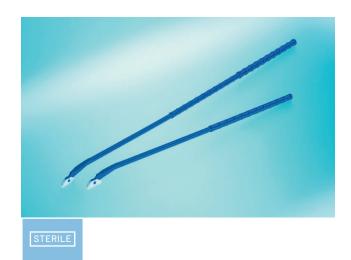

Cell Scraper

- / For gentle mechanical removal of adherent cells
- / Optimised blade design for maximum cell harvest
- / Blade length: 1.8 cm
- / Pivot angle 60°
- / Sterile individual packaging

Description: Cell Scraper, Feature: blade length 1.8 cm

Item no.	Length	Sterile	Qty. inner / outer
541070	28 cm	+	50 / 100
541080	40 cm	+	100 / 100



Media Bottles

- / Made of polyethylene terephthalate (PET)
- / Available in three sizes
- / With graduation
- / Triple-packed for GMP compliant workflow



Flask design: tetragonal, Raw material: PET, Cap design: screw cap

Item no.	Cap colour	Total volume	Sterile	Qty. inner / outer
950700	○white	500 ml	+	50 / 50
951700	○white	100 ml	+	25 / 100
952700	○white	1,000 ml	+	24 / 24



Disposal Bags

Disposal bags made of polypropylene for sterilisation in the steam autoclave. On request they are also available with imprint "Biohazard" if a sufficient number is ordered. For users of hot-air sterilisers, disposal bags made of polyamide are suitable for use up to +160 °C.


Foil thickness: 0.05 mm

Item no.	Length	Width	Nominal capacity	Suitable for steam autoclaves	Suitable for hot air sterilizers	Material	Qty. inner / outer
643201	500 mm	300 mm	10 I	yes		PP	500/500
643401	500 mm	300 mm	10 I		yes	PA	500/500
644201	780 mm	400 mm	301	yes		PP	500/500
644401	780 mm	400 mm	301		yes	PA	500/500
646201	780 mm	600 mm	65 I	yes		PP	500/500
646401	780 mm	600 mm	65 I		yes	PA	300 / 300
649201	1,100 mm	700 mm	130 I	yes		PP	350 / 350
649401	1,100 mm	700 mm	1301		yes	PA	200 / 200



/ For exact dimensions of our cell culture dishes, please refer to the product data sheets on our website.



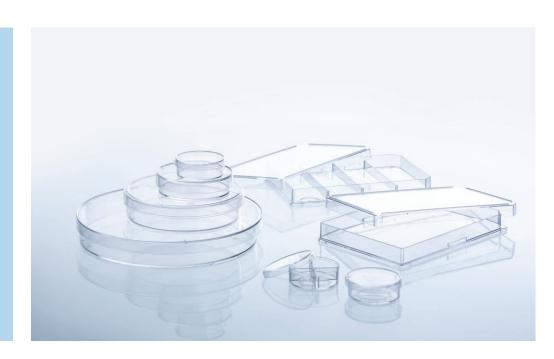

CELLSTAR® CELL CULTURE DISHES

Our CELLSTAR® cell culture dishes are manufactured to meet the highest quality standards at Greiner Bio-One.

This assortment of dishes offers a range of growth sizes from 8.7 to 143 cm². These dimensions are provided in nominal sizes of 35, 60, 100, and 145 mm. For culture dishes associated with growth areas of either 58 or 143 cm², we supply variants with side heights of 20 mm (for special applications). In addition to our standard formats we offer 35 mm cell culture dishes that are equipped with four in-

ner rings. Our standard cell culture dishes have been treated in order to optimise growth and adhesion.

However, we are also capable of providing sterile cell culture trays without any type of coating for suspension culture. Similar to all of our cell culture laboratory products, this range of dishes is packed user-friendly within sterile conditions. This allows technicians to safely handle their samples with a wide variety of cell culture methods without the risk of cell culture contamination.

- / Easy stacking
- Maximal transparency for excellent microscopic analysis
- / Improved cell adhesion through physical surface treatment



Cell Culture Dishes

- / Improved cell adhesion through physical surface treatment
- / Maximal transparency for excellent microscopic analysis
- / Sterile and user-friendly packaging
- / Easy stacking











Surface treatment: TC, Vent nock: yes

Item no.	Compart- ments	Height	Ø nominal size	Growth area	Growth area/unit	Working volume	Total volume	Triple- packed	Sterile	Oty. inner / outer
627160		10 mm	35 mm	8.7 cm ²		≥3 ml	10 ml		+	10 / 740
627170	4	10 mm	35 mm		$0.93\mathrm{cm}^2$	≥0.08 ml	9 ml		+	10 / 740
628160		15 mm	60 mm	21 cm ²		6 ml - 7 ml	28 ml		+	10 / 600
628160-TRI		15 mm	60 mm	21 cm ²		6 ml - 7 ml	28 ml	yes	+	10/300
664160		20 mm	100 mm	58 cm²		16 ml - 17 ml	100 ml		+	15 / 360
664160-TRI		20 mm	100 mm	58 cm ²		16 ml - 17 ml	100 ml	yes	+	15 / 180
639160		20 mm	145 mm	143 cm²		25 ml - 27 ml	240 ml		+	5 / 120
639160-TRI		20 mm	145 mm	143 cm ²		25 ml - 27 ml	240 ml	yes	+	5 / 120

 ${\it I} \quad \hbox{For exact dimensions of our cell culture dishes, please refer to the product data sheets on our website.}$



Cell Culture Dishes - Advanced TC

The Advanced TC surface provides optimal conditions for the cultivation of sensitive and fastidious cells or the usage of restricted growth conditions.

- / Consistent and even cell attachment
- / Homogeneous and optimised cell growth

STERILE FREE OF detectable detectable human DNA RN





Surface treatment: Advanced TC, Vent nock: yes

Item no.	Height	Ø nominal size	Growth area	Working volume	Total volume	Sterile	Qty. inner / outer
627960	10 mm	35 mm	$8.7\mathrm{cm}^2$	≤3 ml	10 ml	+	10 / 740
628960	15 mm	60 mm	21 cm ²	6 ml - 7 ml	28 ml	+	10 / 600

Item no.	Height	Ø nominal size	Growth area	Working volume	Total volume	Sterile	Qty. inner / outer
664960	20 mm	100 mm	58 cm²	16 ml - 17 ml	100 ml	+	15 / 360
639960	20 mm	145 mm	143 cm ²	25 ml - 27 ml	240 ml	+	5 / 120

- / Forum No. 12: Advanced TC: An innovative surface improving cellular assays (F071104)
- $\textit{I} \quad \text{Application Report "Advanced TC for improving the cultivation / differentiation of embryonic stem cells" (F076036) } \\$



Cell Culture Dishes - CELLCOAT®

- / Improved adhesion and cell proliferation
- / Reduced-serum or serum-free cultivation
- / Improved growth of primary cells

Surface treatment: CELLCOAT®, Vent nock: yes

Item no.	Height	Ø nominal size	Growth area	Protein coating	Working volume	Total volume	Qty. inner / outer
628950	15 mm	60 mm	21 cm ²	Collagen Type I	6 ml - 7 ml	28 ml	20 / 100
664950	20 mm	100 mm	58 cm ²	Collagen Type I	16 ml - 17 ml	100 ml	10 / 40
628940	15 mm	60 mm	21 cm ²	Poly-D-Lysine	6 ml - 7 ml	28 ml	20 / 100
664940	20 mm	100 mm	58 cm ²	Poly-D-Lysine	16 ml - 17 ml	100 ml	10 / 40
628930	15 mm	60 mm	21 cm ²	Poly-L-Lysine	6 ml - 7 ml	28 ml	20 / 100
628920	15 mm	60 mm	21 cm ²	Fibronectin	6 ml - 7 ml	28 ml	5/20
664920	20 mm	100 mm	58 cm²	Fibronectin	16 ml - 17 ml	100 ml	5 / 10
628910	15 mm	60 mm	21 cm ²	Laminin	6 ml - 7 ml	28 ml	5/20
664910	20 mm	100 mm	58 cm ²	Laminin	16 ml - 17 ml	100 ml	10 / 10



Cell Culture Dishes Cell-Repellent Surface

Cell culture vessels with cell-repellent surface reliably prevent cell attachment in suspension cultures of semi adherent / adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.










Raw material: PS, Surface treatment: cell-repellent, Vent nock: yes

Item no.	Height	Ø nominal size	Working volume	Total volume	Sterile	Qty. inner / outer
627979	10 mm	35 mm	≤3 mI	10 ml	+	10 / 40
628979	15 mm	60 mm	6 ml - 7 ml	28 ml	+	10 / 20
664970	20 mm	100 mm	16 ml - 17 ml	100 ml	+	1/5

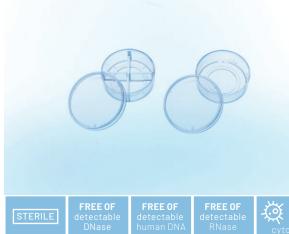

CELLview Dish

The CELLview cell culture dish combines the convenience of a standard size 35 mm disposable plastic cell culture dish with the optical quality of glass, providing superior high-resolution microscopic images of invitro cultivated cultures.

/ Embedded glass bottom for maximal planarity





Height: 10 mm, Ø: 35 mm, Bottom: glass, Vent nock: yes

Item no.	Compart- ments	Growth area	Growth area/unit	Surface treatment	Working volume	Total volume	Total volume (well)	Working volume (well)	Sterile	Oty. inner / outer
627861	1	8.7 cm ²		untreated	2.5 ml - 5 ml	10 ml			+	10 / 40
627860	1	8.7 cm ²		TC	2.5 ml - 5 ml	10 ml			+	10 / 40
627965	1	8.7 cm ²		Advanced TC	2.5 ml - 5 ml	10 ml			+	10 / 40
627871	4		1.9 cm ²	untreated			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40

Item no.	Compart- ments	Growth area	Growth area / unit	Surface treatment	Working volume	Total volume	Total volume (well)	Working volume (well)	Sterile	Oty. inner / outer
627870	4		1.9 cm ²	TC			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40
627975	4		1.9 cm ²	Advanced TC			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40

- / Application Note "Protein localisation using confocal laser scanning microscopy" (F073101)
- $\textit{I} \quad \text{Application Note } \textit{``Live cell imaging on Golgi morpholyy using the CELL view dish"} (F074048)$



OneWell Plate FourWell Plate

- / For the cultivation of a large number of cells the OneWell Plate with ANSI standard dimensions offers the best possible conditions.
- / The FourWell Plates allows users to culture cells on slides while remaining in accordance with an HTS-compatible plate. This four-chambered plate is particularly suitable for parallel experiments.



Height: 14.4 mm, Length: 127.8 mm, Width: 85.5 mm, Lid: yes, Vent nock: yes

Item no.	Well format	Growth area	Surface treatment	Product colour	Total volume (well)	Sterile	Qty. inner / outer
96077307	4		suspension	○ clear	18.6 ml	+	8 / 32
670190	1	95 cm²	suspension	○ clear	113.7 ml	+	8 / 32
670180	1	95 cm²	TC	○ clear	113.7 ml	+	8 / 32

/ Sample packs are available on request.





/ Watch our video:



CELLSTAR®


CELL CULTURE MULTIWELL PLATES

For both adherent and suspension cell culture, choose Greiner Bio-One's multiwell plates in either 6, 12, 24 or 48 well format.

The special chimney wells in conjunction with low evaporation lids significantly reduce the risk of cross-contamination between the wells. The plates are easily stacked, both with and without lids. Due to their ad-

vanced design, the plates still exhibit important features like constant heat and gas exchange even when stacked. Uniform plate dimensions (well depth, diameter and distance between the wells) comply with ANSI standards, enabling easy manual and automated handling. Each plate is sterile, free of detectable DNase, RNase, human DNA, is non-pyrogenic and has a printed lot number for traceability.

- / With hydrophilic surface (TC surface treatment) for improved cell adhesion
- With hydrophobic surface for suspension cultures and hybridoma cells
- / High clarity and low autofluorescence
- / Alphanumeric well coding

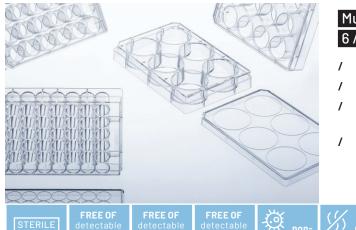


Multiwell Plates 6 / 12 / 24 / 48 Well Format

- / High transparency
- / Alphanumeric well coding
- / With hydrophilic surface (TC surface treatment) for improved cell adhesion
- / With hydrophobic surface for suspension cultures and hybridoma cells

Bottom: solid, Lid: yes, condensation rings

Item no.	Well format	Growth area / unit	Raw material	Surface treatment	Product colour	Total volume (well)	Working volume (well)	Sterile	Qty. inner / outer
657160	6	9.6 cm ²		TC	○ clear	16 mI	2 ml - 5 ml	+	1 / 100
665180	12	3.9 cm ²	PS	TC	○ clear	6.5 ml	2 ml - 4 ml	+	1/100
662160	24	1.9 cm ²		TC	○ clear	3.3 ml	0.5 ml - 1.5 ml	+	1 / 100
677180	48	1 cm²		TC	○ clear	1.7 ml	0.5 ml - 1 ml	+	1/100
657185	6			suspension	○ clear	16 mI		+	1 / 100
665102	12		PS	suspension	○ clear	6.5 ml		+	1 / 100
662102	24			suspension	○ clear	3.3 ml		+	1 / 100
677102	48			suspension	○ clear	1.7 ml		+	1 / 100

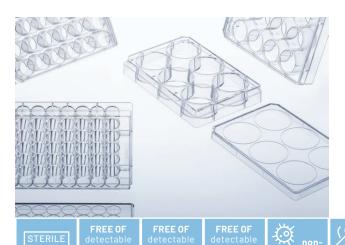

Multiwell Plates Advanced TC

The Advanced TC surface provides optimal conditions for the cultivation of sensitive and fastidious cells or the usage of restricted growth conditions.

/ Consistent and even cell attachment



Item no.	Well format	Growth area / unit	Raw material	Product colour	Total volume (well)	Working volume (well)	Sterile	Qty. inner / outer
657960	6	$9.6\mathrm{cm}^2$		○ clear	16 ml	2 ml - 5 ml	+	1 / 100
665980	12	$3.9\mathrm{cm}^2$	PS	○ clear	6.5 ml	2 ml - 4 ml	+	1/100
662960	24	1.9 cm ²		○ clear	3.3 ml	0.5 ml - 1.5 ml	+	1 / 100
677980	48	1 cm²		○ clear	1.7 ml	0.5 ml - 1 ml	+	1 / 100

- / Forum No. 12: Advanced TC: An innovative surface improving cellular assays (F071104)
- / Application Report "Advanced TC for improving the cultivation / differentiation of embryonic stem cells" (F076036)



Multiwell Plates CELLCOAT®

- / Improved adhesion and cell proliferation
- / Reduced-serum or serum-free cultivation
- / Improved growth of primary cells

Bottom: solid, Surface treatment: CELLCOAT®, Lid: yes, condensation rings

Item no.	Well format	Growth area / unit	Protein coating	Product colour	Total volume (well)	Working volume (well)	Oty. inner / outer
657950	6	9.6 cm²	Collagen Type I	○ clear	16.1 ml	2 ml - 5 ml	5/50
662950	24	1.9 cm²	Collagen Type I	○ clear	3.3 ml	0.5 ml - 1 ml	5/50
657940	6	9.6 cm²	Poly-D-Lysine	○ clear	16.1 ml	2 ml - 5 ml	5/50
662940	24	1.9 cm²	Poly-D-Lysine	○ clear	3.3 ml	0.5 ml - 1 ml	5/50
657930	6	9.6 cm²	Poly-L-Lysine	○ clear	16.1 ml	2 ml - 5 ml	5/50
662930	24	1.9 cm²	Poly-L-Lysine	○ clear	3.3 ml	0.5 ml - 1 ml	5/50



Multiwell Plates Cell-Repellent Surface

Cell culture vessels with cell-repellent surface reliably prevent cell attachment in suspension cultures of semi adherent / adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non- cytotoxic	y py
---------	--------------------------------	------------------------------------	--------------------------------	-------------------	---------

Well profile: F-bottom, Bottom: solid, Raw material: PS, Surface treatment: cell-repellent, Lid: yes, condensation rings

Item no.	Well format	Product colour	Total volume (well)	Working volume (well)	Sterile	Qty. inner / outer
657970	6	○ clear	16.1 ml	2 ml - 5 ml	+	1/5
662970	24	○ clear	3.3 ml	0.5 ml - 1.5 ml	+	1/5
677970	48	○ clear	1.7 ml	0.5 ml - 1 ml	+	1/5

[/] Forum No. 17: CELLSTAR® Cell Culture Vessels with Cell-Repellent Surface (F073777)

 $[\]textit{Application Report "Advantage of CELLSTAR"} \textit{ Cell Culture Vessels with Cell-Repellent Surfaces for 3-D Cell Culture in Hydrogels" (F073792). } \\$





- / Forum No. 8: ThinCert° cell culture products Overview (F073017)
- / Application Note "Immuno cytochemistry"(F073100)
- / For further information and application notes please refer to our website: www.gbo.com

THINCERT® CELL CULTURE INSERTS FOR 6, 12 AND 24 WELL MULTIWELL PLATES

For advanced cell and tissue culture applications,
Greiner Bio-One offers a broad range of membrane supports
- ThinCert®.

Combining 6 different membrane specifications (pore size and density) in geometries to fit 6, 12 and 24 well plates, the ThinCert® cell culture inserts are suitable for a wide range of applications including transport, secretion and diffusion studies, migrational experiments, cytotoxicity testing, co-cultures, trans epithelial electric resist-

ance (TEER) measurements, as well as primary cell culture. ThinCert® cell culture inserts are compatible with standard CELLSTAR® multiwell plates from Greiner Bio-One, and are pre-packed together with the requisite number of plates. The automated production process includes double optical control of each insert produced, ensuring that any biological contamination is avoided. The sterility of the single blisterpacked inserts and multiwell plates is ensured by irradiation.

- / Stable clear polystyrene housing
- / Sealed PET capillary pore membrane
- / Pre-configured multiwell plates with ThinCert® cell culture inserts available on request



ThinCert® Cell Culture Inserts 6 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange



Feature: 4 multiwell plates / box, Height: 16.25 mm, Ø internal: 24.85 mm, Ø external: 27.85 mm, Cultural surface: 452.4 mm², Surface treatment: TC, Working volume (ThinCert®): 1 ml - 3.6 ml, Working volume (well): 2 ml - 4.15 ml

Item no.	Pore density	Ø pore Optical membrane properties		Sterile	Qty. inner / outer
657640	$1 \times 10^{8} \text{cm}^{2}$	0.4 µm	translucent	+	1 / 24
657641	$2 \times 10^{6} \text{ cm}^{2}$	0.4 µm	clear	+	1 / 24
657610	$2 \times 10^{6} \text{ cm}^{2}$	1 µm	clear	+	1 / 24
657630	0,6 x 10 ⁶ cm ²	3 µm	clear	+	1 / 24
657631	$2 \times 10^{6} \text{ cm}^{2}$	3 µm	translucent	+	1 / 24
657638	$0.15 \times 10^{6} \text{cm}^{2}$	8 µm	translucent	+	1 / 24



ThinCert® Cell Culture Inserts 12 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange

Feature: 4 multiwell plates / box, Height: 16.25 mm, Ø internal: 13.85 mm, Ø external: 15.85 mm, Cultural surface: 113.1 mm², Surface treatment: TC, Working volume (ThinCert®): 0.3 ml - 1 ml, Working volume (well): 1 ml - 2 ml

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
665640	1 x 10 ⁸ cm ²	0.4 µm	translucent	+	1 / 48
665641	$2 \times 10^{6} \text{cm}^{2}$	0.4 µm	clear	+	1 / 48
665610	2 x 10 ⁶ cm ²	1 µm	clear	+	1 / 48

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Oty. inner / outer
665630	0,6 x 10 ⁶ cm ²	3 µm	clear	+	1 / 48
665631	$2 \times 10^{6} \text{cm}^{2}$	3 µm	translucent	+	1 / 48
665638	0,15 x 10 ⁶ cm ²	8 µm	translucent	+	1 / 48



ThinCert® Cell Culture Inserts 24 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange









Feature: 2 multiwell plates / box, Height: 16.25 mm, Ø internal: 8.4 mm, Ø external: 10.4 mm, Cultural surface: 33.6 mm², Surface treatment: TC, Working volume (ThinCert®): 0.1 ml - 0.35 ml, Working volume (well): 0.4 ml - 1.2 ml

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
662640	$1 \times 10^{8} \text{cm}^{2}$	0.4 µm	translucent	+	1 / 48
662641	$2 \times 10^{6} \text{cm}^{2}$	0.4 µm	clear	+	1 / 48
662610	$2 \times 10^{6} \text{cm}^{2}$	1 µm	clear	+	1 / 48
662630	0,6 x 10 ⁶ cm ²	3 µm	clear	+	1 / 48
662631	$2 \times 10^6 \text{cm}^2$	3 µm	translucent	+	1 / 48
662638	0,15 x 10 ⁶ cm ²	8 µm	translucent	+	1 / 48



ThinCert® Plate 6 / 12 Well

- / Optimised for use with ThinCert® cell culture inserts
- / Deep wells for an increased volume of medium in air-lift culture
- / Notches for fixed insert position
- / Available in 6 and 12 well format













Height: 39.5 mm, Length: 129.5 mm, Width: 86.6 mm, Lid: yes, yes, condensation rings

Item no.	Well format	Well format Working volume (well)		Qty. inner / outer
657110	6	≤20 ml	+	1/50
665110	12	≤4 ml	+	1 / 60





- / Application Note
 "Establishing a cell
 culture assay based on
 TR-FRET for screening
 G-Protein-coupled
 receptors" (F074058)
- / Application Note "Selection of cell culture surfaces for the adipogenic differentiation of hMSCs" (F010003)

CELLSTAR®

CELL CULTURE MICROPLATES

Cell culture treated microplates are available in the following versions: 96, 384, 1536 well format.

Depending on the application, the well profile is a key feature in a 96 well cell culture microplate. The chimney well cell culture microplate has the same well profile as the standard F-bottom plate. The difference to the standard plate is the chimney-like arrangement of the wells i.e. each well stands on its own. Therefore the risk of contamination from sample material being carried over is minimised. Clear-bottom mi-

croplates have pigmented walls and a transparent thin film bottom, the so-called μ Clear® bottom. In contrast to our standard microplates with a solid polystyrene bottom, they are ideal for cell culture and microscopic applications using fluorescence or luminescence detection methods.

For many applications, a reduction of the sample volume is an important feature. 96 well half area microplates offer an interesting alternative here. They can be pipetted automatically as well as manually without any problem and allow a reduction of the sample volume up to 50 %.

- Available with different cell culture surfaces for optimal cell culture conditions
- / Footprint compatible with automated systems
- / Alphanumeric well coding



Cell Culture Microplates 96 Well

- / With U-bottom, V-bottom or F-bottom
- / Clear / black / white
- / Chimney well design, raised wells and yes, condensation rings in lids prevent cross-contamination
- / Improved cell adhesion through physical surface treatment

STERILE

FREE OF detectable DNase FREE OF detectable numan DNA FREE OF detectable RNase





Well format: 96, Bottom: solid, Raw material: PS, Surface treatment: TC

Item no.	Growth area / unit	Well profile	Product colour	Working volume (well)	Lid	Sterile	Oty. inner / outer
650160	35 mm ²	U-bottom	○ clear	40 μΙ - 280 μΙ	no	+	1 / 100
650180	35 mm²	U-bottom	○ clear	40 μl - 280 μl	yes	+	1 / 100
651160	28 mm ²	V-bottom	○ clear	40 μΙ - 200 μΙ	no	+	1 / 100
651180	28 mm²	V-bottom	○ clear	40 μΙ - 200 μΙ	yes	+	1 / 100
655160	34 mm ²	F-bottom / Chimney Well	○ clear	25 µl - 340 µl	no	+	1 / 100
655162	34 mm ²	F-bottom / Chimney Well	○ clear	25 µl - 340 µl	no	+	5 / 100
655180	34 mm²	F-bottom / Chimney Well	○ clear	25 μΙ - 340 μΙ	yes, yes, condensa- tion rings	+	1 / 100
655182	34 mm²	F-bottom / Chimney Well	○clear	25 μΙ - 340 μΙ	yes, yes, condensa- tion rings	+	10 / 160
655073	34 mm ²	F-bottom / Chimney Well	○white	25 µl - 340 µl	no	+	10 / 40
655083	34 mm²	F-bottom / Chimney Well	○white	25 μΙ - 340 μΙ	yes, yes, condensa- tion rings	+	8 / 32
655079	34 mm²	F-bottom / Chimney Well	● black	25 µl - 340 µl	no	+	10 / 40
655086	34 mm²	F-bottom / Chimney Well	● black	25 μΙ - 340 μΙ	yes, yes, condensa- tion rings	+	8 / 32

[/] For selected products Greiner Bio-One also offers user-friendly bulk packaging.

[/] Barcode labelling on request



Cell Culture Microplates 96 Well - µClear®

Clear-bottom microplates have pigmented walls and a transparent thin film bottom, the so-called µClear® bottom. In contrast to our standard microplates with a solid polystyrene bottom, they are ideal for cell culture and microscopic applications using fluorescence or luminescence detection methods.



Well format: 96, Growth area / unit: 34 mm², Well profile: F-bottom / Chimney Well, Bottom: μClear®, Raw material: PS, Surface treatment: TC, Working volume (well): 25 μl - 340 μl

Item no.	Product colour	Lid	Sterile	Qty. inner / outer
655088	○white	no	+	10 / 40
655098	○white	yes, condensation rings	+	8 / 32
655087	● black	no	+	10 / 40
655090	● black	yes, condensation rings	+	8 / 32



Cell Culture Microplates 96 Well - Half Area

For many applications, a reduction of the sample volume is an important feature. 96 well half area microplates offer an interesting alternative here. They can be pipetted automatically as well as manually without any problem and allow a reduction of the sample volume up to 50 %.



Well format: 96, Growth area / unit: 15 mm², Well profile: F-bottom, Raw material: PS, Surface treatment: TC,

Plate design: half area, Working volume (well): 15 µl - 175 µl, Lid: yes

Item no.	Bottom	Product colour	Sterile	Qty. inner / outer
675180	solid	○ clear	+	8 / 32
675083	solid	○white	+	8 / 32
675086	solid	● black	+	8 / 32
675090	µClear®	● black	+	8 / 32

[/] For selected products Greiner Bio-One also offers user-friendly bulk packaging.



Cell Culture Microplates 96 Well - CELLCOAT®

- / Improved adhesion and cell proliferation
- / Reduced-serum or serum-free cultivation
- / Improved growth of primary cells

Well format: 96, Growth area / unit: 34 mm², Well profile: F-bottom / Chimney Well, Raw material: PS, Surface treatment: CELLCOAT®, Working volume (well): 25 µl - 340 µl


Item no.	Bottom	Protein coating	Product colour	Lid	Qty. inner / outer
655950	solid	Collagen Type I	○clear	yes, condensation rings	5/20
655956	µClear®	Collagen Type I	● black	yes, condensation rings	5/20
655940	solid	Poly-D-Lysine	○clear	yes, condensation rings	5/20
655944	µClear®	Poly-D-Lysine	○white	yes, condensation rings	5/20
655946	µClear®	Poly-D-Lysine	● black	yes	5/20
655948	µClear®	Poly-D-Lysine	● black	yes, condensation rings	20 / 120
655930	solid	Poly-L-Lysine	○clear	yes, condensation rings	5/20
655936	µClear®	Poly-L-Lysine	● black	yes, condensation rings	5/20



Cell Culture Microplates 96 Well - Advanced TC

The Advanced TC surface provides optimal conditions for the cultivation of sensitive and fastidious cells or the usage of restricted growth conditions.

/ Consistent and even cell attachment

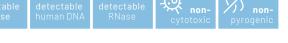

Well format: 96, Growth area / unit: 34 mm², Well profile: F-bottom / Chimney Well, Raw material: PS, Surface treatment: Advanced TC, Working volume (well): 25 μl - 340 μl, Lid: yes, condensation rings

Item no.	Bottom	Product colour	Sterile	Qty. inner / outer
655980	solid	○ clear	+	1 / 100
655983	µClear®	○white	+	8 / 32
655986	µClear®	● black	+	8 / 32



Suspension Culture Microplates 96 Well

- / Hydrophobic surface, ideal for suspension cultures, hybridoma and embryonic stem cells
- / Solid bottom









Well format: 96, Bottom: solid, Raw material: PS, Surface treatment: suspension

Item no.	Well profile	Product colour	Working volume (well)	Lid	Sterile	Qty. inner / outer
650185	U-bottom	○clear	40 µl - 280 µl	yes	+	60 / 240
655185	F-bottom / Chimney Well	○ clear	25 µl - 340 µl	yes, conden- sation rings	+	60 / 240



Cell Culture Microplates 384 Well

- / Clear / black / white
- / Solid bottom or $\mu Clear^{\circ}$ film bottom
- / Barcode labelling on request
- / Alphanumeric well coding

Well format: 384, Growth area / unit: 10 mm², Well profile: F-bottom, Raw material: PS, Working volume (well): 15 µl - 110 µl

Item no.	Bottom	Surface treatment	Protein coating	Product colour	Total volume (well)	Lid	Sterile	Qty. inner / outer
781165	solid	TC		○ clear		no	+	10 / 40
781182	solid	TC		○ clear		yes	+	8 / 32
781073	solid	TC		○white		no	+	10 / 40
781080	solid	TC		○ white		yes	+	8 / 32
781079	solid	TC		● black		no	+	10 / 40
781086	solid	TC		● black		yes	+	8 / 32
781093	μClear®	TC		○white		no	+	10 / 40
781098	μClear®	TC		○ white		yes	+	8 / 32
781092	µClear®	TC		● black		no	+	10 / 40
781091	μClear®	TC		● black		yes	+	8 / 32

Item no.	Bottom	Surface treatment	Protein coating	Product colour	Total volume (well)	Lid	Sterile	Qty. inner / outer
781090	µClear®	TC		● black		yes	+	20 / 120
781950	solid	CELLCOAT®	Collagen Type I	○ clear	131 μΙ	yes		5/20
781956	µClear®	CELLCOAT®	Collagen Type I	● black	131 μΙ	yes		5/20
781940	solid	CELLCOAT®	Poly-D-Lysine	○ clear	131 µI	yes		5/20
781945	solid	CELLCOAT®	Poly-D-Lysine	○white	131 µI	yes		5/20
781944	μClear®	CELLCOAT®	Poly-D-Lysine	○white	131 µI	yes		5/20
781946	µClear®	CELLCOAT®	Poly-D-Lysine	● black	131 µI	yes		5/20
781948	μClear®	CELLCOAT®	Poly-D-Lysine	● black	131 µI	yes		20 / 120
781930	solid	CELLCOAT®	Poly-L-Lysine	○ clear	131 µI	yes		5/20
781936	μClear®	CELLCOAT®	Poly-L-Lysine	● black	131 µI	yes		5/20
781983	µClear®	Advanced TC		○white	131 µI	yes	+	8 / 32
781986	μClear®	Advanced TC		● black	131 µl	yes	+	8 / 32




Cell Culture Microplates 384 Well - Small Volume

HiBase

- / Perfect for top reading even at low working volumes
- / Savings in reagent similar to 1536 well microplates
- / Made of black / white polystyrene for fluorescence or luminescence measurements

Well format: 384, Growth area / unit: 2.7 mm², Well profile: F-bottom, Bottom: solid, Raw material: PS, Plate geometry: HiBase, Plate design: Small Volume, Working volume (well): 4 µl - 25 µl, Lid: yes

Item no.	Surface treatment	Protein coating	Product colour	Sterile	Qty. inner / outer
784080	TC		○white	+	8 / 32
784086	TC		● black	+	8 / 32
784946	CELLCOAT®	Poly-D-Lysine	● black		5/30

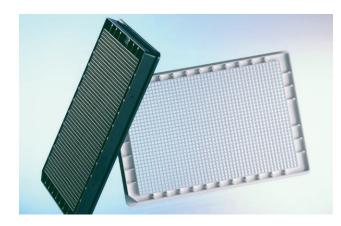


Cell Culture Microplates 1536 Well

- / Clear / black / white
- / Solid bottom or µClear® film bottom
- / Barcode labelling on request

Well format: 1536, Growth area / unit: 2.3 mm², Well profile: F-bottom, Raw material: PS, Plate geometry: HiBase, Working volume (well): 3 µl - 10 µl

Item no.	Bottom	Surface treatment	Protein coating	Product colour	Lid	Sterile	Qty. inner / outer
782180	solid	TC		○ clear	yes	+	1/32
782073	solid	TC		○white	no	+	15 / 60
782080	solid	TC		○white	yes	+	10 / 40
782078	solid	TC		black	no	+	15 / 60
782086	solid	TC		● black	yes	+	10 / 40
782093	μClear®	TC		○white	no	+	15 / 60
782092	µClear®	TC		black	no	+	15 / 60
782946	µClear®	CELLCOAT®	Poly-D-Lysine	black	yes		5/20

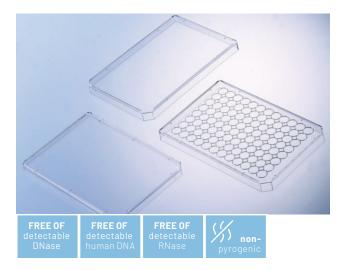

Lids

All sterile lids are non-cytotoxic.

Description: Lid, Raw material: PS

Item no.	Height	Yes, condensation rings	Lid type	Sterile	Qty. inner / outer
656101	9 mm	no	high		1 / 100
656161	9 mm	no	high	+	1 / 100
656170	9 mm	yes	high		1 / 100
656171	9 mm	yes	high	+	1/100

Item no.	Height	Yes, condensation rings	Lid type	Sterile	Qty. inner / outer
656190	6 mm	no	flat		20 / 200
656191	6 mm	no	flat	+	20 / 200
691101	4.8 mm	no	ultra low		25 / 100
691161	4.8 mm	no	ultra low	+	5 / 100





- / Application Note: Cultivation of Suspension and Hybridoma Cells in CELLSTAR® CELLreactor Tubes (F073918)
- / Application Note:
 Superior protein yields
 in suspension CHO cells
 using FectoPRO™-mediated transient transfection in CELLSTAR®
 CELLreactor (F073926)

CELLSTAR® CELL CULTURE TUBES

CELLSTAR® CELLreactor tube can be used as small bioreactor for suspension and spheroid cell culture, facilitating miniaturisation of large-scale setups and maximising the number of parallel experiments. Each CELLreactor tube cap features several holes and a membrane with a pore size of 0.2 µm to guarantee maximal sterility while providing excellent gas exchange. In case the aeration has to be reduced, individual openings can be sealed.

Agitation of internal liquids is achieved with standard shaking lab equipment minimising foam formation and shearing forces induced by integrated mixing devices. Compared to cell culture and spinner flask as well as other cultivation disposables, no transfer for cell harvest is reguired. Based on the conical design, the tubes fit in standard 15 ml / 50 ml centrifuge rotors and cells can be spun down in the same tube. In addition to cell culture applications, the CELLSTAR® CELLreactor tube can also be applied for the expansion of aerobic bacteria, yeast or other microorganisms in shaken cultures as well as storage of components and liquids requiring gas exchange.

- / Bioreactor for suspension and spheroid cells
- Expansion of aerobic bacteria, yeast and microorganisms
- / Storage of components and liquids requiring gas exchange



CELLreactor

15 ml and 50 ml polypropylene tube with filter screw cap

- / For cultivation of suspension cells and expansion of aerobic microorganisms
- / Facilitates a high number of parallel experiments
- / Maximal sterility and excellent gas exchange
- Conical tube design and in-tube harvest

STERILE FREE OF detectable DNase hu

FREE OF FRE detectable LIMB RN.





Graduation: yes, Writing field: yes, Raw material: PP, Bottom shape: conical, Cap design: filter screw cap

Item no.	Height	Ø	Cap colour	Working volume	Nominal volume	Sterile	Qty. inner / outer
188241	120 mm	17 mm	blue	1 ml - 5 ml	15 mI	+	20 / 500
227245	115 mm	30 mm	blue	1 ml - 35 ml	50 ml	+	20 / 500

- $\textit{I} \quad \text{Application Note: Cultivation of Suspension and Hybridoma Cells in CELLSTAR} \text{ } \textit{CELL} \textit{reactor Tubes} (F073918) \\$
- / Application Note: Superior protein yields in suspension CHO cells using FectoPRO™-mediated transient transfection in CELLSTAR® CELLreactor(F073926)




CELLSTAR® Cell Culture Tubes

- / Manufactured from crystal clear polystyrene
- / Improved cell adhesion through physical surface treatment
- / Available with screw cap, bayonet cap or two-position cap

Raw material: PS, Surface treatment: TC

Item no.	Height	Ø	Support skirt	Cap colour	Working volume	Nominal volume	Cap design	Sterile	Oty. inner / outer
120160	75 mm	12.4 mm	no		≤4 ml	4.5 ml		+	1 / 1,000
120190	75 mm	12.4 mm	no		≤4 ml	4.5 ml		+	25 / 2,000
163160	100 mm	17 mm	no	● red	≤12 ml	12 ml	screw cap	+	5 / 1,000
164160	100 mm	16.8 mm	yes	● red	≤12 ml	12 ml	bayonet cap	+	5 / 1,000
191160	95 mm	18 mm	no		≤12.5 ml	14 mI		+	1 / 750

⁷ The dimensions and volumes of our tubes are only nominal values. For exact dimensions and volumes, please refer to the product data sheets on our website: www.gbo.com.

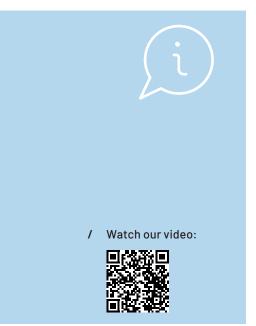



EASYSTRAINER CELL STRAINERS

With EASYstrainer cell strainers, Greiner Bio-One offers an innovative and user-friendly solution for the filtration of cell suspensions.

EASYstrainer can be used, for example, after enzymatic tissue digestion for primary cell isolation or for cell preparation prior to flow cytometry. The large cell strainers fit on all conical 50 ml tubes and are available with mesh sizes of 40, 70 and 100 μm. EASYstrainer Small fits 15 ml tubes as well as smaller tubes and reaction vessels. It is available with mesh sizes of 20, 40, 70 and 100 μm.

Thestackability of EASY strainers with different mesh sizes allows for the separation of cells with different sizes in one step. In addition, the upper part of EASYstrainer Small can be inverted. This allows retained cells to be rinsed out for further use. EASYstrainer makes working aseptically much easier: It can be held via a surrounding rim or a handle in order to avoid accidental contact with the sterile filter material. Additional safety is provided by the transparent blister packaging from which the strainers can be conveniently and aseptically removed.

- / Flexible fit, suitable for tubes from 1.5 ml to 50 ml
- / Available Mesh Sizes: 20, 40, 70 und 100 μm
- / No liquid overspill



EASYstrainer

- / Flexible fit, suitable for tubes from 1.5 ml to 50 ml
- / Handle and ridged skirt for improved aseptic handling
- / Venting slot for fast filtration
- / No liquid overspill
- / Single packaging reduces risk of contamination

















Item no.	Description	Mesh size	Product colour	Sterile	Qty. inner / outer
542040	EASYstrainer for 50 ml tubes	40 µm	● green	+	1/50
542070	EASYstrainer for 50 ml tubes	70 µm	blue	+	1/50
542000	EASYstrainer for 50 ml tubes	100 µm	yellow	+	1/50
542120	EASYstrainer Small for tubes: 1.5 / 5 / 15 ml and 12x75 mm	20 μm	● red	+	1/50
542140	EASYstrainer Small for tubes: 1.5 / 5 / 15 ml and 12x75 mm	40 µm	● green	+	1/50
542170	EASYstrainer Small for tubes: 1.5 / 5 / 15 ml and 12x75 mm	70 µm	blue	+	1/50
542100	EASYstrainer Small for tubes: 1.5 / 5 / 15 ml and 12x75 mm	100 μm	yellow	+	1/50

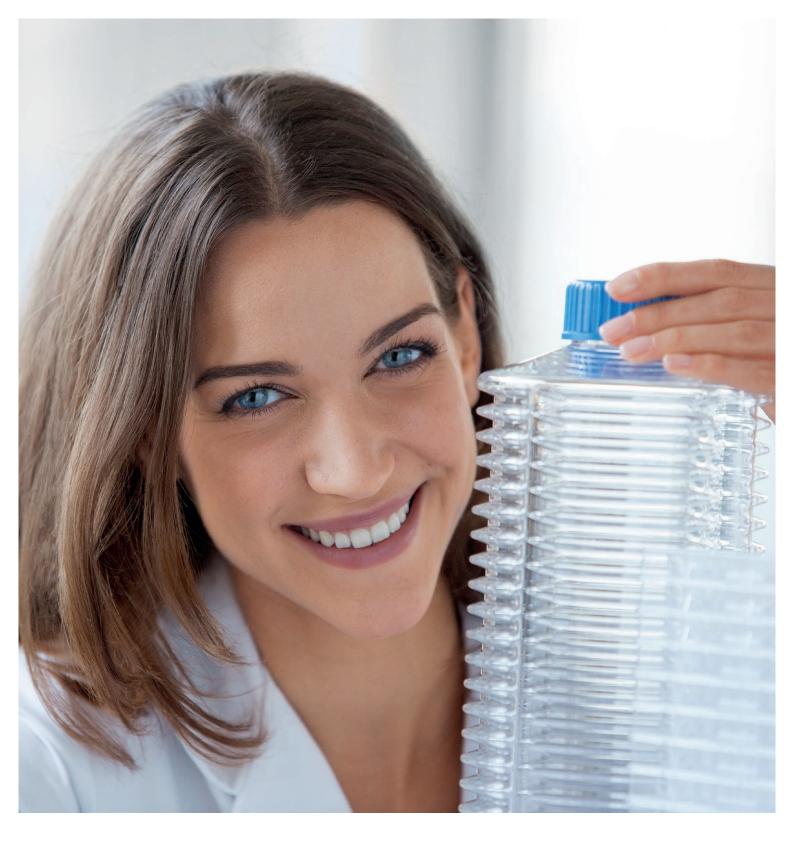

The cultivation of cells as mass cultures has become increasingly important over the past few decades. Mass cell cultures are mainly used for the production of vaccines or recombinant proteins for therapeutic approaches.

MASS CELL CULTURE

/	CELLMASTER Cell Culture Roller Bottles	3
		.74
	Roller Bottles Polystyrene	.75
	Roller Bottles Polyethylene	
	Terephthalate(PET)	.76
/	CELLdisc Cell Culture Device	.77
	CELLdisc 1 / 4 / 8 / 12 / 16 / 24 / 40 Layers	.78
	CELLdisc External Filter	.79
	CELLdisc Closed Filling Caps	.80
	CELLhandle	. 8
	CELLstage Filling Accessory	. 8
	CELLevator	.82
	CELL ring	.82

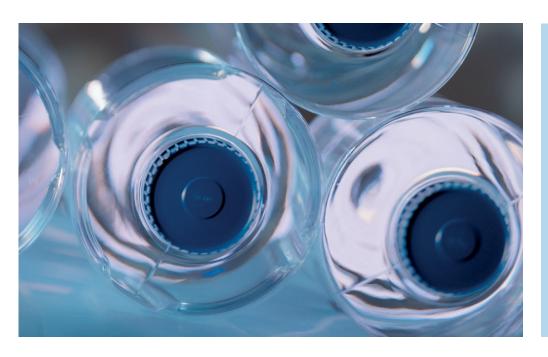



/ For further information, please refer to our website: www.gbo.com



CELLMASTERCELL CULTURE ROLLER BOTTLES

Roller bottles made of polystyrene or PET range in size from $850\,\mathrm{cm^2}$ to $4250\,\mathrm{cm^2}$ growth area and are available either with standard or filter screw caps. The filter membrane with a pore size of $0.2\,\mu\mathrm{m}$ guarantees optimal protection against contamination. It provides a high gas exchange rate which enables a mass cultivation of cells using e.g. bicarbonate buffer and $\mathrm{CO_2}$. The safety screw cap enables a tight closing and contamination-free cultivation. All roller

bottles are sterilised by irradiation according to validated procedures (ISO 11137). Endotoxin testing is conducted in accordance with USP 85 with a tolerance level of 0.03 EU/ml.

The product range includes a short and a long form, that are labelled as X and XL accordingly. Both sizes are available with a smooth or radially ribbed surface. The ribbed design increases the growth area of the roller bottle without changing space requirements.

- / PS or PET roller bottles
- / Different sizes with or without radially ribbed surface
- / Graduations from 200 to 2000 ml
- / Seamless production technique rules out leaking seams



Roller Bottles Polystyrene

- / For adherent cell culture
- / Manufactured from crystal clear polystyrene
- / Certified USP Class VI end product testing
- / Lot number and expiry date on each bottle
- / Filter screw cap with 0.2 µm pore size















Ø: 122 mm, Graduation: yes, Raw material: PS, Surface treatment: TC

Item no.	Flask design	Height	Growth area	Cap colour	Surface	Total volume	Cap design	Sterile	Oty. inner / outer
680660	1 X	271 mm	850 cm ²	blue	smooth	2,520 ml	screw cap	+	2 / 24
680665	1 X	271 mm	850 cm ²	blue	smooth	2,520 ml	screw cap	+	24 / 24
680648	1 X	271 mm	850 cm ²	blue	smooth	2,520 ml	screw cap	+	24 / 48
680658	1 X	271 mm	850 cm ²	blue	smooth	2,520 ml	filter screw cap	+	2 / 24
680668	1 X	271 mm	850 cm ²	blue	smooth	2,520 ml	filter screw cap	+	24 / 24
680645	1 X	271 mm	850 cm ²	blue	smooth	2,520 ml	filter screw cap	+	24 / 48
681670	2.5 X	271 mm	2,125 cm²	blue	ribbed	2,300 ml	screw cap	+	2 / 24
681675	2.5 X	271 mm	2,125 cm ²	blue	ribbed	2,300 ml	screw cap	+	24 / 24
681672	2.5 X	271 mm	2,125 cm²	blue	ribbed	2,300 ml	filter screw cap	+	2 / 24
682660	1XL	500 mm	1,700 cm ²	blue	smooth	4,970 ml	screw cap	+	1 / 12
682612	1XL	500 mm	1,700 cm ²	blue	smooth	4,970 ml	screw cap	+	12 / 12
682624	1XL	500 mm	1,700 cm ²	blue	smooth	4,970 ml	screw cap	+	12 / 24
682625	1XL	500 mm	1,700 cm ²	blue	smooth	4,970 ml	filter screw cap	+	12 / 24
682615	1 XL	500 mm	1,700 cm ²	blue	smooth	4,970 ml	filter screw cap	+	12 / 12
682670	5 XL	500 mm	4,250 cm ²	blue	ribbed	4,640 ml	screw cap	+	1 / 12
682672	5 XL	500 mm	4,250 cm ²	blue	ribbed	4,640 ml	screw cap	+	12 / 24
682678	5 XL	500 mm	4,250 cm ²	blue	ribbed	4,640 ml	filter screw cap	+	1 / 12

 $[\]emph{I} \quad \text{Thread enables quick opening with a 2/3 turn. Screw caps with larger knurls for improved grip and ease of opening \textit{I} closing.}$

 $[\]emph{I} \quad \textbf{Sterile, individually packed screw caps: standard (item no. 383361) and filter (item no. 383382)}.$



Roller Bottles

Polyethylene Terephthalate (PET)

- / For adherent cell culture
- / Certified USP Class VI end product testing
- / With standard screw cap

















Flask design: 1 X, Height: 271 mm, Ø: 116 mm, Growth area: 850 cm², Graduation: yes, Raw material: PET, Surface: smooth, Total volume: 2,300 ml, Cap design: screw cap

Item no.	Cap colour	Sterile	Oty. inner / outer
680190	blue	+	-/18
680195	blue	+	-/30

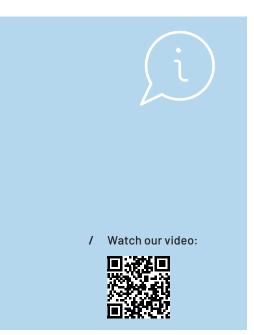

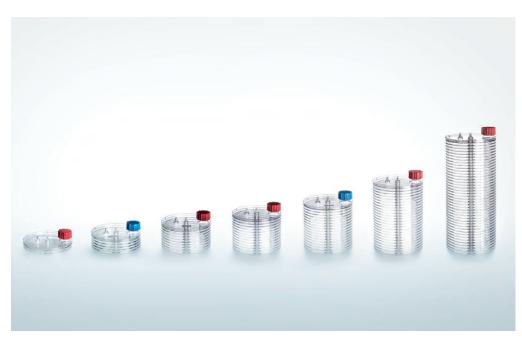

CELL CULTURE DEVICE

CELLdisc is a multilayer cell culture device covering a range of 250 cm² to 10,000 cm² growth area. The innovative ergonomic design provides a versatile system from research scale to industrial batches with a 40 % higher surface/volume ratio then conventional multilayer systems.

A centrally located ventilation channel assures uniform distribution of gas throughout the device whereas the wide opening port simplifies manual filling. The compact and robust cylindrical device is ideally suited for automation and upscaling of cell

cultures. For the connection of the individual layers a proprietary, particle, and adhesive free assembly technique is used and the complete end product is USP Class VI certified.

To guarantee an ideal cell culture environment CELLdisc is provided with two surface treatments: Beside the TC surface for standard cells and applications CELLdisc is also available with the Advanced TC surface, improving cultivation of sensitive cells, even under restricted growth conditions and increasing cellular adhesion, proliferation and transfection rates.

- / Mass Cell Culture
- / Antibody, virus and vaccine production
- / Production of recombinant or therapeutic proteins



CELLdisc

1 / 4 / 8 / 12 / 16 / 24 / 40 Layers

- / 40 % higher surface/volume ratio than comparable systems
- / Easy operation and minimal space occupation
- / Media exchange without contact to cell layers





FREE OF detectable human DNA FREE OF detectable RNase





Ø: 200 mm, Raw material: PS, Cap design: screw cap

Item no.	Layers	Height	Growth area	Surface treatment	Cap colour	Working volume	Sterile	Qty. inner / outer
678101	1	61 mm	250 cm ²	TC	● red	15 ml - 50 ml	+	1/8
678104	4	93 mm	1,000 cm ²	TC	● red	60 ml - 200 ml	+	1/4
678108	8	135 mm	2,000 cm ²	TC	● red	120 ml - 400 ml	+	1/3
678112	12	177 mm	3,000 cm ²	TC	● red	180 ml - 600 ml	+	1/2
678116	16	220 mm	4,000 cm ²	TC	● red	240 ml - 800 ml	+	1/2
678124	24	304 mm	6,000 cm ²	TC	● red	360 ml - 1,200 ml	+	1/2
678140	40	474 mm	10,000 cm ²	TC	● red	600 ml - 2,000 ml	+	1/1
678904	4	93 mm	1,000 cm ²	Advanced TC	blue	60 ml - 200 ml	+	1/4
678908	8	135 mm	2,000 cm ²	Advanced TC	blue	120 ml - 400 ml	+	1/3
678912	12	177 mm	3,000 cm ²	Advanced TC	blue	180 ml - 600 ml	+	1/2
678916	16	220 mm	4,000 cm ²	Advanced TC	blue	240 ml - 800 ml	+	1/2
678924	24	304 mm	6,000 cm ²	Advanced TC	blue	360 ml - 1,200 ml	+	1/2
678940	40	474 mm	10,000 cm ²	Advanced TC	blue	600 ml - 2,000 ml	+	1/1



CELLdisc External Filter

Feature: External Filter, Ø: 200 mm, Raw material: PS, Surface treatment: TC, Cap design: screw cap

- / Triple-packed CELLdiscs, pre-assembled with external filters
- / USP VI certified venting filters, manufactured with reinforced hydrophobic PTFE membranes
- / A defined pore size of 0.2 µm prevents contamination and ensures sufficient airflow to the cells



Item no.	Layers	Height	Growth area	Cap colour	Working volume	Sterile	Qty. inner / outer
678101-EXF	1	61 mm	250 cm ²	● red	15 ml - 50 ml	+	1 / 4
678104-EXF	4	93 mm	1,000 cm ²	● red	60 ml - 200 ml	+	1/3
678108-EXF	8	135 mm	$2,000 \text{cm}^2$	●red	120 ml - 400 ml	+	1/2
678112-EXF	12	177 mm	$3,000 \text{cm}^2$	● red	180 ml - 600 ml	+	1/2
678116-EXF	16	220 mm	$4,000 \text{cm}^2$	● red	240 ml - 800 ml	+	1/2
678124-EXF	24	304 mm	6,000 cm ²	● red	360 ml - 1,200 ml	+	1/2
678140-EXF	40	474 mm	10,000 cm ²	●red	600 ml - 2,000 ml	+	1/1

[/] Assembled product is sterilized (SAL $10^{\text{-}6}$) to reduce risk of contamination



CELLdisc Closed Filling Caps

- / Triple-packed CELLdiscs, pre-assembled with external filters and closed filling cap
- / Choice of single tubing (CF1) or double tubing (CF2) depending on desired emptying process
- / Tubing with MPC type connector for quick and safe connection / disconnection

Ø: 200 mm, Raw material: PS, Surface treatment: TC, Cap design: screw cap

Item no.	Feature	Layers	Height	Growth area	Cap colour	Working volume	Sterile	Qty. inner / outer
678101-CF1	hose connector without Dip-In	1	61 mm	250 cm ²	● red	15 ml - 50 ml	+	1/4
678104-CF1	hose connector without Dip-In	4	93 mm	1,000 cm ²	● red	60 ml - 200 ml	+	1/3
678104-CF2	hose connector with Dip-In	4	93 mm	1,000 cm ²	● red	60 ml - 200 ml	+	1/3
678108-CF1	hose connector without Dip-In	8	135 mm	2,000 cm ²	● red	120 ml - 400 ml	+	1/2
678108-CF2	hose connector with Dip-In	8	135 mm	2,000 cm ²	● red	120 ml - 400 ml	+	1/2
678112-CF1	hose connector without Dip-In	12	177 mm	3,000 cm ²	● red	180 ml - 600 ml	+	1/2
678112-CF2	hose connector with Dip-In	12	177 mm	3,000 cm ²	● red	180 ml - 600 ml	+	1/2
678116-CF1	hose connector without Dip-In	16	220 mm	4,000 cm ²	● red	240 ml - 800 ml	+	1/2
678116-CF2	hose connector with Dip-In	16	220 mm	4,000 cm ²	● red	240 ml - 800 ml	+	1/2
678124-CF1	hose connector without Dip-In	24	304 mm	6,000 cm ²	● red	360 ml - 1,200 ml	+	1/2
678124-CF2	hose connector with Dip-In	24	304 mm	6,000 cm ²	● red	360 ml - 1,200 ml	+	1/2
678140-CF1	hose connector without Dip-In	40	474 mm	10,000 cm ²	● red	600 ml - 2,000 ml	+	1/1
678140-CF2	hose connector with Dip-In	40	474 mm	10,000 cm ²	● red	600 ml - 2,000 ml	+	1/1

 $[\]emph{\emph{I}}$ Assembled product is sterilized (SAL 10 $^{\text{-6}}$) to reduce risk of contamination



CELLhandle

- Gripping device for easy lifting and emptying of large-sized CELLdisc
- / Enables single-hand usage

Item no.	0ty. inner / outer
878074	1/1



CELLstage

Filling Accessory

- / Available for CELLdisc 4-24 layers and CELLdisc 40 layers
- Creates the optimum angle and position for CELLdisc filling
- / Stainless steel allows multiple sterilization methods
- / Suitable for left- and right handed users

Item no.	Description	Qty. inner / outer
878072	for CD4 - CD24	1/1
878073	for CD40	1/1



CELLevator

Easy and secured CELLdisc stacking

- / Maximum loading capacity 8 kg
- / Space saving storage
- / Autoclavable (120 °C, 2 bar)

Item no.	Qty. inner / outer
878071	1/9

/ For more information please see instruction for use No. F073253 on www.gbo.com



CELLring

Levelling ring

/ Ensures exact planar positioning of CELLdisc

Item no.	Qty. inner / outer
878075	-/3

TECHNICAL NOTE



CELLdisc

COMPARISON OF INCUBATOR SPACE USAGE BETWEEN GREINER BIO-ONE **CELLdisc** AND RECTANGULAR MULTILAYER DEVICES

The aim of mass cell culture is to produce large quantities of cells mostly used for industrial or clinical applications. Users can choose from a variety of different disposables and concepts. To reduce the cost of the scale up process to a minimum, different factors need to be considered and evaluated.

If multilayer devices such as the Greiner Bio-One CELL disc are to be used, not only growth area and the number of cells per device should be taken into account but also how effectively incubator space is utilised by them.

The innovative CELL disc with its round and ergonomic design stands out from existing rectangular devices. The cuboid nature of these alternatives would, at first glance, imply an optimal use of space within an incubator that is itself cuboid in design. However, in reality, the length and width of such systems do not correspond ideally to the dimensions of standard incubators and leave a lot of space unused. Additionally such devices cannot be positioned next to each other as this could compromise consistent thermal distribution, because when units touch each other so access to the temperature environment of the incubator is compromised.

CELLdisc however, thanks to the cylindrical nature of the design, can be placed directly next to one another and still receive uniform exposure to the optimum temperatue of the incubator. Due to this fact and the 40% higher surface volume/ratio CELLdisc offers over standard competitive systems, utilisation of incubator space is far more efficient when using CELLdisc than rectangular devices. (See Table 1)

TABLE 1

Utilisation of incubator space - exemplary calculation of multilayer devices with comparible dimesions

	CELLdisc 16 layers	Competitor A 10 layers, rectangular form	Competitor B 10 layers, rectangular form (high density version)
Dimensions (L x W x H incl. cap)	ø 20 cm, H 22 cm	33 cm x 20 cm x 20.5 cm	33 cm x 23 cm x 21 cm
Growth surface (device)	4.000 cm ²	6.360 cm ²	8.216 cm²
Max. devices / shelf	6	2	2
Max. devices / incubator	12	4	4
Growth surface (incubator load)	12 x 4.000 = 48.000 cm ²	$4 \times 6.360 = 25.440 \text{ cm}^2$	$4 \times 8.216 =$ 32.864 cm^2
Utilisation	100 %	53 %	68.47 %

FIG. 1.

Standard incubator, eg: HeraCell 240 (Thermo Fisher Scientific) Incubator size internal dimensions: W 56 cm, H: 67 cm



Fig. 1a: Front view of the incubator:
1) Top shelf: Cuboid multilayer devices
2) Lower shelf: CELLdisc

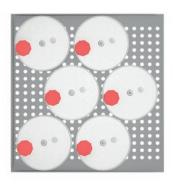


Fig. 1b: Cross section of the incubator. Optimal positioning of CELL disc and utilisation of incubator space.



In 3D cell culture, cells form in a spatial orientation - similar to the body. This enables, for example, the replacement of animal experiments and - usually as high-throughput screening - research on drugs, stem cells and tumor development.

3D CELL CULTURE

/	CELLSTAR* Cell Culture Vessels
	Cell-Repellent Surface
	Cell Culture Dishes Cell-Repellent Surface87
	Cell Culture Flasks Cell-Repellent Surface87
	Multiwell Plates / Microplates
	Cell-Repellent Surface88
/	Magnetic 3D Cell Culture 89
	Spheroid Bioprinting 96 Well90
	Spheroid Bioprinting 384 Well90
	Magnetic Levitation 6 / 24 Well91
	Screening & Imaging 96 / 384 Well92
	MagPen Single / 24 Well / 96 Well93
	Consumables / Accessories Magnetic
	3D Cell Culture93
/	ThinCert® Cell Culture Inserts for 6, 12
	and 24 Well Multiwell Plates94
	ThinCert® Cell Culture Inserts 6 Well95
	ThinCert® Cell Culture Inserts 12 Well95
	ThinCert® Cell Culture Inserts 24 Well96
	ThinCert® Plate 6 / 12 Well 97





- / Forum No. 17: CELLSTAR® Cell Culture Vessels with Cell-Repellent Surface (F073777)
- / Application Report "Advantage of CELLSTAR®
 Cell Culture Vessels
 with Cell-Repellent
 Surface for 3-D Cell
 Culture in Hydrogels"
 (F073792)

CELLSTAR° CELL CULTURE VESSELS CELL-REPELLENT SURFACE

Cell culture is an essential tool in drug discovery, tissue engineering, toxicology testing, stem cell research, as well as in basic research.

Beside conventional two-dimensional (2D) monolayer cell culture, 3D cell culture models are becoming a routine tool which enables the expression of extracellular matrix (ECM) components as well as the formation of cell-cell and cell-matrix interactions. These characteristics are important for replicating

in-vivo cell differentiation, proliferation, and function in vitro. Greiner Bio-One developed CELLSTAR® cell culture vessels with cell-repellent surface specifically for culturing in 3D. The cell-repellent surface effectively prevents cell adherence and therefore can promote the spontaneous formation of three-dimensional spheroids. Cell culture vessels with cell-repellent surface are also ideal platforms for long-term cultivation in hydrogels.

- / Spheroid and organoid culture
- / Aggregation of stem cells
- / Suspension culture of semi-adherent/ adherent cell lines
- / 3D culture in hydrogels



Cell Culture Dishes Cell-Repellent Surface

Cell culture vessels with cell-repellent surface reliably prevent cell attachment in suspension cultures of semi adherent / adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.





FREE OF detectable human DNA

FREE OF detectable RNase





Raw material: PS, Surface treatment: cell-repellent, Vent nock: yes

Item no.	Height	Ø nominal size	Working volume	Total volume	Sterile	Qty. inner / outer
627979	10 mm	35 mm	≤3 ml	10 ml	+	10 / 40
628979	15 mm	60 mm	6 ml - 7 ml	28 ml	+	10 / 20
664970	20 mm	100 mm	16 ml - 17 ml	100 ml	+	1/5



Cell Culture Flasks

Cell-Repellent Surface

A cell-repellent surface reliably prevent cell attachment in suspension cultures of semi-adherent and adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.













Raw material: PS, Surface treatment: cell-repellent

Item no.	Flask design	Cap colour	Total volume	Cap design	Sterile	Qty. inner / outer
690980	flat	○ white	50 ml	screw cap	+	10 / 20
690985	flat	○ white	50 ml	filter screw cap	+	10 / 20
658980	flat	○white	250 ml	screw cap	+	5 / 15
658985	flat	○ white	250 ml	filter screw cap	+	5 / 15
660980	flat	○white	550 ml	screw cap	+	5/5
660985	flat	○ white	550 ml	filter screw cap	+	5/5
661980	high	○white	650 ml	screw cap	+	4/4
661985	high	○ white	650 ml	filter screw cap	+	4/4



Multiwell Plates / Microplates Cell-Repellent Surface

- / 6 / 12 / 24 / 48 well multiwell plates available
- / 96 / 384 well plates with various well geometries and optional µClear® film bottom

STERILE

FREE OF detectable DNase FREE OF detectable numan DNA

FREE OF detectable RNase





Raw material: PS, Surface treatment: cell-repellent

Item no.	Well format	Well profile	Bottom	Product colour	Total volume (well)	Working volume (well)	Lid	Sterile	Oty. inner / outer
657970	6	F-bottom	solid	○ clear	16.1 ml	2 ml - 5 ml	yes, conden- sation rings	+	1/5
665970	12	F-bottom	solid	○ clear	6.5 ml	2 ml - 4 ml	yes, conden- sation rings	+	1/5
662970	24	F-bottom	solid	○clear	3.3 ml	0.5 ml - 1.5 ml	yes, conden- sation rings	+	1/5
677970	48	F-bottom	solid	○ clear	1.7 ml	0.5 ml - 1 ml	yes, conden- sation rings	+	1/5
650970	96	U-bottom	solid	○ clear	323 µI	40 μl - 280 μl	yes, conden- sation rings	+	1/6
650979	96	U-bottom	solid	○ clear	323 µI	40 μl - 280 μl	yes, conden- sation rings	+	8 / 32
651970	96	V-bottom	solid	○ clear	234 μΙ	40 μl - 200 μl	yes, conden- sation rings	+	1/6
655970	96	F-bottom / Chimney Well	solid	○ clear	392 μΙ	25 μl - 340 μl	yes, conden- sation rings	+	1/6
655976	96	F-bottom / Chimney Well	µClear®	● black	392 µI	25 μl - 340 μl	yes, conden- sation rings	+	8 / 32
655976-SIN	96	F-bottom / Chimney Well	μClear®	● black	392 μΙ	25 μl - 340 μl	yes, conden- sation rings	+	1/32
781970	384	F-bottom	solid	○ clear	131 μΙ	15 μl - 110 μl	yes	+	1/60
781974	384	F-bottom	µClear®	○white	131 μΙ	15 μl - 110 μl	yes	+	8 / 32
781976	384	F-bottom	µClear®	● black	131 μΙ	15 μl - 110 μl	yes	+	8 / 32
781976-SIN	384	F-bottom	µClear®	● black	131 μΙ	15 μl - 110 μl	yes	+	1/32
787979	384	U-bottom	solid	○ clear	122 μΙ	10 µl - 90 µl	yes	+	8 / 32



- / Brochure "3D Cell Culture" (F071076)
- / For further information please refer to our microsite:






MAGNETIC 3D CELL CULTURE

The core technology of Greiner Bio-One's magnetic 3D cell culture is the magnetisation of cells with NanoShuttle-PL. The cells can be aggregated with magnetic forces, either by levitation or Bioprinting, to form structurally and biologically representative 3D models in vitro. NanoShuttle-PL consists of gold, iron oxide, and Poly-L-Lysine. These nanoparticles (\emptyset < 50 nm) magnetise cells by electrostatically attaching to cell membranes during an overnight static incubation. Magnetised cells will appear peppered with dark nanoparticles after incubation.

NanoShuttle-PL is biocompatible, having no effect on metabolism, proliferation and inflammatory stress. Additionally, it does not interfere with experimental techniques, such as fluorescence or Western blotting. With magnetised spheroids, solution addition and removal are made easy by using magnetic force to hold them in a stationary position during aspiration, thereby limiting spheroid loss.

- / 3D in a 2D workflow
- / Easy handling
- / Fast 3D tissue assembly
- / No sample loss
- / Scalable
- / Ready for automation

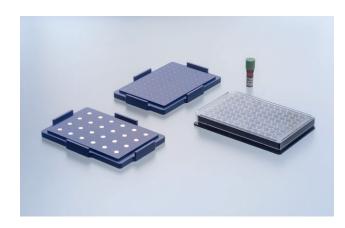

Spheroid Bioprinting 96 Well

Magnetized cells are brought together to form spheroids at the well bottom using weak magnetic forces.

Item no.	Description	Content kit	Qty. inner / outer
655840	96 Well Bioprinting Kit, clear	NanoShuttle-PL (3 vials), Spheroid Drive, Holding Drive, 96 well cell culture microplates (clear) with cell-repellent surface (2 x 655970)	1/1
655841	96 Well Bioprinting Kit, black, µClear®	NanoShuttle-PL (3 vials), Spheroid Drive, Holding Drive, 96 well cell culture microplates (black, µclear®) with cell-repellent surface (2 x 655976-SIN)	1/1
655850	96 Well Ring Drive	96 Well Ring Drive for the formation of 3D ring structures	1/1
655830	96 Well Spheroid and Holding Drive	Spheroid Drive, Holding Drive	-/1



Spheroid Bioprinting 384 Well

Magnetized cells are brought together to form spheroids at the well bottom using weak magnetic forces.

Item no.	Description	Content kit	Qty. inner / outer
781840	384 Well Bioprinting Kit, clear	NanoShuttle-PL (2 vials), Spheroid Drive, Holding Drive, 384 well cell culture microplates (clear) with cell-repellent surface (2 x 781970)	1/1
781841	384 Well Bioprinting Kit, black, μClear®	NanoShuttle-PL (2 vials), Spheroid Drive, Holding Drive, 384 well cell culture microplates (black, µClear®) with cell-repellent surface (2x 781976- SIN)	1/1
781850	384 Well Ring Drive	384 Well Ring Drive for the formation of 3D ring structures	1/1

Item no.	Description	Content kit	Qty. inner / outer
781830	384 well spheroid and holding drive	Spheroid Drive, Holding Drive	-/1



Magnetic Levitation 6 / 24 Well

Magnetised cells are levitated off the plate bottom by a magnet and rapidly form aggregates.


Item no.	Description	Content kit	Sterile	Qty. inner / outer
657840	6 Well Bio-Assembler Kit	Levitation Drive, Holding Drive, NanoShuttle-PL (2 Vials), 6 well cell culture multiwell plates (2 x 657970) and 6 Well Intermediate lid (2 x 657825) with cell-repellent surface		1/1
657825	6 Well Intermediate lid	Intermediate lid with cell-repellent surface	+	2 / 10
657830	6 Well Levitation and Holding Drive	Levitation Drive (1), Holding Drive (1)		-/1
662840	24 Well Bio-Assembler Kit	Levitation Drive, Holding Drive, NanoShuttle-PL (2 Vials), 24 well cell culture multiwell plates (2 x		1/1
662825	24 Well Intermediate lid	Intermediate lid with cell-repellent surface	+	1 / 10
662830	24 well Levitation and Holding Drive	Levitation Drive (1), Holding Drive (1)		-/1



Screening & Imaging 96 / 384 Well

Ideal kits for combining applications and examination methods.

Item no.	Description	Content kit	Qty. inner / outer
655846	96 Well BiO Assay Kit	NanoShuttle-PL (3 vials), 6 Well Levitation Drive, 96 Well Spheroid, Holding and Ring Drive, 96 Well Deep Well Plate, 6 Well cell culture multiwell plates with cell-repellent surface (2 x 657970), 96 Well cell culture microplates (clear) with cell-repellent surface (2 x 655970)	1/1
655849	96 Well BiO Assay Kit & Imaging System	NanoShuttle-PL (3 vials), 6 Well Levitation Drive, 96 Well Spheroid, Holding and Ring Drive, 96 Well Deep Well Plate, 6 Well cell culture multiwell plates with cell-repellent surface (2 x 657970), 96 Well cell culture microplates (transparent) with cell-repellent surface (2 x 655970), Imaging Kit (657860)	1/1
781846	384 Well BiO Assay Kit	NanoShuttle-PL (2 vials), 6 Well Levitation Drive, 384 Well Spheroid, Holding and Ring Drive, 96 Well Deep Well plate, 6 Well cell culture multiwell plates with cell-repellent surface (2 x 657970), 384 Well cell culture microplates (clear) with cell-repellent surface (2 x 781970)	1/1
781849	384 Well BiOAssay Kit & Imaging System	NanoShuttle-PL (2 vials), 6 Well Levitation Drive, 384 Well Spheroid, Holding and Ring Drive, 96 Well Deep Well plate, 6 Well cell culture multiwell plates with cell-repellent surface (2 x 657970), 384 Well cell culture microplates (clear) with cell-repellent surface (2 x 781970), Imaging Kit (657860)	1/1
657860	Imaging Kit	Imaging System, Light Pad, Cooling Fan	-/1

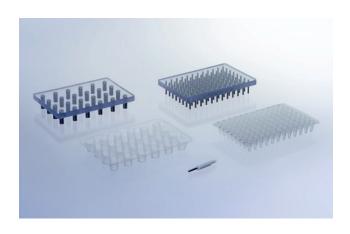

MagPen Single / 24 Well / 96 Well

MagPen is a smart assistant for easy and fast transfer and collection of magnetized cell cultures by a simple "pick up-and-drop"-step.

Item no.	Description	Content kit	Sterile	Qty. inner / outer
657850	MagPen 3-pack	Teflon caps (3), magnets (3)		-/3
657824	24 Well Multi-MagPen Kit	24 Well Multi-MagPen Drive and 24 Well Multi-MagPen Sleeve (2 x 651524) with cell-repellent surface		-/1
651524	24 Well Multi-MagPen Sleeve	Multi-MagPen Sleeve with cell-repellent surface	+	1 / 10
657896	96 Well Multi-MagPen Kit	96 Well Multi-MagPen Drive und 96 Well Mul- ti-MagPen Sleeve (2 x 61596) with cell-repellent surface		-/1
651596	96 Well Multi-MagPen Sleeve	Multi-MagPen Sleeve with cell-repellent surface	+	1/10



Consumables / Accessories Magnetic 3D Cell Culture

- / NanoShuttle-PL consists of gold, iron oxide and Poly-L-Lysine
- / MagPen is a magnetic tool to pick up the spheroids

Item no.	Description	Content kit	Qty. inner / outer
657841	NanoShuttle-PL	600 µl vials of NanoShuttle-PL (1)	-/1
657843	NanoShuttle-PL 3-pack	600 µl vials of NanoShuttle-PL (3)	-/3
657846	NanoShuttle-PL 6-pack	600 µl vials of NanoShuttle-PL (6)	-/6
657852	NanoShuttle-PL 12-pack	600 µl vials of NanoShuttle-PL (12)	-/12
657847	NanoShuttle-PL 6-pack with free iPod	NanoShuttle-PL (6), free iPod (with purchase of either 655849 or 781849)	-/1





- / Forum No. 8: ThinCert®
 cell culture products Overview (F073017)
- / Application Note "Immuno cytochemistry"(F073100)
- / For further information and application notes please refer to our website: www.gbo.com

THINCERT® CELL CULTURE INSERTS FOR 6, 12 AND 24 WELL MULTIWELL PLATES

For advanced cell and tissue culture applications,
Greiner Bio-One offers a broad range of membrane supports
- ThinCert®.

Combining 6 different membrane specifications (pore size and density) in geometries to fit 6, 12 and 24 well plates, the ThinCert® cell culture inserts are suitable for a wide range of applications including transport, secretion and diffusion studies, migrational experiments, cytotoxicity testing, co-cultures, trans epithelial electric resist-

well as primary cell culture. ThinCert® cell culture inserts are compatible with standard CELLSTAR® multiwell plates from Greiner Bio-One, and are pre-packed together with the requisite number of plates. The automated production process includes double optical control of each insert produced, ensuring that any biological contamination is avoided. The sterility of the single blisterpacked inserts and multiwell plates is ensured by irradiation.

ance (TEER) measurements, as

- / Stable clear polystyrene housing
- / Sealed PET capillary pore membrane
- / Pre-configured multiwell plates with ThinCert® cell culture inserts available on request



ThinCert® Cell Culture Inserts 6 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange



Feature: 4 multiwell plates / box, Height: 16.25 mm, Ø internal: 24.85 mm, Ø external: 27.85 mm, Cultural surface: 452.4 mm², Surface treatment: TC, Working volume (ThinCert®): 1 ml - 3.6 ml, Working volume (well): 2 ml - 4.15 ml

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
657640	$1 \times 10^{8} \text{cm}^{2}$	0.4 µm	translucent	+	1 / 24
657641	$2 \times 10^{6} \text{ cm}^{2}$	0.4 µm	clear	+	1 / 24
657610	$2 \times 10^{6} \text{ cm}^{2}$	1 µm	clear	+	1 / 24
657630	0,6 x 10 ⁶ cm ²	3 µm	clear	+	1 / 24
657631	$2 \times 10^{6} \text{ cm}^{2}$	3 µm	translucent	+	1 / 24
657638	$0.15 \times 10^{6} \text{cm}^{2}$	8 µm	translucent	+	1 / 24



ThinCert® Cell Culture Inserts 12 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange

Feature: 4 multiwell plates / box, Height: 16.25 mm, Ø internal: 13.85 mm, Ø external: 15.85 mm, Cultural surface: 113.1 mm², Surface treatment: TC, Working volume (ThinCert®): 0.3 ml - 1 ml, Working volume (well): 1 ml - 2 ml

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
665640	1 x 10 ⁸ cm ²	0.4 µm	translucent	+	1 / 48
665641	$2 \times 10^{6} \text{cm}^{2}$	0.4 µm	clear	+	1 / 48
665610	2 x 10 ⁶ cm ²	1 µm	clear	+	1 / 48

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Oty. inner / outer
665630	0,6 x 10 ⁶ cm ²	3 µm	clear	+	1 / 48
665631	$2 \times 10^{6} \text{cm}^{2}$	3 µm	translucent	+	1 / 48
665638	$0.15 \times 10^6 \text{cm}^2$	8 µm	translucent	+	1 / 48



ThinCert® Cell Culture Inserts 24 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange











Feature: 2 multiwell plates / box, Height: 16.25 mm, Ø internal: 8.4 mm, Ø external: 10.4 mm, Cultural surface: 33.6 mm², Surface treatment: TC, Working volume (ThinCert®): 0.1 ml - 0.35 ml, Working volume (well): 0.4 ml - 1.2 ml

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
662640	1 x 10 ⁸ cm ²	0.4 μm	translucent	+	1 / 48
662641	$2 \times 10^{6} \text{cm}^{2}$	0.4 μm	clear	+	1 / 48
662610	$2 \times 10^{6} \text{cm}^{2}$	1 µm	clear	+	1 / 48
662630	0,6 x 10 ⁶ cm ²	3 µm	clear	+	1 / 48
662631	$2 \times 10^{6} \text{cm}^{2}$	3 µm	translucent	+	1 / 48
662638	0,15 x 10 ⁶ cm ²	8 µm	translucent	+	1 / 48



ThinCert® Plate 6 / 12 Well

- / Optimised for use with ThinCert® cell culture inserts
- / Deep wells for an increased volume of medium in air-lift culture
- / Notches for fixed insert position
- / Available in 6 and 12 well format














Height: 39.5 mm, Length: 129.5 mm, Width: 86.6 mm, Lid: yes, condensation rings

Item no.	Well format	Working volume (well)	Sterile	Qty. inner / outer
657110	6	≤20 ml	+	1/50
665110	12	≤4 ml	+	1 / 60

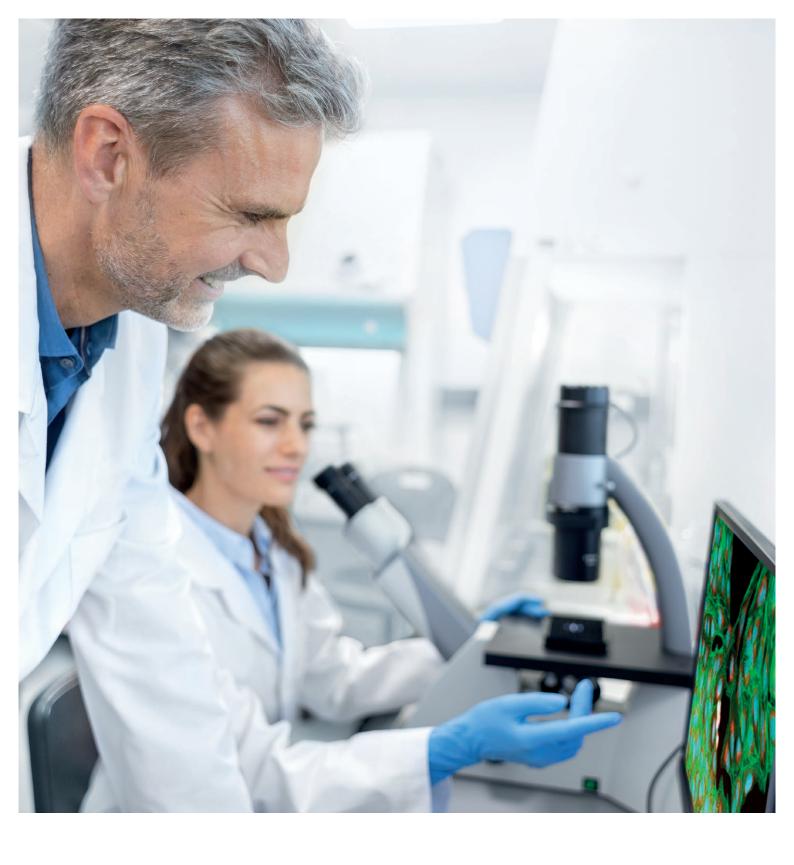

Microscopy is a basic and important method, which is often used in research as well as in medical diagnostics. For this purpose, Greiner Bio-One offers a variety of tailor-made solutions, which guarantee optimal basic conditions for microscopic experiments.


PRODUCTS FOR MICROSCOPY

/	CELLview Dish Cell Culture Dish with Glass Bottom
/	CELLview Slide Cell Culture Slide with Glass Bottom
/	CELLview Plate Cell Culture Plate with Glass Bottom
/	SCREENSTAR Microplates110 SCREENSTAR Microplates 96 / 384 / 1536 Well.111
/	SensoPlate Glass Bottom Microplates112 SensoPlate Glass Bottom Plates 24/96/384/ 1536 Well113

PRODUCTS FOR MICROSCOPY

Well distance equal to a 96 well microplate

Positioning notch for automated microscopy in well A1

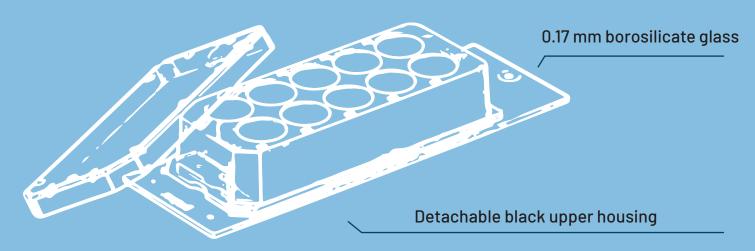

10 wells with alphanumeric coding



PRODUCTS FOR MICROSCOPY

Technological progression in confocal microscopy, optical systems and emerging technologies continues to elevate microscopy as one of the most powerful tools in cell biology. With its advantages for molecular selectivity and capability of live observation, fluorescence microscopy currently is among the most widely used approaches for high-resolution, non-in-

vasive imaging of living cells. Depending on the complexity of live cell imaging experiments and the requirements of the corresponding microscope, the requirements for the utilised disposables can be as comprehensive. Greiner Bio One's imaging consumables are tailored solutions to provide optimal basic settings for your microscopic experiment.

FEATURES:

- / Maximal light transmission
- / Innovative design for maximal planarity
- / Reduced meniscus effect
- Optimal cell attachment and viability



CELLview Products
Tailored solutions
to provide optimal
basic settings for
your microscopic
experiment

For live cell imaging experiments and high-resolution microscopic applications high numerical aperture (N.A.) objectives are required. Microplates with a standard bottom thickness of Immare much too thick to image when using N.A. objectives. Using such microplates or thicker coverslips will lead to spherical aberration resulting in a loss of contrast and image sharpness.

CELLview products combine the convenience of a plastic disposable with the high optical quality of a 0.17 mm thin cover glass bottom, providing superior high-resolution microscopic images of in-vitro cultures. The specific design with the embedded cover glass bottom guarantees a single-plane, flat bottom with a consistent working distance, maximal planarity and optimal thermal conductivity in heated platforms.





µCLEAR® MICROPLATES

The 190 µm polystyrene film bottom is ideal for standard imaging applications with low to medium magnification

SCREENSTAR microplates with a 0.19 mm cycloolefin film bottom are suitable for sophisticated microscopic applications, in high-content screening (HCS) or high-resolution microscopy with water and oil immersion objectives. Cycloolefins possess

excellent optical features and display a low background in the lower UV, with a refractive index and focus background comparable to glass.



CELLview Products

Dishes, slides and plates with 0.17 mm thin cover glass bottom for superior high-resolution microscopic images of in-vitro cultures.



SCREENSTAR Microplates

96 / 384 / 1536 well plates with a 0.19 mm thin cycloolefin film bottom for sophisticated microscopic applications, in high content screening or high-resolution microcopy.



SensoPlate Glass Bottom Plates

Glass bottom plates without surface treatment for fluorescence correlation spectroscopy and microscopic applications.





- / Application Note "Protein localisation using confocal laser scanning microscopy" (F073101)
- / Application Note "Live cell imaging on Golgi morpholgy using the CELLview dish" (F074048)

CELLVIEW DISH

CELL CULTURE DISH WITH GLASS BOTTOM

CELLview Dish combines the convenience of a standard size 35 mm disposable plastic cell culture dish with the optical quality of glass, providing superior high-resolution microscopic images of in-vitro cultivated cultures.

It is made from high-grade polystyrene combined with an integrated glass bottom. The innovative design of the dish provides a single-plane, flat bottom with a consistent working distance and maximal planarity. Moreover, the dish bottom configuration facilitates optimal thermal conductivity and avoids thermal variations in heated platforms

The subdivided version of CELLview Dish enables simultaneous multiplex analyses of different cell lines, various stimulations or diverse transfections. Ouartering the cell culture dish provides four individual compartments with a growth area of approximately 1.9 cm², allowing minimisation of cells and reagents required per individual assay. In addition to the untreated glass surface, Greiner Bio-One provides a tissue culture surface as well as the Advanced TC surface modification to enhance the attachment of adherent cells, thus eliminating the need for protein coating.

used for live cell imaging.

- / TC and Advanced TC surface available
- / Maximal spectral transmission
- No autofluorescence, no depolarisation of light



CELLview Dish

The CELLview cell culture dish combines the convenience of a standard size 35 mm disposable plastic cell culture dish with the optical quality of glass, providing superior high-resolution microscopic images of *in-vitro* cultivated cultures.

/ Embedded glass bottom for maximal planarity

STERILE FREE OF detectab DNase

FREE OF detectable human DNA FREE OF detectable RNase





Height: 10 mm, Ø: 35 mm, Bottom: glass, Vent nock: yes

Item no.	Compart- ments	Growth area	Growth area / unit	Surface treatment	Working volume	Total volume	Total volume (well)	Working volume (well)	Sterile	Oty. inner / outer
627861	1	8.7 cm ²		untreated	2.5 ml - 5 ml	10 ml			+	10 / 40
627860	1	8.7 cm ²		TC	2.5 ml - 5 ml	10 ml			+	10 / 40
627965	1	8.7 cm ²		Advanced TC	2.5 ml - 5 ml	10 ml			+	10 / 40
627871	4		1.9 cm ²	untreated			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40
627870	4		1.9 cm ²	TC			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40
627975	4		1.9 cm ²	Advanced TC			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40

[/] Application Note "Protein localisation using confocal laser scanning microscopy" (F073101)

 $[\]emph{I} \quad \textbf{Application Note "Live cell imaging on Golgi morpholgy using the CELL view dish" (F074048)}$





/ For further information, please refer to our website: www.gbo.com



CELLVIEW SLIDE

CELL CULTURE SLIDE WITH GLASS BOTTOM

CELLview Slide consists of a transparent slide with a black upper housing that effectively subdivides the slide into 10 compartments, which have been designed to mimic the size and layout of a standard 96 well microplate. Because of this standard layout, the slide is compatible with multichannel pipettes making it simple and efficient to use. Furthermore, the round well design helps to reduce meniscus effects for optimum results in cell culture and microscopic analysis. The slide has a 0.17 mm thin cover glass embedded

in its bottom for improved optical clarity and imaging.

Embedding the cover glass guarantees an even focal plane which is a prerequisite for all high-speed and high-resolution microscopy applications. Furthermore, the black upper housing reduces cross talk between adjacent wells during fluorescence microscopy and the optical glass, which exhibits virtually no autofluorescence, allows for maximum spectral transmission without depolarisation of transmitted light.

- / 10 wells with alphanumeric coding
- / Black detachable compartmentalization
- / Well distance is equal to a 96 well microplate
- / Positioning notch for automated microscopy
- / Highly transparent achromatic borosilicate glass bottom



CELLview Slide

- / 10 wells with alphanumeric coding
- / Positioning notch for automated microscopy
- / Black detachable compartmentalization
- / Reduced meniscus effect due to round well design
- / Glass thickness: 0.17 mm







FREE OF detectable RNase





Well format: 10, Height: 12.6 mm, Length: 75 mm, Width: 25 mm, Growth area / unit: 34 mm², Bottom: glass, Total volume (well): 440 µl

Item no.	Surface treatment	Sterile	Qty. inner / outer
543078	TC	+	1/20
543079	TC	+	5 / 45
543978	Advanced TC	+	1/20
543979	Advanced TC	+	5 / 45

[/] Sample packs are available on request.





/ For plates with an optical cycloolefin film bottom, please refer to our SCREENSTAR plates in this chapter.

CELLVIEW PLATE

CELL CULTURE PLATE WITH GLASS BOTTOM

CELLview glass bottom microplates are designed for demanding and high-resolution microscopic applications.

They consist of a cycloole-fin-based black frame with a 0.17 mm thin borosilicate glass bottom providing superior images of in-vitro cultures. The optimised microplate geometry and

the recessed bottom facilitate imaging of all peripheral wells even with immersion objectives. The round conical well design reduces the meniscus effect in order to assure equal cellular distribution and constant imaging results. An appropriate surface treatment improves cellular attachment and growth.

- / Outstanding image quality and resolution
- / Cycloolefin-based frame with 0.17 mm glass bottom
- / Excellent optical transparency
- / Recessed well bottom facilitating the use of lenses with low working distance and high aperture
- / Compatible with advanced confocal microscopic systems



CELLview Plate

- / For outstanding image quality and resolution
- / Recessed well bottom facilitating the use of objectives with low working distance
- / Ditch at the perimeter can be filled with liquid to minimise edge effects and evaporation
- / Compatible with advanced automated microscopic systems

Well format: 96, Growth area / unit: 33 mm², Well profile: F-bottom / Chimney Well, Bottom: glass, Raw material: COP, Working volume (well): 25 µl - 440 µl, Lid: yes

Item no.	Surface treatment	Product colour	Sterile	Qty. inner / outer
655891	TC	● black	+	1 / 16
655981	Advanced TC	● black	+	1 / 16





/ Forum No. 15: SCREENSTAR: A 1536 Well Microplate for High-Content and High-Throughput Screening (F073120)

SCREENSTAR MICROPLATES

SCREENSTAR are specialised microplates for sophisticated microscopic applications, in high content screening (HCS) or high-resolution microscopy with water and oil immersion objectives.

They combine outstanding glass-like optical properties with an excellent surface for adherent cell culture.

Moreover, the plates display ex-

cellent optical properties with reduced autofluorescence in the lower UV range, low birefringence and a refractive index of 1.53 comparable to glass. SCREENSTAR microplates enable complete periphery access for high magnification objectives. They are entirely manufactured out of cycloolefin with a black pigmented cycloolefin frame and a 190 μ m ultraclear cycloolefin film bottom.

- / 96 / 384 / 1536 well format
- For sophisticated microscopic applications and high content screening
- / Highly transparent cycloolefin film bottom
- / Adherent TC surface treatment



SCREENSTAR Microplates 96 / 384 / 1536 Well

- / For sophisticated microscopic applications and high content screening
- / Highly transparent cycloolefin film bottom
- / Adherent TC surface treatment













Bottom: Cycloolefin film, Raw material: COP, Surface treatment: TC

Item no.	Well format	Growth area / unit	Well profile	Product colour	Working volume (well)	Lid	Sterile	Qty. inner / outer
655866	96	33 mm²	F-bottom / Chimney Well	● black	25 μΙ - 440 μΙ	yes	+	1 / 16
781866	384	8.1 mm ²	F-bottom	black	10 μΙ - 110 μΙ	yes	+	8 / 32
789866	1536	2.1 mm ²	F-bottom	black	3 μΙ - 15 μΙ	no	+	17 / 68

/ Forum No. 15: SCREENSTAR: A 1536 Well Microplate for High-Content and High-Throughput Screening (F073120)





/ For further information, please refer to our website: www.gbo.com



SENSOPLATE

GLASS BOTTOM MICROPLATES

The research of biomolecular processes on the level of single molecules and in volume ranges equivalent to the size of a single bacterium is of immense importance, both in basic research and in industrial high-throughput screening. The combination of modern confocal optics, new fluorescent dyes, sensitive photomultipliers and improved data processing has revolutionised the technique of fluorescence correlation spectroscopy (FCS). Over the past few years this has led to its widespread application, and alongside the technological advances in hardware development, Greiner Bio-One worked hand-in-hand with customers and instrument suppliers to develop the glass bottom microplates. These better satisfy the requirements of fluorescence correlation spectroscopy with regards to optical clarity and deformation when compared to standard polystyrene plates.

The SensoPlate family was developed in a complete product line consisting of 24, 96, 384 and 1536 well glass bottom formats.

- / For fluorescence correlation spectroscopy and microscopic applications
- / 24 / 96 / 384 / 1536 well format
- / Black frame with highly transparent glass bottom
- / Glass bottom thickness of 175 µm is equivalent to the light path of standard coverslips
- / Sterile, with lid and single-packed



SensoPlate Glass Bottom Plates 24 / 96 / 384 / 1536 Well

- / For fluorescence correlation spectroscopy and microscopic applications
- / 24 / 96 / 384 / 1536 well format
- / Black frame with highly transparent glass bottom
- / Glass bottom thickness of 175 μm is equivalent to the light path of standard coverslips
- / Sterile, with lid and single-packed

Well profile: F-bottom, Bottom: glass, Raw material: PS, Surface treatment: untreated, Lid: yes


Item no.	Well format	Product colour	Plate geometry	Working volume (well)	Sterile	Qty. inner / outer
662892	24	● black		0.5 ml - 1.5 ml	+	1 / 12
655892	96	black		25 μΙ - 340 μΙ	+	1 / 16
781892	384	black		10 μΙ - 130 μΙ	+	1 / 16
783892	1536	● black	LoBase	3 μΙ - 10 μΙ	+	1 / 16



Greiner Bio-One offers a comprehensive range of tubes and multipurpose beakers for various applications. The vessels are available in different materials, with and without caps and with labelling options.

TUBES / MULTIPURPOSE BEAKERS

/	Tubes without closures116
	Tube Polystyrene117
	Tube Polypropylene117
	Grip Stopper118
	Tube Polypropylene for Storage Box118
/	Tubes with closures119
	Centrifuge Tube Polypropylene - 15 ml 120
	Centrifuge Tube Polypropylene - 50 ml121
	CELLreactor122
	Tubes with closures Polystyrene 122
	Tube Two-position cap123
/	Separation Tubes124
	Leucosep 12 ml and 50 ml 125
	OncoQuick® 125
/	Multipurpose Containers / Beakers 126
	Multipurpose Containers Polystyrene127
	Multipurpose Beaker Polypropylene 128
	Containers for Plant Cultures Polystyrene 128
	Drosophila Containers
	Ceaprene stopper

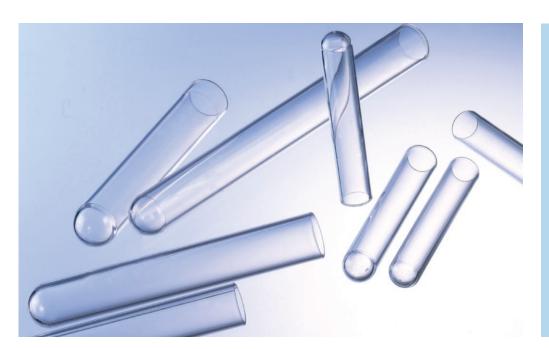



/ For the max. relative centrifugal force (RCF) and the chemical/thermal resistance of our tubes please refer to the technical appendix.

TUBES WITHOUT CLOSURES

The range of Greiner Bio-One tubes is very versatile and meets a wide variety of different demands.

Our tubes are made of different materials: **Polystyrene (PS)** is ideally suited for optical measurements as a result of its high clarity.

Polypropylene (PP) is recommended for the storage of chemical and biological samples due to its thermal, mechanical and chemical resistance. Greiner Bio-One tubes without closure are available with round bottom. In addition, it is possible to order appropriate closures in the form of grip stoppers and screw caps.

- / With round or conical bottom
- / With or without skirt
- / Available in polystyrene or polypropylene



Tube Polystyrene

- / High transparency
- / Available in different sizes and packaging units

Raw material: PS, Bottom shape: round

Item no.	Height	Ø	Working volume	Nominal volume	Oty. inner / outer
112101	55 mm	12 mm	≤3 ml	4 ml	240 / 3,600
115101	75 mm	12 mm	≤4 mI	5 ml	250 / 2,000
136101	100 mm	14 mm	≤8 ml	10 ml	1,400 / 1,400
160101	100 mm	16 mm	≤10.5 ml	12 ml	1,600 / 1,600
187101	100 mm	17 mm	≤12 ml	14 mI	1,500 / 1,500
169101	152 mm	16 mm	≤18 ml	20 ml	1,500 / 1,500



Tube Polypropylene

- / Suited for sample storage
- / Thermal, chemical and mechanical stability
- / Available in different sizes and packaging units

Raw material: PP, Bottom shape: round

Item no.	Height	Ø	Working volume	Nominal volume	Qty. inner / outer
112201	55 mm	12 mm	≤3 ml	4 ml	240 / 3,600
115201	75 mm	12 mm	≤4 mI	5 ml	250 / 2,000
160201	100 mm	16 mm	≤10 mI	12 ml	-/1,600
187201	100 mm	17 mm	≤12 mI	14 ml	1,500 / 1,500



Grip Stopper

/ Made of high-grade polyethylene

Raw material: PE

Item no.	Description	Product colour	Suitable for tubes Ø	Qty. inner / outer
303321	grip stopper	○ natural	12 mm	500/20,000
307321	grip stopper	natural	14 mm	100 / 10,000
310321	grip stopper	○ natural	16 mm	100 / 8,000
318321	grip stopper	natural	17 mm	250 / 8,000
724170	screw cap	○ natural	61 mm	720 / 720



Tube Polypropylene for Storage Box

/ Tubes for storage box item no. 975502

Height: 44 mm, Ø: 8.5 mm, Raw material: PP, Bottom shape: round, Working volume: ≤1 ml, Nominal volume: 1.3 ml

Item no.	Feature	Sterile	Qty. inner / outer
102201	tube		1,000 / 1,000
102261	tube chain with attached strip cap	+	-/1,000
102270	tube chain with attached strip cap		-/1,000



/ For the max. relative centrifugal force (RCF) and the chemical/thermal resistance of our tubes please refer to the technical appendix.



TUBES WITH CLOSURES

Greiner Bio-One tubes with closures are made of different materials: Polystyrene (PS) is ideally suited for optical measurements as a result of its high clarity. Polypropylene (PP) is recommended for the storage of chemical and biological samples due to its thermal, mechanical and chemical resistance. Greiner Bio-One tubes without

closure are available with round or with conical bottom as well as with or without skirt. They are available in sterile or non-sterile versions. The product range is completed by tubes with a two-position cap. This special capenables ventilation or an airtight closure of the tube, depending on the position of the stopper.



Centrifuge Tube Polypropylene - 15 ml

- / Optimal mechanical, thermal and chemical stability
- / Available with blue or white cap
- / Triple-packed options available
- / Light protection tubes for light-sensitive materials and reactions



FREE OF detectable DNase FREE OF detectable numan DNA

FREE OF detectable RNase





Height: 120 mm, Ø: 17 mm, Support skirt: no, Graduation: yes, Writing field: yes, Raw material: PP, Bottom shape: conical, Working volume: ≤14 ml, Nominal volume: 15 ml, Cap design: screw cap

Item no.	Feature	Product colour	Cap colour	Colour writing field	Graduation colour	Triple- packed	Sterile	Qty. inner / outer
188261-N		○natural	blue	○white	blue		+	50/500
188271-N		natural	blue	○ white	blue		+	50 / 1.000
188261		○natural	blue	○white	blue		+	50 / 500
188271		natural	blue	○ white	blue		+	100 / 1,000
188271-TRI		○natural	blue	○white	blue	yes	+	5/500
188283	light protection tube	brown	blue	○white	blue		+	50 / 500
188281		○ natural	○ white	○white	blue		+	50 / 500
188285		natural	O white	○white	blue		+	100 / 700
188280	light protection tube	brown	blue	○white	blue		+	100 / 1.000

^{/ 188261-}N &188721-N currently only available in Germany!



Centrifuge Tube Polypropylene - 50 ml

- / Optimal mechanical, thermal and chemical stability
- / Available with blue or white cap
- / Triple-packed options available
- / Light protection tubes for light-sensitive materials and reactions



FREE OF detectable DNase FREE OF detectable human DNA

FREE OF detectable RNase





Height: 115 mm, Ø: 30 mm, Graduation: yes, Writing field: yes, Raw material: PP, Bottom shape: conical, Working volume: ≤50 ml, Nominal volume: 50 ml, Cap design: screw cap

Item no.	Feature	Support skirt	Product colour	Cap colour	Colour writing field	Graduation colour	Triple- packed	Sterile	Oty. inner/ outer
210261		yes	○ natural	blue	○white	blue		+	25 / 450
210270		yes	onatural	blue	○white	blue		+	25 / 300
227261		no	○ natural	blue	○white	blue		+	20 / 500
227270		no	onatural	blue	○white	blue		+	25 / 300
227261-TRI		no	○ natural	blue	○white	blue	yes	+	5 / 250
227285		no	onatural	O white	○white	blue		+	20 / 500
227281		no	○ natural	○ white	○white	blue		+	25 / 300
227280	light protection tube	no	brown	blue	○white	blue		+	20 / 500
227283	light protection tube	no	brown	blue	○white	blue		+	25 / 300



CELLreactor

15 ml and 50 ml polypropylene tube with filter screw cap

- / For cultivation of suspension cells and expansion of aerobic microorganisms
- / Facilitates a high number of parallel experiments
- / Maximal sterility and excellent gas exchange
- Conical tube design and in-tube harvest

STERILE FREE OF detectable DNase

FREE OF detectable human DNA

FREE OF detectable RNase





Graduation: yes, Writing field: yes, Raw material: PP, Bottom shape: conical, Cap design: filter screw cap

Item no.	Height	Ø	Cap colour	Working volume	Nominal volume	Sterile	Qty. inner / outer
188241	120 mm	17 mm	blue	1 ml - 5 ml	15 mI	+	20 / 500
227245	115 mm	30 mm	blue	1 ml - 35 ml	50 ml	+	20 / 500

- $\textit{I} \quad \text{Application Note: Cultivation of Suspension and Hybridoma Cells in CELLSTAR} \text{ } \textit{CELL} \textit{reactor Tubes} (F073918) \\$
- / Application Note: Superior protein yields in suspension CHO cells using FectoPRO™-mediated transient transfection in CELLSTAR® CELLreactor(F073926)



Tubes with closures Polystyrene

- / High transparency
- / Available with screw cap or bayonet cap in blue / black / white
- / With or without skirt
- / In bulk or rack packaging

Raw material: PS

Item no.	Height	Ø	Sup- port skirt	Grad- uation	Writing field	Bottom shape	Cap colour	Working volume	Nomi- nal volume	Cap design	Sterile	Oty. inner / outer
164180	100 mm	16.8 mm	yes	_	no	coni- cal	O white	≤12 ml	12 ml	bayonet cap		1,300 / 1,300
164161	100 mm	16.8 mm	yes	-	no	coni- cal	blue	≤12 ml	12 ml	bayonet cap	+	25 / 1,000
163177	100 mm	17 mm	no	-	no	round	• black	≤12 ml	12 ml	screw cap		900 / 900

Item no.	Height	Ø	Sup- port skirt	Grad- uation	Writing field	Bottom shape	Cap colour	Working volume	Nomi- nal volume	Cap design	Sterile	Oty. inner / outer
188161	120 mm	17 mm	no	yes	yes	coni- cal	blue	≤14 ml	15 ml	screw cap	+	50 / 500
188171	120 mm	17 mm	no	yes	yes	coni- cal	blue	≤14 ml	15 ml	screw cap	+	100 / 1,000
186171	120 mm	17 mm	no	-	no	round	blue	≤15 ml	15 ml	screw cap	+	100 / 1,000

 $\textit{I} \quad \text{Item no. 163177 is also available as polypropylene version with white screw cap (item no. 163270)}.$



Tube

Two-position cap

The two-position cap enables ventilation or an airtight closure of the tube, depending on the position of the stopper.

Bottom shape: round, Cap design: Two-position cap

Item no.	Height	Ø	Gradua- tion	Writing field	Raw material	Graduation colour	Working volume	Nominal volume	Sterile	Qty. inner / outer
115261	75 mm	12 mm	yes	no	PP		≤4 ml	5 ml	+	1 / 1,000
115262	75 mm	12 mm	yes	no	PP		≤4 ml	5 ml	+	25 / 2,000
120161	75 mm	12.4 mm	-	no	PS	blue	≤4 ml	4.5 ml	+	1 / 1,000
120180	75 mm	12.4 mm	-	no	PS	blue	≤4 ml	4.5 ml	+	25 / 2,000
187261	95 mm	18 mm	yes	yes	PP		≤12 ml	14 ml	+	1/800
187262	95 mm	18 mm	yes	yes	PP		≤12 ml	14 ml	+	25 / 1,000
191161	95 mm	18 mm	-	no	PS		≤12.5 ml	14 ml	+	1 / 750
191180	95 mm	18 mm	-	no	PS		≤12.5 ml	14 ml	+	25 / 1,000





/ For further information, please refer to our website: www.gbo.com



SEPARATION TUBES

Different separation techniques can be used to enrich certain particles (DNA, RNA, proteins, cells etc.) specifically from complex biological mixtures such as cell and tissue homogenates, blood or urine, so that they can then be selectively investigated. The separation of cells by density gradient centrifugation has proven to be the most often used method.

Leucosep has been developed for optimal separation of lymphocytes and peripheral mononuclear cells (so-called PB-MCs) from human whole blood and bone marrow by means of density gradient centrifugation.

The key feature of Leucosep is the porous barrier incorporated into the centrifuge tube made of highly translucent polypropylene.

OncoQuick® is a simple-to-use, rapid and efficient system for the enrichment of circulating tumour cells that are released into the blood by a solid epithelial tumour or malignant melanoma. OncoQuick® combines the advantages of cell separation by density gradient centrifugation with recovery rates that are comparable with immunobead methods.

- / Leucosep for the separation of lymphocytes/ mononuclear cells from blood/bone marrow
- / OncoQuick® for the enrichment of disseminated, circulating tumour cells from peripheral blood



Leucosep 12 ml and 50 ml

Efficient separation of lymphocytes and mononuclear cells from peripheral blood and bone marrow.

- / Easy to fill
- / No re-contamination with unwanted red blood cells
- / PBMC isolation from whole blood in 15 min

Type of sample: blood

Item no.	Separation medium	Bottom shape	Sample volume	Sterile	Qty. inner / outer
163288	pre-filled with Leucosep separation medium	round	3 - 8 ml		50 / 500
163289		round	3 - 8 mI		50/500
163290		round	3 - 8 mI	+	50/500
227288	pre-filled with Leucosep separation medium	conical	15 - 30 ml		25 / 250
227289		conical	15 - 30 ml		25 / 300
227290		conical	15 - 30 ml	+	25 / 300

/ For IFU please refer to our website: www.gbo.com



OncoQuick®

Enrichment of disseminated, circulating tumour cells from peripheral blood.

- / Time request approx.45 minutes
- / Reproducible recovery: > 70 %
- / No additional laboratory equipment required
- / Tubes are pre-filled with separation medium

Sample volume: 15 - 30 ml, Type of sample: blood

Item no.	Cap colour	Qty. inner / outer
227255	blue	1/1
227250	blue	10 / 10

/ For IFU please refer to our website: www.gbo.com





MULTIPURPOSE CONTAINERS / BEAKERS

Multipurpose containers or beakers can be used for a variety of different applications. They can be applied universally for academic and non-human medical purposes. Multipurpose containers are made of polystyrene or polypropylene and are available in different sizes. They feature either plastic or metal caps and can be supplied with or without plain or printed label.

Containers for plant culture are made of a very clear material which ensures maximum light transmission and thus rapid

and successful growth of plant cultures. These culture containers are not only suitable for the proliferation of plant cultures but can also be used as transport containers and are available with or without lids. Our drosophila containers are ideally suited for the cultivation of Drosophila melanogaster. The bottom part of the container is made of polystyrene. Drosophila containers can be supplied with a separately orderable ceaprene stopper. This stopper is gas-permeable and made of water-repellent material.

- / Multipurpose containers in polystyrene or polypropylene
- / Polystyrene containers for plant culture
- / Polystyrene drosophila containers



Multipurpose Containers Polystyrene

- / Manufactured from crystal clear polystyrene
- / Available in different sizes
- Can be used universally for academic and non-human medical purposes
- / With or without plain or printed label
- / Plastic or metal cap

Raw material: PS, Cap design: screw cap

Item no.	Feature	Height	Ø	Support skirt	Bottom shape	Cap colour	Working volume	Nominal volume	Cap material	Sterile	Oty. inner/ outer
201150		91 mm	25 mm	yes	conical	○white	≤23 ml	30 ml	plastics		400 / 400
201170		91 mm	25 mm	yes	conical	○white	≤23 ml	30 ml	plastics	+	400 / 400
201152	with plain label	91 mm	25 mm	yes	conical	○white	≤23 ml	30 ml	plastics		400 / 400
201172	with plain label	91 mm	25 mm	yes	conical	○white	≤23 ml	30 ml	plastics	+	400 / 400
224170		77 mm	49 mm			⊚silver	≤80 ml	100 ml	metal		20 / 160
225170		115 mm	49 mm			○silver	≤150 ml	150 ml	metal		20 / 120
225180	with plain label	115 mm	49 mm			○silver	≤150 ml	150 ml	metal		20 / 120



Multipurpose Beaker Polypropylene

- / Can be used universally for non human medical purposes
- / Optimal mechanical, thermal and chemical stability



Height: 70 mm, Ø: 55 mm, Graduation: yes, Working volume: ≤120 ml, Nominal volume: 120 ml, Cap design: screw cap

Item no.	Feature	Product colour	Cap colour	Sterile	Oty. inner / outer
724411		○ natural	○white		-/300
724412		natural	○white	+	-/300
724413	Tamper proof label	○ natural	○white	+	-/300
724414		natural	● red	+	-/300

/ Item No., lot number and expiry date are printed on the screw cap of the beaker to provide additional information and improve traceability



Containers for Plant Cultures Polystyrene

The use of a very clear material ensures maximum light transmission and thus rapid growth. These culture containers are not only suitable for the proliferation of plant cultures but can also be used as transport containers.

Height: 100 mm, Raw material: PS

Item no.	Description	Ø	Working volume	Nominal Cap design		Sterile	Qty. inner / outer
960161	container	53 mm	≤150 mI	175 ml	pull cap	+	4/300
968177	container, bottom part	68 mm	≤300 ml	330 ml			192 / 192
968162	container	68 mm	≤300 ml	330 ml	pull cap	+	1 / 168

/ Also available with lid as sterile version.



Drosophila Containers

- / Ideally suited for the cultivation of Drosophila melanogaster
- / Choice of different sizes
- / Bottom part of container made of polystyrene

Description: Drosophila Containers, Raw material: PS

Item no.	Height	Ø	Working volume	Nominal volume	Qty. inner / outer
205101	64 mm	27 mm		28 ml	1,500 / 1,500
217101	82 mm	36 mm		68 ml	605 / 605
960177	100 mm	53 mm	≤150 ml	175 ml	315 / 315

/ Can be supplied with a separately orderable ceaprene stopper. This stopper is gas-permeable and made of water-repellent material.



Ceaprene stopper

- / gas-permeable
- / made of water-repellent material

Item no.	Suitable for tubes Ø	Oty. inner / outer
323070	22 mm	- /1,500
354070	27 mm	- /1,500
330070	36 mm	-/605
332070	53 mm	-/315



For sample freezing and storage, Greiner Bio-One offers a comprehensive product portfolio including Cryo.s tubes, racks as well as electronic equipment for tube closure and sample tracking. With the brand Cryo.s Greiner Bio-One combines more than 30 years of experience in the manufacturing of cryogenic tubes with latest technology and innovation.

CRYOTECHNICS

/	Cryo.s Cryo Tubes	132
	Cryo.s - 1 ml	133
	Cryo.s - 2 ml	133
	Cryo.s - 4 ml	135
	Cryo.s-5 ml	135
	Support Rack	136
	Cryo storage box	136
/	Cryo.s with Barcode	137
	Cryo.s with Barcode	
	— Preproduced unique codes	138
	Cryo.s with Barcode	
	— Customised Codes	138
	48-way Datamatrix Cryo Rack	139
	81-way Datamatrix Cryo Rack	140
/	Cryo.s Biobanking Tubes	14
	Cryo.s Biobanking Tubes 300 µl	142
	Cryo.s Biobanking Tubes 600 µl	143
	Cryo.s Biobanking Tubes 1000 µl	144
	Screw Caps Biobanking Tubes	145
	96-way Datamatrix Cryo Rack	145
	Cryo.s Decapper	146
	Cryo's Rack Scanner	14 F





/ For a freezing and thawing protocol as well as safety advisory for Cryo.s please refer to the technical appendix.

CRYO.S CRYO TUBES

Cryo.s, Cryo.s with datamatrix and linear barcode and Cryo.s biobanking tubes are for storage of tissue, cells, fungi, bacteria, spores, cellular extracts or body fluids at ultra-low temperature for research and development purposes as well as in-vitro diagnostics. They must not be stored in the liquid phase of liquid nitrogen, but only in the gas phase above. Cryo.s tubes are not intended for any application in the context of reproductive medicine.

An outstanding feature of all Cryo.s is the utilisation of a USP class VI certified, medical grade polypropylene for tube manufacturing. This material does not release significant amounts of leachables, hence stored samples stay clean and free of contamination. Sterile product versions are sterilised applying an ISO 11137 validated irradiation procedure yielding an SAL level of 10^{-6} .

The portfolio of Cryo.s tubes includes cryogenic tubes with different volumes, different base forms as well as several cap colours. Cryo.s offer several features for the identification and labelling of individual samples, to include coloured screw caps, a white, scratch-resistant writing area and barcoding options.

- / CE-marked
- USP class VI certified medical grade polypropylene
- Sterilised applying an ISO 11137 validated irradiation procedure yielding an SAL level of 10 -6
- / Suited for storage in liquid nitrogen vapour phase (-196 °C)



Cryo.s - 1 ml

- / High thermal resistance
- / Cryo.s with internal thread have a screw cap with silicone gasket
- / Cap inserts item no. 304134 (50 pieces per bag)
- / USP class VI certified medical grade polypropylene







FREE OF detectable DNase FREE OF detectable human DN







Height: 42 mm, Ø: 12.5 mm, Starfoot: yes, Raw material: PP, Bottom shape: conical, Working volume: ≤1.2 ml, Cap design: screw cap

Item no.	Barcode	Graduation	Writing field	Cap colour	Thread type	Triple-packed	Sterile	Qty. inner / outer
123261	no	no	no	○ natural	internal		+	100 / 500
123263	no	yes	yes	natural	internal		+	100 / 500
123263-TRI	no	yes	yes	○ natural	internal	yes	+	10 / 200
123277	no	yes	yes	● green	internal		+	100 / 500
123278	no	yes	yes	yellow	internal		+	100 / 500
123279	no	yes	yes	blue	internal		+	100 / 500
123280	no	yes	yes	● red	internal		+	100 / 500

/ with 150 inserts per case



Cryo.s - 2 ml

- / High thermal resistance
- / Cryo.s with internal thread have a screw cap with silicone gasket
- / Cap inserts item no. 304134 (50 pieces per bag)
- / USP class VI certified medical grade polypropylene

FREE OF detectable human DNA RNase STERILE DNase FREE OF detectable RNase cytotoxic pyrogeni

Raw material: PP, Bottom shape: round, Cap design: screw cap

Item no.	Height	Ø	Star- foot	Bar- code	Gradu- ation	Writ- ing field	Cap colour	Work- ing volume	Thread type	Tri- ple-packed	Ster- ile	Oty. inner/ outer
121261	48 mm	12.5 mm	no	no	-	no	○ natural	≤2 ml	internal		+	100 / 500

Item no.	Height	Ø	Star- foot	Bar- code	Gradu- ation	Writ- ing field	Cap colour	Work- ing volume	Thread type	Tri- ple-packed	Ster- ile	Oty. inner/ outer
121263	48 mm	12.5 mm	no	no	yes	yes	onatural	≤2 ml	internal		+	100 / 500
121277	48 mm	12.5 mm	no	no	yes	yes	green	≤2 ml	internal		+	100 / 500
121278	48 mm	12.5 mm	no	no	yes	yes	yellow	≤2 ml	internal		+	100 / 500
121279	48 mm	12.5 mm	no	no	yes	yes	blue	≤2 ml	internal		+	100 / 500
121280	48 mm	12.5 mm	no	no	yes	yes	● red	≤2 ml	internal		+	100 / 500
122261	48 mm	12.5 mm	yes	no	-	no	○ natural	≤2 ml	internal		+	100 / 500
122263	48 mm	12.5 mm	yes	no	yes	yes	onatural	≤2 ml	internal		+	100 / 500
122263-TRI	48 mm	12.5 mm	yes	no	yes	yes	○ natural	≤2 ml	internal	yes	+	10 / 200
122277	48 mm	12.5 mm	yes	no	yes	yes	green	≤2 ml	internal		+	100 / 500
122278	48 mm	12.5 mm	yes	no	yes	yes	yellow	≤2 ml	internal		+	100 / 500
122279	48 mm	12.5 mm	yes	no	yes	yes	blue	≤2 ml	internal		+	100 / 500
122280	48 mm	12.5 mm	yes	no	yes	yes	● red	≤2 ml	internal		+	100 / 500
126261	47 mm	12.4 mm	yes	no	-	no	onatural	≤2.2 ml	exter- nal		+	100 / 500
126263	47 mm	12.4 mm	yes	no	yes	yes	○ natural	≤2.2 ml	exter- nal		+	100 / 500
126263-TRI	47 mm	12.4 mm	yes	no	yes	yes	natural	≤2.2 ml	exter- nal	yes	+	10 / 200
126277	47 mm	12.4 mm	yes	no	yes	yes	green	≤2.2 ml	exter- nal		+	100 / 500
126278	47 mm	12.4 mm	yes	no	yes	yes	yellow	≤2.2 ml	exter- nal		+	100 / 500
126279	47 mm	12.4 mm	yes	no	yes	yes	blue	≤2.2 ml	exter- nal		+	100 / 500
126280	47 mm	12.4 mm	yes	no	yes	yes	● red	≤2.2 ml	exter- nal		+	100 / 500

/ with 150 inserts per case



Cryo.s - 4 ml

- / High thermal resistance
- / Cap inserts item no. 304134 (50 pieces per bag)
- / USP class VI certified medical grade polypropylene







Description: Cryo.s 4 ml, Height: 83 mm, Ø: 12.4 mm, Starfoot: yes, Raw material: PP, Bottom shape: round, Working volume: ≤4 ml, Cap design: screw cap

Item no.	Barcode	Graduation	Writing field	Cap colour	Thread type	Sterile	Oty. inner / outer
127261	no	no	no	○ natural	external	+	50 / 300
127263	no	yes	yes	natural	external	+	50/300
127277	no	yes	yes	● green	external	+	50/300
127278	no	yes	yes	yellow	external	+	50/300
127279	no	yes	yes	blue	external	+	50/300
127280	no	yes	yes	● red	external	+	50/300

/ with 100 inserts per case



Cryo.s - 5 ml

- / High thermal resistance
- / Cryo.s with internal thread have a screw cap with silicone gasket
- / Cap inserts item no. 304134 (50 pieces per bag)
- / USP class VI certified medical grade polypropylene



Description: Cryo.s 5 ml, Height: 86 mm, Ø: 12.5 mm, Starfoot: no, Raw material: PP, Bottom shape: round, Working volume: ≤4.5 ml, Cap design: screw cap

Item no.	Barcode	Graduation	Writing field	Cap colour	Thread type	Sterile	Qty. inner / outer
124261	no	-	no	○ natural	internal	+	50/300
124263	no	yes	yes	natural	internal	+	50/300

Item no.	Barcode	Graduation	Writing field	Cap colour	Thread type	Sterile	Oty. inner / outer
124273	no	yes	yes	● red	internal	+	50/300
124275	no	yes	yes	● green	internal	+	50/300
124276	no	yes	yes	yellow	internal	+	50/300
124274	no	yes	yes	blue	internal	+	50/300

/ with 100 inserts per case



Support Rack

- / Suitable for Cryo.s with starfoot base (Item No. 122XXX, 123XXX, 126XXX, 127XXX)
- / Improved handling since the tubes can be opened with one hand
- / Rubber base to prevent slipping
- / Offers space for up to 40 Cryo.s

Description: support rack for one-hand operation, Height: 22 mm, Length: 200 mm, Width: 100 mm, Raw material: PC

Item no.	Rack colour	Oty. inner / outer
802501	blue	1/50



Cryo storage box

The box material polypropylene is very temperature and shock resistant, thus allowing storage temperatures as low as -90 °C.

- / Holds 81 Cryo.s sample tubes
- / Transparent lid for optimum visibility of box content

Description: Cryo storage box, Height: 51 mm, Length: 126.5 mm, Width: 126.5 mm, Raw material: PP

Item no.	Rack colour	Oty. inner / outer
802202	○ natural	5 / 20
802203	● red	5 / 20
802204	blue	5 / 20
802206	yellow	5 / 20
802225	● green	5 / 20

 $[\]emph{I} \quad \text{The indicated height refers to the Cryo Storage Box filled with 2 ml Cryo.s and covered with a lid.}$





/ For further information, please refer to our website: www.gbo.com



CRYO.S WITH BARCODE

Cryo.s, Cryo.s with datamatrix and linear barcode and Cryo.s biobanking tubes are for storage of tissue, cells, fungi, bacteria, spores, cellular extracts or body fluids at ultra-low temperature for research and development purposes as well as in-vitro diagnostics. They must not be stored in the liquid phase of liquid nitrogen, but only in the

gas phase above. Cryo.s tubes are not intended for any application in the context of reproductive medicine.

With Cryo.s with Datamatrix and the accessory racks in 48-way and 81-way formats, Greiner Bio-One expands its portfolioby optimum solutions for semi-automated and automated sample handling and storage.

- / Datamatrix code ECC200 on the tube bottom with reliable Reed-Solomon error correction
- / Linear barcode type
 128 with plain text on the tube side
- Application of barcodes and datamatrix codes by means of innovative laser technology
- / 100 % controlled code readability
- / Suited for storage in liquid nitrogen vapour phase (-196 °C)
- / Coded tubes are suitable for airfreight (IATA conformity)



Cryo.s with Barcode

- Preproduced unique codes

- / Pre-produced unique datamatrix code ECC 200 with 14 x 14 dots on tube bottom, additional linear barcode type 128 with human readable text on tube side
- / Highest code resistance against chemicals, mechanical stress and extremely low storage temperatures (down to -196 °C)









FREE OF detectable human DNA FREE OF detectable RNase





Starfoot: yes, Barcode type: linear barcode and datamatrix code, Raw material: PP, Cap design: screw cap

Item no.	Height	Ø	Barcode	Bottom shape	Cap colour	Working volume	Thread type	Sterile	Qty. inner / outer
123263-2DG	42 mm	12.5 mm	yes	conical	O natural	≤1.2 ml	internal	+	100 / 500
122263-2DG	48 mm	12.5 mm	yes	round	o natural	≤2 ml	internal	+	100 / 500
126263-2DG	47 mm	12.4 mm	yes	round	O natural	≤2.2 ml	external	+	100 / 500
127263-2DG	83 mm	12.4 mm	yes	round	o natural	≤4 ml	external	+	50 / 300

 ${\it I} \quad \hbox{Other screw cap colours on request.}$



Cryo.s with Barcode

— Customised Codes

- / Customised datamatrix code ECC 200 with 14 x 14 dots on tube bottom, additional customised linear barcode type 128 with human readable text on tube side
- / Highest code resistance against chemicals, mechanical stress and extremely low storage temperatures (down to -196 °C)







Raw material: PP, Cap design: screw cap

Item no.	Height	Ø	Star- foot	Bar- code	Barcode type	Bottom shape	Cap colour	Work- ing vol- ume	Nomi- nal volume	Thread type	Ster- ile	Oty. inner/ outer
123263-2D3	42 mm	12.5 mm	yes	yes	linear bar- code and datamatrix code	coni- cal	○ natural	≤1.2 ml	1 ml	inter- nal	+	100 / 500

Item no.	Height	Ø	Star- foot	Bar- code	Barcode type	Bottom shape	Cap colour	Work- ing vol- ume	Nomi- nal volume	Thread type	Ster- ile	Oty. inner/ outer
122263-2D3	48 mm	12.5 mm	yes	yes	linear bar- code and datamatrix code	round	○natural	≤2 ml	2 ml	inter- nal	+	100 / 500
126263-2D1	47 mm	12.4 mm	yes	yes	linear bar- code and datamatrix code	round	○ natural	≤2.2 ml	2 ml	ex- ter- nal	+	100 / 500
121263-128	48 mm	12.5 mm	no	yes	linear barcode	round	natural	≤2 ml	2 ml	inter- nal	+	100 / 500
127263-2D1	83 mm	12.4 mm	yes	yes	linear bar- code and datamatrix code	round	○natural	≤4 ml	4 ml	ex- ter- nal	+	50 / 300
124263-128	86 mm	12.5 mm	no	yes	linear barcode	round	onatural	≤4.5 ml	5 ml	inter- nal	+	50 / 300

- / Other screw cap colours on request.
- / Barcodes to be specified in order form F071004.



48-way Datamatrix Cryo Rack

- / SLAS-standard format
- / Automation compatible
- / Suitable for storage in vapour phase above liquid nitrogen
- / Barcoded versions on request

Length: 85.5 mm, Width: 127.8 mm, Raw material: PP

Item no.	Description	Height	Rack colour	Lid type	Suitable for Cryo.s	Qty. inner / outer
803277	low rack	30 mm	● black	without lid	1/2/4/5 ml	5 / 20
803202	low rack	52.5 mm	● black	with lid	1 / 2 ml	-/20
803270	high rack	88.5 mm	● black	with lid	4 ml	- /15



81-way Datamatrix Cryo Rack

- / Shock-proof and temperature-resistant polycarbonate
- / Stacking feature and rotation stoppers
- / 133 x 133 mm footprint for storage in classic freezer setups
- / Suitable for storage in vapour phase above liquid nitrogen
- / Barcoded versions on request

Length: 132.4 mm, Width: 132.4 mm, Raw material: PC, Lid type: with lid

Item no.	Description	Height	Rack colour	Suitable for Cryo.s	Qty. inner / outer
802576	low rack	52 mm	black	1 / 2 ml	1 / 10
802506	high rack	88.1 mm	black	4 ml	1/10





- / Forum No. 21: Sample Storage Tubes as Quality-Critical Components in Biobanking (F073072)
- / Watch our video:



CRYO.S BIOBANKING TUBES

With Cryo.s biobanking tubes, Greiner Bio-One offers an ideal solution for the efficient storage of biological samples in large-scale biorepositories.

Cryo.s biobanking tubes are available with working volumes of 235 μ l, 580 μ l and 975 μ l and offered in highly automation-friendly 96-way racks. The innovative design of tubes and racks allows for a very space-efficient storage with up to 30 % better utilisation of storage space in freezers or liquid nitrogen tanks.

Cryo.s, Cryo.s with datamatrix and linear barcode and Cryo.s biobanking tubes are for storage of tissue, cells, fungi, bacteria, spores, cellular extracts or body fluids at ultra-low temperature for research and development purposes as well as in-vitro diagnostics. They must not be stored in the liquid phase of liquid nitrogen, but only in the gas phase above. Cryo.s tubes are not intended for any application in the context of reproductive medicine.

- Application of barcodes and datamatrix codes by means of innovative laser technology
- / 100 % controlled code readability
- / Coded tubes are suitable for airfreight (IATA conformity)
- / Suited for storage at -20 °C, -80 °C and in vapour phase above liquid nitrogen



Cryo.s Biobanking Tubes 300 µl

- / Pre-produced unique datamatrix code on tube, datamatrix code and linear barcode on rack
- / Height-reduced screw cap conserving up to 30 % freezer space
- Customised code sequences on request (order form F071003)















Height: 18.7 mm, Total rack height: 21.6 mm, Ø: 8.8 mm, Barcode type: datamatrix code, Raw material: PP, Working volume: ≤235 μl

Item no.	Barcode	Cap colour	Cap design	Thread type	Type of packaging	Sterile	Qty. inner / outer
976570	yes		without screw cap	internal	rack		480 / 960
976580	yes	onatural	screw cap	internal	rack		480 / 960
976561	yes	○ natural	screw cap	internal	rack	+	480 / 960
976586	yes	green	screw cap	internal	rack		480 / 960
976566	yes	● green	screw cap	internal	rack	+	480 / 960
976585	yes	yellow	screw cap	internal	rack		480 / 960
976565	yes	yellow	screw cap	internal	rack	+	480 / 960
976584	yes	blue	screw cap	internal	rack		480 / 960
976564	yes	blue	screw cap	internal	rack	+	480 / 960
976583	yes	● red	screw cap	internal	rack		480 / 960
976563	yes	● red	screw cap	internal	rack	+	480 / 960
976588	yes	pink	screw cap	internal	rack		480 / 960
976568	yes	pink	screw cap	internal	rack	+	480 / 960
976589	yes	brown	screw cap	internal	rack		480/960
976569	yes	brown	screw cap	internal	rack	+	480 / 960
976587	yes	black	screw cap	internal	rack		480/960
976567	yes	● black	screw cap	internal	rack	+	480 / 960
131202	yes	natural	screw cap	internal	bulk		480 / 960
131263	yes	○natural	screw cap	internal	bulk	+	480 / 960

[/] Follow the instructions of use provided within each box.

[/] USP class VI certified medical grade polypropylene



Cryo.s Biobanking Tubes 600 µl

- / Pre-produced unique datamatrix code on tube, datamatrix code and linear barcode on rack
- / Height-reduced screw cap conserving up to 30 % freezer space
- Customised code sequences on request (order form F071003)







FREE OF detectable human DNA

FREE OF detectable RNase





Height: 33.3 mm, Total rack height: 36.2 mm, Ø: 8.8 mm, Barcode type: datamatrix code, Raw material: PP, Working volume: ≤580 µl

Item no.	Barcode	Cap colour	Cap design	Thread type	Type of packaging	Sterile	Qty. inner / outer
977570	yes		without screw cap	internal	rack		192 / 960
977580	yes	natural	screw cap	internal	rack		192 / 960
977561	yes	○natural	screw cap	internal	rack	+	192 / 960
977586	yes	● green	screw cap	internal	rack		192 / 960
977566	yes	■ green	screw cap	internal	rack	+	192 / 960
977585	yes	yellow	screw cap	internal	rack		192 / 960
977565	yes	yellow	screw cap	internal	rack	+	192 / 960
977584	yes	blue	screw cap	internal	rack		192 / 960
977564	yes	blue	screw cap	internal	rack	+	192 / 960
977583	yes	● red	screw cap	internal	rack		192 / 960
977563	yes	● red	screw cap	internal	rack	+	192 / 960
977588	yes	pink	screw cap	internal	rack		192 / 960
977568	yes	pink	screw cap	internal	rack	+	192 / 960
977589	yes	brown	screw cap	internal	rack		192 / 960
977569	yes	brown	screw cap	internal	rack	+	192 / 960
977587	yes	black	screw cap	internal	rack		192 / 960
977567	yes	● black	screw cap	internal	rack	+	192 / 960
132202	yes	natural	screw cap	internal	bulk		192 / 960
132263	yes	natural	screw cap	internal	bulk	+	192 / 960

[/] Follow the instructions of use provided within each box.

[/] USP class VI certified medical grade polypropylene



Cryo.s Biobanking Tubes 1000 µl

- / Pre-produced unique datamatrix code on tube, datamatrix code and linear barcode on rack
- Height-reduced screw cap conserving up to 30 %freezer space
- Customised code sequences on request (order form F071003)
- USP class VI certified medical grade polypropylene











Height: 50.8 mm, Total rack height: 53.9 mm, Ø: 8.8 mm, Barcode type: datamatrix code, Raw material: PP, Working volume: ≤975 μl

Item no.	Barcode	Cap colour	Cap design	Thread type	Type of packaging	Sterile	Qty. inner / outer
978570	yes		without screw cap	internal	rack		192 / 960
978580	yes	onatural	screw cap	internal	rack		192 / 960
978561	yes	○ natural	screw cap	internal	rack	+	192 / 960
978586	yes	● green	screw cap	internal	rack		192 / 960
978566	yes	● green	screw cap	internal	rack	+	192 / 960
978585	yes	yellow	screw cap	internal	rack		192 / 960
978565	yes	yellow	screw cap	internal	rack	+	192 / 960
978584	yes	blue	screw cap	internal	rack		192 / 960
978564	yes	blue	screw cap	internal	rack	+	192 / 960
978583	yes	● red	screw cap	internal	rack		192 / 960
978563	yes	● red	screw cap	internal	rack	+	192 / 960
978588	yes	pink	screw cap	internal	rack		192 / 960
978568	yes	pink	screw cap	internal	rack	+	192 / 960
978589	yes	brown	screw cap	internal	rack		192 / 960
978569	yes	brown	screw cap	internal	rack	+	192 / 960
978587	yes	black	screw cap	internal	rack		192 / 960
978567	yes	● black	screw cap	internal	rack	+	192 / 960
133202	yes	natural	screw cap	internal	bulk		192 / 960
133263	yes	○ natural	screw cap	internal	bulk	+	192 / 960

[/] Follow the instructions of use provided within each box.

[/] USP class VI certified medical grade polypropylene



Screw Caps Biobanking Tubes

- / 96 screw caps in cap carrier
- / Non-sterile







FREE OF detectable human DN*A* FREE OF detectable RNase





Description: 96 screw caps in cap carrier, Raw material: PP, Cap design: screw cap

Item no.	Barcode	Cap colour	Qty. inner / outer
385270	no	○ natural	960 / 960
385276	no	● green	960 / 960
385275	no	yellow	960 / 960
385274	no	blue	960 / 960
385273	no	● red	960 / 960
385278	no	pink	960 / 960
385279	no	● brown	960 / 960
385277	no	● black	960 / 960



96-way Datamatrix Cryo Rack

- / Pre-produced unique datamatrix code ECC 200 and linear barcode type 128
- / Made from polycarbonate















Barcode type: linear barcode and datamatrix code, Raw material: PC

Item no.	Height	Barcode	Rack colour	Suitable for Cryo.s	Qty. inner / outer
976501	19.1 mm	yes	● black	300 µl	5 / 10
977501	33.7 mm	yes	● black	600 µI	2 / 10
978501	51.4 mm	yes	● black	1000 μΙ	2 / 10



Cryo.s Decapper

The new Cryo.s Decapper enables the precise simultaneous de- and recapping of cryotubes.

- / Individual motors for precise application of defined torque values
- / 6 channel version for 1, 2 and 4 ml Cryo.s in 48 way cryo racks
- / 8 channel version for Cryo.s Biobanking Tubes
- / Docking station with charging function included

Description: Cryo.s 8-channel handheld decapper, Content kit: 1 Decapper, Power supply unit, Docking station with charging function, Quick start guide

Item no.	Qty. inner / outer
852070	1/1



Cryo.s Rack Scanner


With the Cryo.s rack scanner, Greiner Bio-One provides an optimum solution for scanning Cryo.s with datamatrix and Cryo.s biobanking tubes in SLAS-standard formatted racks.

- / Single tube scan feature
- / Small footprint
- / 5 year warranty

Height: 7 cm, Length: 31 cm, Width: 19.5 cm, Content kit: 1 scanner, power cable, USB cable, driver, user manual

Item no.	Description	Qty. inner / outer
849070	rack scanner EU/UK/Asia edition	1/1
849050	rack scanner US edition	1/1

CryotechnicsCryo.s Biobanking Tubes

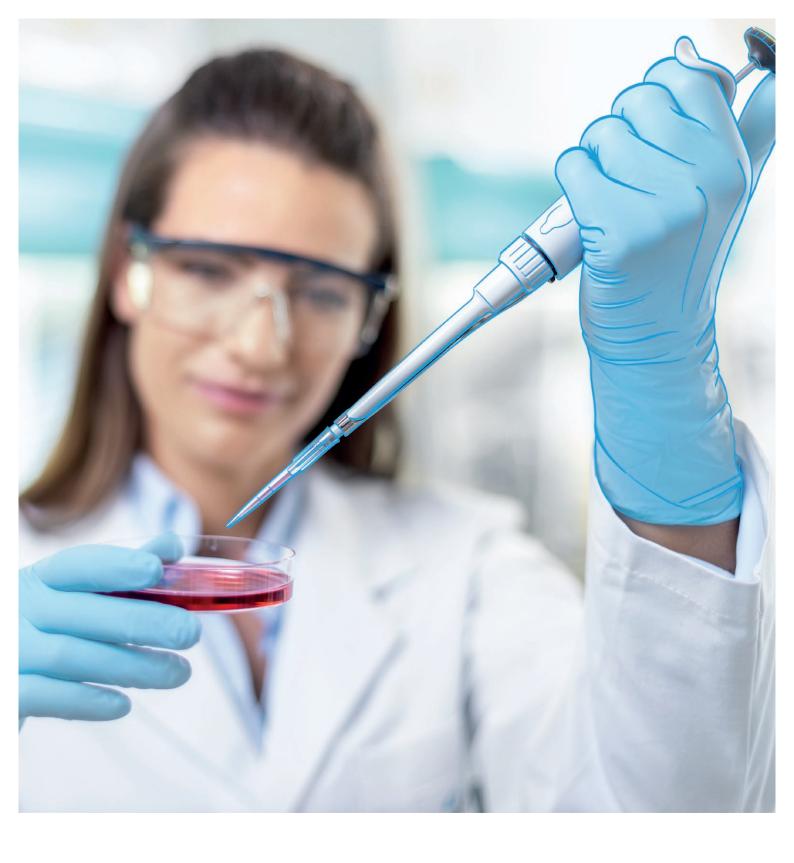

To achieve consistent, reproducible results when handling liquids in the laboratory, reliable dispensers and pipette tips are essential. With high quality materials and pipetting solutions, maximum precision can be achieved while reducing the consumption of reagents and time. Comfortable and fatigue-free working is also crucial for everyday laboratory life.

LIQUID HANDLING

/	Sapphire Pipettes	150
	Single-Channel Pipettes	15
	Multi-Channel Pipettes	15
	Pipette Carrousel	152
	Pipette Carrousel	152
/	Sapphire Pipette Tips	153
	Pipette Tips 10 μl	154
	Pipette Tips 10 μl extended	155
	Filter Tips 20 µl	156
	Filter Tips 100 µl	
	Pipette Tips 200 μl	157
	Pipette Tips 300 μl	158
	Pipette Tips 1000 μl	
	Pipette Tips 1250 µl	160
	Racks (unfilled)	16
	Macro Pipette Tip 1-5 ml	16
/	CELLSTAR® Serological Pipettes	162
	Serological Pipettes	163
	Serological Pipettes Triple-packed	164
	Serological Pipettes Special Models	
	MaxiPette	
	Pastaur / Sarum Pinattas	





/ Watch our video:



SAPPHIRE PIPETTES

Sapphire pipettes are a variable volume range of air-displacement pipettes designed and built to deliver optimal performance, ergonomics, and robustness in day-to-day use.

The lightweight design coupled with a state-of-the-art mechanism reduces pipetting forces and makes our pipettes both comfortable and easy-to-use. The contoured body shape allows for a relaxed grip whatever size your hand, and the 3-position tip ejector button makes

pipetting comfortable for both left-handed and right-handed users. The colour-coded push button allows for easier identification of the relevant tips. Sapphire pipettes provide an excellent performance within ISO 8655 which is considered the basis for determining the accuracy and precision of air-displacement pipettes. In combination with Sapphire pipette tips they show one of the lowest error tolerances on the market. All Sapphire pipettes are fully autoclavable.

- / Outstanding accuracy and precision
- / Ergonomic design and minimal pipetting forces for a new level of comfort
- Lightweight, robust and fully autoclavable
- / Digital volume setting
- / Colour-coded for tip identification



Single-Channel Pipettes

- / Light and comfortable design
- / Low pipetting forces for ease of use
- / Digital volume setting
- / Colour-coded for tip identification
- / 5 ml & 10 ml pipettes without ejector lever
- / Fully autoclavable

Description: single-channel pipette

Item no.	Colour code	Volume range	Qty. inner / outer
89000002	orange	0.2 - 2 µl	1/1
89000010	● red	1 - 10 μΙ	1/1
89000020	light yellow	2 - 20 µl	1/1
89000100	pink	10 - 100 μΙ	1/1
89000200	yellow	20 - 200 µl	1/1
89001000	blue	100 - 1,000 µl	1/1
89000500	● violet	500 - 5,000 μl	1/1
89010000	○ light blue	1,000 - 10,000 μΙ	1/1

[/] Spare parts and other accessories available on request



Multi-Channel Pipettes

- / Light and comfortable design
- / Low pipetting forces for ease of use
- / Digital volume setting
- / Colour-coded for tip identification
- / Optimised for use with Greiner Bio-One tips
- / Fully autoclavable

Item no.	Description	Colour code	Volume range	Qty. inner / outer
89000810	8-channel pipette	● red	0.5 - 10 μΙ	1/1
89000820	8-channel pipette	olight yellow	2 - 20 µl	1/1
89008200	8-channel pipette	yellow	20 - 200 µl	1/1
89008300	8-channel pipette	● green	20 - 300 µl	1/1

Item no.	Description	Colour code	Volume range	Oty. inner / outer
89001210	12-channel pipette	● red	0.5 - 10 µl	1/1
89001220	12-channel pipette	olight yellow	2 - 20 µI	1/1
89012200	12-channel pipette	yellow	20 - 200 µl	1/1
89012300	12-channel pipette	● green	20 - 300 µl	1/1

[/] Spare parts and other accessories available on request



Pipette Carrousel


- / Holds up to 7 pipettes
- / Robust and space-saving
- / Holds single and multi-channel pipettes

Item no.	Qty. inner / outer
89000099	1/1



Watch our video:



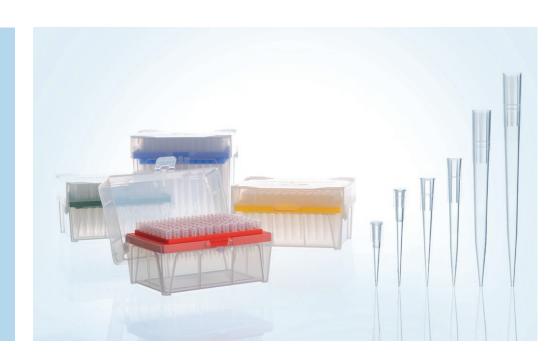

SAPPHIREPIPETTE TIPS

- Fully automated production and packaging
- / Universal fit for all common pipettes
- / Tips and racks are autoclavable
- Coloured box inserts for easy volume identification
- Low retention surface properties for high recovery rate and maximum precision

We understand that flexibility is critical in terms of modern liquid handling solutions. This is why the Sapphire tips are manufactured in eight different sizes in a volume range from 10 µl to 1250 µl, including an extended length 10 µl tip for the recovery of small sample volumes.

All our tips are made of medi-

cal grade polypropylene and are available as standard, low-retention, filter or low-retention filter tips.

Depending on the format, Sapphire pipette tips are packed in racks, bulk or refill units. Sapphire tips and pipettes when used together provide a fully optimised and harmonised liquid handling solution.



Pipette Tips

10 µl

- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification















Raw material: PP, Volume range: 0.2 - 10 µl, Nominal volume: 10 µl, Matching rack: 770310

Item no.	Feature	Product colour	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
771350	standard	○natural	● red	bag		1,000 / 10,000
771351	standard	natural	● red	rack		960 / 5,760
771352	standard	○natural	● red	rack	+	960 / 5,760
771354	standard	natural	● red	refill		960 / 4,800
771355	standard	○natural	● red	refill	+	960 / 4,800
771361	low retention	natural	● red	rack		960 / 5,760
771362	low retention	○ natural	● red	rack	+	960 / 5,760
771364	low retention	natural	● red	refill		960 / 4,800
771365	low retention	○ natural	● red	refill	+	960 / 4,800
771353	filter	natural	● red	rack	+	960 / 5,760
771363	low retention filter	natural	● red	rack	+	960 / 5,760



Pipette Tips 10 µl extended

- / Extended 10 µl tip for better recovery of small sample volumes
- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation















Raw material: PP, Volume range: 0.1 - 10 μl, Nominal volume: 10 μl, Matching rack: 770310

Item no.	Feature	Product colour	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
772350	standard	○ natural	● red	bag		1.000 / 5.000
772351	standard	natural	● red	rack		960 / 5.760
772352	standard	○ natural	● red	rack	+	960 / 5.760
772354	standard	natural	● red	refill		960 / 4.800
772355	standard	○ natural	● red	refill	+	960 / 4.800
772361	low retention	onatural on the same of the sa	● red	rack		960 / 5.760
772362	low retention	○ natural	● red	rack	+	960 / 5.760
772364	low retention	natural	● red	refill		960 / 4.800
772365	low retention	natural	●red	refill	+	960 / 4.800
772353	filter	natural	● red	rack	+	960 / 5.760
772363	low retention filter	○natural	● red	rack	+	960 / 5.760





- / Prevent contamination with liquids and aerosols during pipetting
- / filters manufactured without additives
- / Easy-to-read graduation
- / Colour-coded for tip identification

















Raw material: PP, Volume range: 2 - 20 μl, Nominal volume: 20 μl, Type of packaging: rack

Item no.	Feature	Box insert colour	Sterile	Qty. inner / outer
773353	filter	orange	+	960 / 5,760
773363	low retention filter	orange	+	960 / 5,760





- / Prevent contamination with liquids and aerosols during pipetting
- / filters manufactured without additives
- / Easy-to-read graduation
- / Colour-coded for tip identification

















Raw material: PP, Volume range: 5 - 100 μl, Nominal volume: 100 μl, Type of packaging: rack

Item no.	Feature	Box insert colour	Sterile	Qty. inner / outer
774353	filter	dark yellow	+	960 / 5,760
774363	low retention filter	dark yellow	+	960 / 5,760





- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification

















Raw material: PP, Volume range: 5 - 200 μl, Nominal volume: 200 μl, Matching rack: 770320

Item no.	Feature	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
775350	standard	yellow	bag		1,000 / 15,000
775351	standard	yellow	rack		960 / 5,760
775352	standard	yellow	rack	+	960 / 5,760
775354	standard	yellow	refill		960 / 4,800
775355	standard	yellow	refill	+	960 / 4,800
775361	low retention	yellow	rack		960 / 5,760
775362	low retention	yellow	rack	+	960 / 5,760
775364	low retention	yellow	refill		960 / 4,800
775365	low retention	yellow	refill	+	960 / 4,800
775353	filter	yellow	rack	+	960 / 5,760
775363	low retention filter	vellow	rack	+	960 / 5,760



Pipette Tips 300 µl

- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification















Raw material: PP, Volume range: 10 - 300 μl, Nominal volume: 300 μl, Matching rack: 770330

Item no.	Feature	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
776350	standard	● green	bag		1,000 / 10,000
776351	standard	● green	rack		960 / 5,760
776352	standard	● green	rack	+	960 / 5,760
776354	standard	● green	refill		960 / 4,800
776355	standard	● green	refill	+	960 / 4,800
776361	low retention	● green	rack		960 / 5,760
776362	low retention	● green	rack	+	960 / 5,760
776364	low retention	● green	refill		960 / 4,800
776365	low retention	● green	refill	+	960 / 4,800
776353	filter	● green	rack	+	960 / 5,760
776363	low retention filter	• green	rack	+	960 / 5,760



Pipette Tips 1000 µl

- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification















Raw material: PP, Volume range: 50 - 1.000 μl, Nominal volume: 1,000 μl, Matching rack: 770340

Item no.	Feature	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
777350	standard	blue	bag		1,000 / 5,000
777351	standard	blue	rack		960 / 3,840
777352	standard	blue	rack	+	960 / 3,840
777354	standard	blue	refill		960 / 4,800
777355	standard	blue	refill	+	960 / 4,800
777361	low retention	blue	rack		960/3,840
777362	low retention	blue	rack	+	960/3,840
777364	low retention	blue	refill		960 / 4,800
777365	low retention	blue	refill	+	960 / 4,800



Pipette Tips 1250 µI

- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification















Raw material: PP, Volume range: 50 - 1.250 μl, Nominal volume: 1,250 μl, Matching rack: 770340

Item no.	Feature	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
778350	standard	blue	bag		1,000 / 5,000
778351	standard	blue	rack		960 / 3,840
778352	standard	blue	rack	+	960/3,840
778354	standard	blue	refill		960 / 4,800
778355	standard	blue	refill	+	960 / 4,800
778361	low retention	blue	rack		960/3,840
778362	low retention	blue	rack	+	960 / 3,840
778364	low retention	blue	refill		960 / 4,800
778365	low retention	blue	refill	+	960 / 4,800
778353	filter	blue	rack	+	960/3,840
778363	low retention filter	blue	rack	+	960 / 3,840



Racks (unfilled)

- / Highly transparent and stable
- / Can be autoclaved and reused up to $50\,\mathrm{times}$
- / Stackable and automation-friendly
- / Colour-coded rack inserts for quick volume identification



Item no.	Description	Feature	Box insert colour	Rack colour	Qty. inner / outer
770310	suitable for 10 μl / 10 μl ext.		● red	O natural	10 / 60
770320	suitable for 200 μl		yellow	o natural	10 / 60
770330	suitable for 300 μl		● green	O natural	10 / 60
770340	suitable for 1000 μl / 1250 μl		blue	o natural	10 / 40
770370	suitable for 10 - 300 μl	without rack insert		O natural	10 / 60
770380	suitable for 1000 - 1250 μl	without rack insert		o natural	10 / 40



Macro Pipette Tip

- / Volume range from 1 5 ml
- / Suitable for Gilson® P5000

Feature: Macro, Graduation: -, Volume range: 1 - 5 ml

Item no.	Product colour	Oty. inner / outer
745290	○ natural	250 / 2,500



/ Serological pipettes are also available triple-packed.



CELLSTAR® SEROLOGICAL PIPETTES

- / Tip design guarantees a drop-free pipetting
- / Maximum accuracy
- / High optical clarity
- / Clear, easily legible graduation
- / A coloured vertical Schellbach stripe on 1, 2, 5, 10 and 25 ml pipettes makes it considerably easier to read off the volume
- / Colour code according to international standards

Greiner Bio-One offers a wide range of different serological pipettes and transfer pipettes.

The volume capacity of the serological pipettes is increased by a negative graduation. The pipettes are provided with colour codes according to international standards. The use of highgrade polystyrene guarantees maximum clarity. Expiry date and LOT number are printed on each single packaging unit. All pipettes are supplied with a filter for protection against the suction of liquid into the pipetting device.

CELLSTAR® serological pipettes are available in three different packaging options: in plastic bulk packs as well as individually wrapped in plastic or paper/plastic, both with peel-off function and additional breakthrough function.



Serological Pipettes

- / Sterile
- / Increased volume range with negative graduations
- / Lot number and expiry date on each bag
- / Filter for protection against suction of liquids into the pipetting device
- / Various packaging options





FREE OF detectable human DNA

FREE OF detectable RNase





Raw material: PS

Item no.	Description	Graduation	Nominal volume	Packaging	Type of packaging	Sterile	Qty. inner / outer
604107	1 ml pipette	1/100	1 ml		bulk	+	25 / 1,000
604160	1 ml pipette	1/100	1 ml	plastic / plastic	single-packed	+	100 / 1,000
604181	1 ml pipette	1/100	1 ml	paper/plastic	single-packed	+	100 / 1,000
710107	2 ml pipette	1/100	2 ml		bulk	+	25 / 1,000
710160	2 ml pipette	1/100	2 ml	plastic / plastic	single-packed	+	100 / 1,000
710180	2 ml pipette	1/100	2 ml	paper/plastic	single-packed	+	100 / 1,000
606107	5 ml pipette	1/10	5 ml		bulk	+	25 / 500
606160	5 ml pipette	1/10	5 ml	plastic / plastic	single-packed	+	50 / 200
606180	5 ml pipette	1/10	5 ml	paper/plastic	single-packed	+	50 / 200
607107	10 ml pipette	1/10	10 ml		bulk	+	25 / 500
607160	10 ml pipette	1/10	10 ml	plastic / plastic	single-packed	+	50 / 200
607180	10 ml pipette	1/10	10 ml	paper/plastic	single-packed	+	50 / 200
760107	25 ml pipette	2/10	25 ml		bulk	+	25 / 200
760160	25 ml pipette	2/10	25 ml	plastic / plastic	single-packed	+	50 / 200
760180	25 ml pipette	2/10	25 ml	paper/plastic	single-packed	+	50 / 200
768160	50 ml pipette	1/2	50 ml	plastic / plastic	single-packed	+	20 / 100
768180	50 ml pipette	1/2	50 ml	paper / plastic	single-packed	+	20 / 100



Serological Pipettes Triple-packed

- / Sterile
- / Increased volume range with negative graduations
- / Lot number and expiry date on each bag
- / Filter for protection against suction of liquids into the pipetting device
- / Triple-packed


STERILE DNase detectable human DNA RNase cytotoxic pyrogeni


Raw material: PS, Packaging: plastic / plastic, Triple-packed: yes

Item no.	Description	Graduation	Nominal volume	Sterile	Oty. inner / outer
604160-TRI	1 ml pipette	1/100	1 ml	+	10 / 100
710160-TRI	2 ml pipette	1/100	2 ml	+	10 / 100
606160-TRI	5 ml pipette	1/10	5 ml	+	10 / 100
607160-TRI	10 ml pipette	1/10	10 ml	+	10 / 100
760160-TRI	25 ml pipette	2/10	25 ml	+	10 / 100
768160-TRI	50 ml pipette	1/2	50 ml	+	10 / 100

Serological Pipettes

Special Models

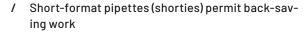

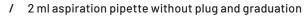

/ Single-packed







FREE OF detectable human DNA







Raw material: PS

Item no.	Description	Feature	Graduation	Nominal volume	Packaging	Sterile	Qty. inner / outer
710183	2 ml aspiration pipette	no plug	_	2 ml	paper/plastic	+	100 / 1,000
606190	5 ml pipette	shorty	1/10	5 ml	plastic / plastic	+	50 / 200
607190	10 ml pipette	shorty	2/10	10 ml	plastic / plastic	+	50 / 200



MaxiPette

- / Ergonomic design for easy pipetting
- / Variable speed control
- / Cordless and rechargeable
- / Up to eight hours of continuous operation
- / Compatible with Greiner Bio-One serological pipettes

Description: MaxiPette, Content kit: pipette controller charger benchtop stand extra filter replaceable lithium battery

Item no.	Qty. inner / outer
847070	1/1

/ Device with US or UK plug available on request




Pasteur / Serum Pipettes

- / Ideally suited for the rapid transfer of liquids
- / Sterile / non-sterile

Length: 153 mm

Item no.	Description	Graduation	Working volume	Sterile	Qty. inner / outer
700370	pasteur pipette	-	≤0,1 ml		500 / 1.500
700361	pasteur pipette	-	≤0,1 ml	+	25 / 1.000
612301	serum pipette	yes	≤2,5 ml		500 / 1.500
612361	serum pipette	yes	≤2,5 ml	+	1/800
612362	serum pipette	yes	≤2,5 ml	+	25 / 1.000

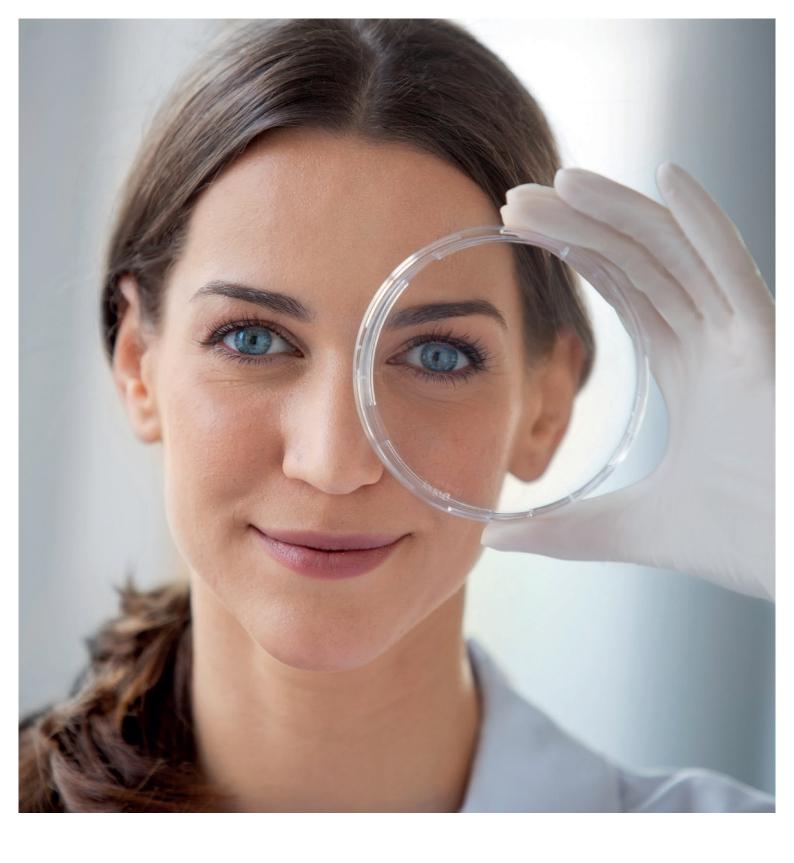

The study of microorganisms, their metabolism and their effects on other (macro-) organisms generally refers to the field of microbiology. The various detection methods are used in many research areas and in industry. For example, tests for the detection of microorganisms and for quality assurance are also applied in the production of pharmaceuticals, cosmetics, food and beverages.

MICROBIOLOGY

/	Dishes / Plates / Others168
	Petri Dishes
	Contact Dishes Divided Petri Dish
	Macroplate / Square Petri Dish
	/ OneWell Plate170
	Swab Tube Cotton Swab
	Disposable Inoculation Loops / Needles 171





/ For exact dimensions of our petri dishes, please refer to the product data sheets on our website.

DISHES / PLATES / OTHERS

Greiner Bio-One offers petri dishes with diameters of 35, 60, 94, 100 and 145 mm. Manufactured from highly transparent polystyrene (PS), they are heat-resistant up to 60 °C for use with hot agar. Depending on the application vented and non-vented design can be selected. Besides the standard round design, a square-profile dish (120 mm x 120 mm x 17 mm) and the OneWell Plate in SBS format are available.

Furthermore, Greiner Bio-One offers contact dishes which are used with hygiene monitoring including detection of contaminations as well as testing the effectiveness of cleaning and disinfection on plane surfaces by means of contact cultures. The product range is completed by a special model with two compartments which enables parallel experiments with different samples or parameters in one dish.



Petri Dishes

- / Available in different sizes
- / Easy stacking
- / With or without vents
- / Manufactured from crystal clear polystyrene

Item no.	Height	Ø nominal size	Vent nock	Qty. inner / outer
627102	10 mm	35 mm	yes	10 / 740
628102	15 mm	60 mm	yes	20 / 600
632180	16 mm	94 mm	no	20 / 480
633180	16 mm	94 mm	yes	20 / 480
664102	20 mm	100 mm	yes	15 / 360
639102	20 mm	145 mm	yes	15 / 120



Contact Dishes Divided Petri Dish

- / With or without vents
- / Manufactured from crystal clear polystyrene
- / Contact dish (sterile) with graduation to enable quick and easy analyses

Height: 15 mm

Item no.	Feature	Compartments	Ø nominal size	Total volume (well)	Vent nock	Sterile	Qty. inner / outer
629102			65 mm		no		20 / 600
629161	graduated		65 mm		no	+	20 / 600
629180	graduated		65 mm		yes	+	20 / 600
635102	2 compartments	2	94 mm	20 ml	yes		20 / 480



Macroplate / Square Petri Dish / OneWell Plate

- / With vents
- / Manufactured from crystal clear polystyrene
- / Optimised space requirement compared to round dishes


Item no.	Feature	Well format	Length	Width	Product colour	Total volume (well)	Lid	Vent nock	Oty. inner / outer
657102	microplate format	6	127.8 mm	85.5 mm	○ clear	16 mI	conden- sation rings		2 / 100
670102	microplate format	1	127.8 mm	85.5 mm	O clear	113.7 ml	yes	yes	8 / 32
688102	square		120 mm	120 mm	O clear		yes	yes	10 / 240



Swab Tube Cotton Swab

- / Suitable for taking non-human bacteriological, serological or cytological samples in veterinary research
- / For hygienic controls in food industry as well as in environmental sampling

Item no.	Description	Feature	Height	Ø	Sterile	Oty. inner / outer
420161	Swab Tube	swab tube, polystyrol	110 mm	16 mm	+	1,300 / 1,300
420180	Swab Tube	swab tube, polystyrol	110 mm	16 mm	+	1 / 1,000
421161	Swab Tube	swab tube, polystyrol	152 mm	16 mm	+	-/1,000
421180	Swab Tube	swab tube, polystyrol	152 mm	16 mm	+	1 / 700
421084	Cotton Swab		145 mm		+	1 / 1,400

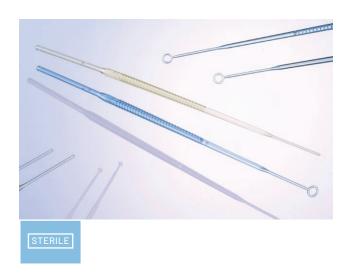


Disposable Inoculation Loops / Needles

- / Needles ideal for picking single colonies
- / Flexible loops for easier collection and inoculation
- / Colour coding for volume differentiation

Length: 200 mm

Item no.	Description	Product colour	Volume	Sterile	Qty. inner / outer
731101	inoculation loop	○white	1μΙ	+	50/2,000
731161	inoculation loop	○white	1 μΙ	+	1/600
731165	inoculation loop	○white	1 μΙ	+	10 / 3,000
731170	inoculation loop	blue	10 μΙ	+	50/2,000
731171	inoculation loop	blue	10 μΙ	+	1/600
731175	inoculation loop	blue	10 μΙ	+	10 / 3,000
731180	inoculation needle	yellow		+	50/2,000
731181	inoculation needle	yellow		+	1/600
731185	inoculation needle	yellow		+	1/3,000

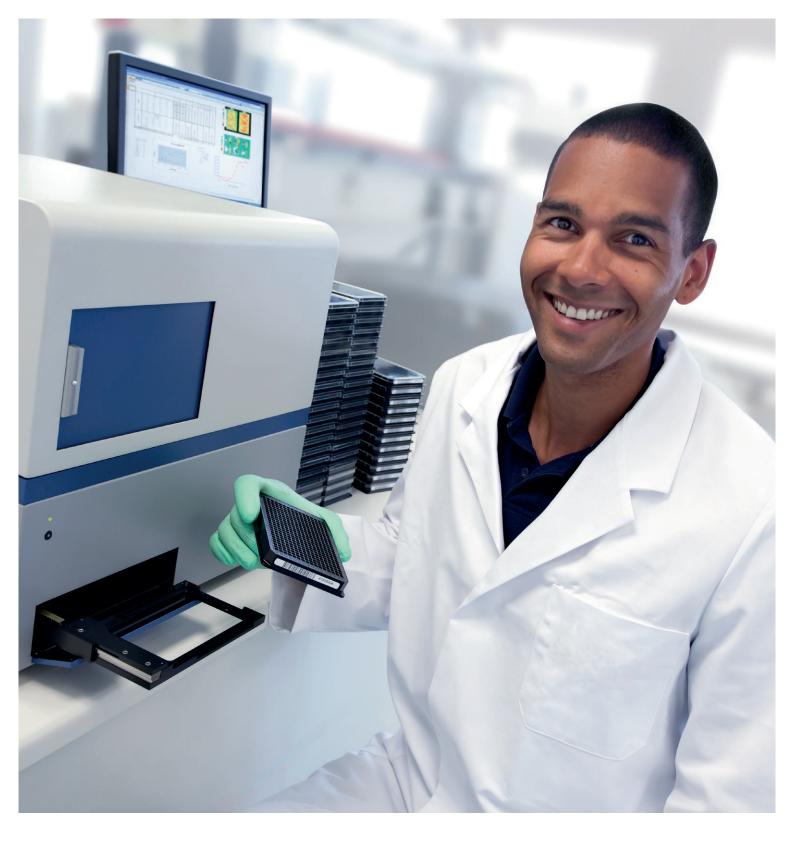

Continued progress in research and related technologies, such as microscopy, imaging, detection and liquid handling, has given rise to a wide variety of platforms used in basic science, biotechnology and pharmaceutical drug development. Today, researchers can select application-specific microplates among a broad range of products that differ in format, design, base material, colour and surface properties.


HTS-MICROPLATES

/	96 Well Microplates	. 178
	96 Well Microplates Polystyrene	179
	96 Well Microplates Polystyrene - Half Area.	180
	96 Well Microplates Polypropylene	180
/	384 Well Microplates	. 182
	384 Well Microplates Polystyrene	183
	384 Well Microplates Polystyrene	
	- Small Volume	184
	384 Well Microplates Polypropylene	184
/	1536 Well Microplates	. 185
	1536 Well Microplates	186
/	Polypropylene Storage Plates	. 187
	96 Well MASTERBLOCK® Polypropylene	188
	96 Well Storage Box	188
	384 Deep Well MASTERBLOCK®	
	Polypropylene	189
	1536 Deep Well Microplates Polypropylene	189
/	Microplates for Compound Storage	. 190
	Microplates for Compound Storage 384	
	and 1536 Well	191
/	Non-binding Microplates	. 192
	Non-binding Microplates 96 Well	193
	Non-binding Microplates 384 and 1536 Well .	193
/	Streptavidin-coated Microplates	. 195
	Streptavidin-coated 96 and 384 Well	196

/	UV-Star® Microplates	197
	UV-Star® Microplates	198
/	Lids / Sealers / CapMats	199
	Lids	200
	Sealers	200
	CapMats	20
/	Protein Crystallisation Plates .	202
	96 Well CrystalQuick	
	/ CrystalQuick Plus Plates	20
	24 Well ComboPlate CrystalBridge	
	/ Coverslips	203
	Tarasaki Platas 60 and 72 Wall	20/

MICROPLATES

- / Comprehensive microplate portfolio
- / 96 / 384 / 1536 Well
- / High-quality materials for each application
- / Different well geometries
- / Various surface modifications

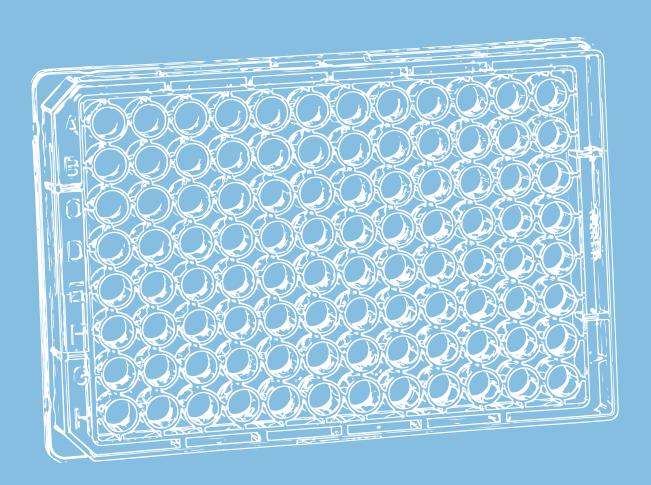

HIGH-THROUGHPUT SCREENING MICROPLATES FROM GREINER BIO-ONE


Polypropylene (PP) and polystyrene (PS) are the standard materials used to manufacture the majority of microplates. Polystyrene is a highly clear polymer with excellent optical properties which makes it ideal for precise optical measurements. Polystyrene is also characterised by its ability to bind biomolecules, such as proteins, and it is therefore often used for manufacturing immunological products. Polystyrene is suitable for work with cell cultures.

Polypropylene is characterised by its excellent chemical

and thermal stability. It is the ideal polymer for storage vessels or microplates. In addition, Greiner Bio-One manufactures microplates with special requirement profiles, such as the UV-Star® or SCREENSTAR microplates made from different cycloolefins. They are characterised by their low level of autofluorescence, exceptionally high clarity, especially in the UV range, and greater chemical stability when compared with polystyrene.

GREINER BIO-ONE MICROPLATES:

- / Manufactured under DIN ISO 9001 quidelines
- / Full traceability
- Footprint compatible with automated systems
- / Free of detectable endotoxins
- / Free of detectable DNase, RNase and human DNA
- Available with barcode on request

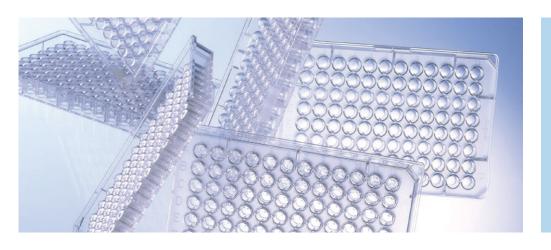

OUR PORTFOLIO
OFFERS THE
RIGHT PLATFORM
AND COATING
FOR EVERY
APPLICATION

μClear® and UV-Star® Film Bottom Microplates

The development of a new and patented processing technique has made it possible for us to produce microplates with ultra-thin films, without the use of adhesives or solvents – the μ Clear® and UV-Star® products. This special method eradicates the risk of leaking wells.

SensoPlate Glass Bottom Microplates

SensoPlate glass bottom microplates consist of a black pigmented polystyrene frame on to which a 175 µm thick borosilicate glass bottom is bonded. Thanks to the high optical quality of the glass bottom as well as the minimal bending, SensoPlate are

especially recommended for fluorescence correlation spectroscopy and sophisticated microscopic applications.

Black or White?

White microplates are usually used for luminescence measurements (e. g. Luciferase Reporter Assays) and black microplates for fluorescence. measurements (e.g. Green Fluorescence Protein). The critical properties in these methods, such as background, autofluorescence or crosstalk are considerably improved by the use of blackor white pigmented microplates.





Further information on microplates:

Forum No. 19: Base Material and Surface Modifications of Greiner Bio-One Multiwell Plates and Microplates - An Overview (F073793)

Microplate Selection Guide (F073048)

Microplate Dimensions Guide (F073027).

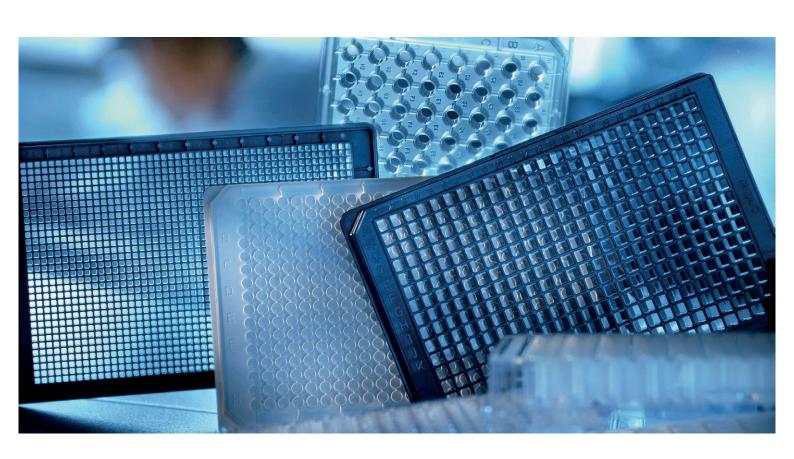

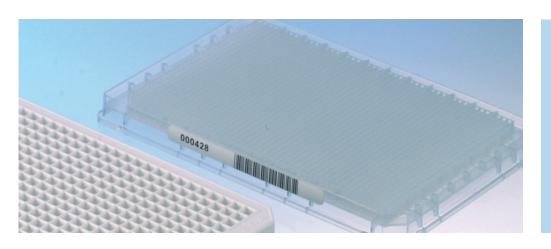

BARCODE:

You can download the microplate order form directly from our website www.gbo.com (Item no. F073010)

MICROLON, FLUOTRAC, LUMITRAC

These brands stand for the quality of our immunology products. MICROLON are clear microplates for transmission measurements. FLUOTRAC are black microplates for fluorescence measurements. LUMITRAC are white microplates for luminescence measurements.

Non-binding Surfaces

Non-binding surfaces are characterised by their low binding capacity for biomolecules such as DNA, RNA, peptides and proteins. The repellent property of the non-binding surfaces for biomoleculescan be advantageous in biochemical assays by increasing the sensitivity, reducing the background and improving the signal-to-noise ratio.

MICROLON, FLUOTRAC, LUMITRAC

Quality brands for immunological Plates:

MICROLON

clear microplates for transmission measurements

FLUOTRAC

black microplates for fluorescence measurements

LUMITRAC

white microplates for luminescence measurements


Non-binding surfaces

Low binding capacity for biomolecules such as DNA, RNA, peptides and proteins for sensitive biochemical assays

Cell Culture Surface

Significantly improves the adhesion of cells and the binding of proteins to the plastic surface

(see chapter cell culture)

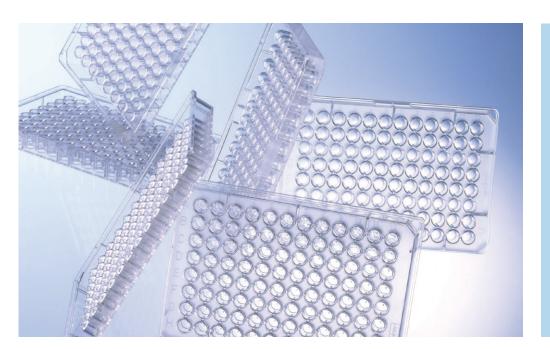



/ For details regarding the different well profiles, please refer to the technical appendix.

96 WELL MICROPLATES

Since its introduction in the 1960's applications for the 96 well microplate have continually increased to the extent that it is impossible to envisage modern research and industry without it today. Greiner Bio-One has been manufacturing microplates and strip microplates for diagnostics

and immunological research for over 40 years. A large number of different 96 well microplates is available in a wide variety of surface treatments. The spectrum ranges from clear bottom microplates and completely black or white microplates to UV-Star® products.

- Available in polystyrene, polypropylene and cycloolefin
- / Clear / black / white
- / With U-bottom, V-bottom or F-bottom
- / Sterile / non-sterile
- / Non-treated or in high-binding quality
- / In non-binding quality



96 Well Microplates Polystyrene

- / With U-bottom, V-bottom or F-bottom
- / Sterile / non-sterile
- / Non-treated or in high binding quality
- / With solid bottom or µClear® film bottom

Well format: 96, Raw material: PS, Lid: no

Item no.	Well profile	Bottom	Binding character- isitc	Binding brand name	Surface treatment	Product colour	Working volume (well)	Sterile	Oty. inner/ outer
650101	U-bottom	solid			untreated	○clear	40 μl - 280 μl		10 / 100
650161	U-bottom	solid			untreated	○ clear	40 μl - 280 μl	+	2 / 100
651101	V-bottom	solid			untreated	○clear	40 μl - 200 μl		10 / 100
651161	V-bottom	solid			untreated	○ clear	40 μl - 200 μl	+	2 / 100
655101	F-bottom	solid			untreated	○ clear	25 μl - 340 μl		10 / 100
655161	F-bottom	solid			untreated	○ clear	25 μl - 340 μl	+	2 / 100
655075	F-bottom / Chimney Well	solid			untreated	○white	25 μl - 340 μl		10 / 40
655074	F-bottom / Chimney Well	solid	high-bind- ing	LUMITRAC 600		○white	25 μl - 340 μl	+	10 / 40
655077	F-bottom / Chimney Well	solid	high-bind- ing	FLUOTRAC 600		● black	25 μl - 340 μl	+	10 / 40
655076	F-bottom / Chimney Well	solid			untreated	● black	25 μl - 340 μl		10 / 40
655095	F-bottom / Chimney Well	µClear®			untreated	○white	25 μl - 340 μl		10 / 40
655094	F-bottom / Chimney Well	μClear®	high-bind- ing			○white	25 μl - 340 μl	+	10 / 40
655097	F-bottom / Chimney Well	μClear®	high-bind- ing			● black	25 μl - 340 μl	+	10 / 40
655096	F-bottom / Chimney Well	µClear®			untreated	● black	25 µl - 340 µl		10 / 40



96 Well Microplates Polystyrene - Half Area

- / Sterile / non-sterile
- / Reduction of sample volume by up to 50 %
- / Standardised pathlength (1 cm=170 µl, 0.5 cm=80 µl)
- / Non-treated or in high-binding quality
- / With solid bottom or µClear® film bottom

Well format: 96, Well profile: F-bottom, Raw material: PS, Plate design: half area, Working volume (well): 15 μl - 175 μl, Lid: no

Item no.	Bottom	Binding characterisitc	Surface treatment	Product colour	Sterile	Qty. inner / outer
675161	solid		untreated	○ clear	+	10 / 40
675101	solid		untreated	○clear		10 / 40
675074	solid	high-binding		○white	+	10 / 40
675075	solid		untreated	○white		10 / 40
675077	solid	high-binding		● black	+	10 / 40
675076	solid		untreated	● black		10 / 40
675096	μClear®		untreated	● black		10 / 40

[/] Further information on Half Area Microplates: Forum No. 16: 96 Well Half Area Microplates and their Application in Fluorescence, Luminescence and Transmission Measurements (F073121)




96 Well Microplates Polypropylene

- / Ideally suited for the storage of active agents, DNA/RNA or stock cultures
- / High chemical resistance and temperature tolerance
- / Black plates for fluorescence polarisation

Well format: 96, Bottom: solid, Raw material: PP, Surface treatment: untreated, Lid: no

Item no.	Well profile	Product colour	Working volume (well)	Sterile	Qty. inner / outer
650201	U-bottom / Chimney Well	○ natural	50 μΙ - 300 μΙ		10 / 100

Item no.	Well profile Product cold		Working volume (well)	Sterile	Oty. inner / outer
650261	U-bottom / Chimney Well	onatural	50 μΙ - 300 μΙ	+	10 / 100
650209	U-bottom / Chimney Well	● black	50 μΙ - 300 μΙ		10 / 100
655201	F-bottom / Chimney Well	onatural	25 μΙ - 370 μΙ		10 / 100
655209	F-bottom / Chimney Well	● black	25 μΙ - 370 μΙ		10 / 100
651201	V-bottom/Chimney Well	onatural	50 μΙ - 335 μΙ		10 / 100
651209	V-bottom/Chimney Well	● black	50 μΙ - 335 μΙ		10 / 100

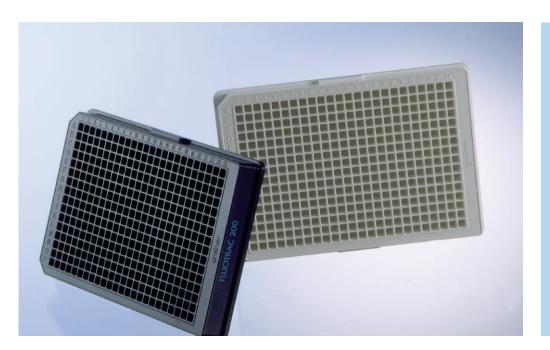



/ For cell culture treated microplates, please refer to chapter cell culture

384 WELL MICROPLATES

Drug screening has undergone rapid development over the past years.

The number of tests with new targets and the number of active agents to be tested is constantly increasing. Volume reduction, simple testing and cost savings are some of the highest priorities and high-format microplates with a low well volume are one of the most important tools in achieving this. One of the

first higher-format microplates was the 384 well plate, launched by Greiner Bio-One in 1994/1995. Compared with the 96 well microplate, the number of wells is quadrupled in this microplate, combined with a volume reduction from 382 μ l to 131 μ l. The well-to-well spacing is 4.5 mm (96 well plate: 9 mm). The external dimensions of the 384 well microplates are compatible with automated systems.

- Available in polystyrene, polypropylene and cycloolefin
- / Clear / black / white
- / With F-bottom, V-bottom or as Small Volume option
- / Solid bottom or µClear° film bottom
- / Non-treated or in high-binding quality
- / In non-binding quality




384 Well Microplates Polystyrene

- / Solid bottom or µClear® film bottom
- / Sterile / non-sterile
- / Non-treated or in high-binding quality

Well format: 384, Well profile: F-bottom, Raw material: PS, Working volume (well): $15\,\mu$ l - $110\,\mu$ l

Item no.	Bottom	Binding characterisitc	Binding brand name	Surface treatment			Sterile	Qty. inner / outer
781101	solid			untreated	○ clear	no		10 / 100
781061	solid	high-binding			○ clear	no	+	10 / 40
781162	solid			untreated	○ clear	no	+	10 / 100
781185	solid			untreated	○ clear	yes	+	1/32
781186	solid			untreated	○ clear	yes	+	8 / 32
781074	solid	high-binding	LUMITRAC 600		○ white	no	+	10 / 40
781075	solid			untreated	○white	no		10 / 40
781077	solid	high-binding	FLUOTRAC 600		● black	no	+	10 / 40
781076	solid			untreated	● black	no		10 / 40
781095	µClear®			untreated	○ white	no		10 / 40
781097	µClear®	high-binding			● black	no	+	10 / 40
781096	μClear®			untreated	black	no		10 / 40

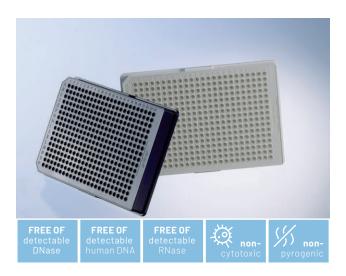

384 Well Microplates

Polystyrene - Small Volume

HiBase

- Perfect for top reading even at low working volumes
- / Savings in reagent similar to 1536 well microplates
- / Made of clear / black / white polystyrene for transmission, fluorescence or luminescence measurements

Well format: 384, Well profile: F-bottom, Bottom: solid, Raw material: PS, Surface treatment: untreated, Plate geometry: HiBase, Plate design: Small Volume, Working volume (well): 4 μl - 25 μl, Lid: no

Item no.	Product colour	Qty. inner / outer
784101	○ clear	10 / 40
784075	○ white	10 / 40
784075-25	○white	25 / 150
784076	● black	10 / 40
784076-25	● black	25 / 150



384 Well Microplates

Polypropylene


- / Ideally suited for the storage of active agents, DNA/RNA or stock cultures
- / High chemical resistance and temperature tolerance
- / Available natural or black
- / Item No. 781201-906 for Acoustic Liquid Handling

Well format: 384, Bottom: solid, Raw material: PP, Surface treatment: untreated, Lid: no

Item no.	Well profile	Product colour	Plate geometry	Plate design	Working volume (well)	Qty. inner / outer
781201	F-bottom	○ natural			15 μΙ - 145 μΙ	10 / 100
784201	V-bottom	natural	Deep Well	Small Volume	1 μΙ - 90 μΙ	10 / 100
781201-906	F-bottom	○ natural			15 µl - 145 µl	10 / 100
781209	F-bottom	black			15 μΙ - 145 μΙ	10 / 100
781280	V-bottom	○ natural			13 μΙ - 120 μΙ	10 / 100



/ For cell culture treated microplates, please refer to chapter cell culture

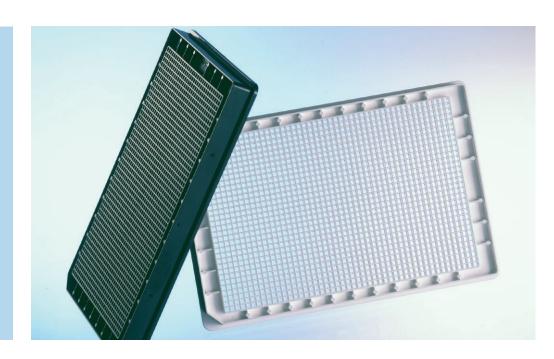


1536 WELL MICROPLATES

- / Available in polystyrene, polypropylene and cycloolefin
- / Clear / black / white
- HiBase for top-reading applications
- / Solid bottom or µClear° film bottom
- / Sterile / non-sterile
- / Non-treated or in high-binding quality

The highest possible degree of automation, optimal performance and cost savings continue to be the requirements placed on microplates for high-throughput screening.

In 1997, shortly after the launch of the 384 well microplates, Greiner Bio-One was the first manufacturer to introduce another innovative microplate format - the 1536 well microplate. The external dimensions were the same as those used in the 96 well and 384 well microplates. However, to utilise the available space most efficiently, the number of wells was increased fourfold from 384 to 1536.

Close cooperation with numerous users has now led to the development of a broad product range, and the constant drive towards improvements in quality has, for example, led to a reduction in curvature of the plates to $<100\ \mu m.$

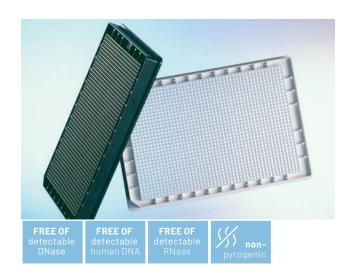

1536 Well Microplates


- / Solid bottom or $\mu Clear^{\circ}$ film bottom
- / Sterile / non-sterile
- / Non-treated or in high-binding quality

Well format: 1536, Well profile: F-bottom, Raw material: PS, Plate geometry: HiBase, Working volume (well): 3 μl - 10 μl, Lid: no

Item no.	Bottom	Binding characterisitc	Binding brand name	Surface treatment	Product colour	Sterile	Qty. inner / outer
782101	solid			untreated	○ clear		15 / 60
782061	solid	high-binding	MICROLON 600		○ clear	+	15 / 60
782075	solid			untreated	○white		15 / 60
782074	solid	high-binding	LUMITRAC 600		○white	+	15 / 60
782076	solid			untreated	● black		15 / 60
782077	solid	high-binding	FLUOTRAC 600		● black	+	15 / 60
782095	µClear®			untreated	○white		15 / 60
782097	µClear®	high-binding			● black	+	15 / 60
782096	µClear®			untreated	● black		15 / 60



/ Compound storage plates for acoustic liquid handling can also be found in this chapter.

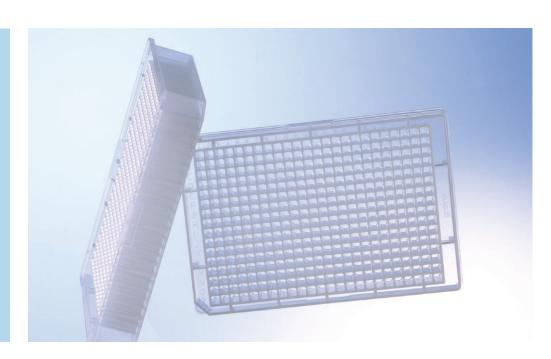

POLYPROPYLENE STORAGE PLATES

Greiner Bio-One polypropylene microplates are perfect storage plates for active agents, patient samples or biomolecules.

Their most important properties are biological inertness, resistance to numerous solvents commonly used in the laboratory, such as DMSO and high temperature resistance. The footprint

is compatible with automated systems.

The microplates are also characterised by elevated well walls which make it possible to easily seal them. The portfolio ranges from the 96 well MASTERBLOCK® with volumes of 0.5 ml, 1 ml or 2 ml over a MASTERBLOCK® with 384 wells or 1536 wells up to the 96 well storage box.

- / Ideally suited for the storage of active agents, DNA/RNA or stock cultures
- / High chemical resistance and temperature tolerance




96 Well MASTERBLOCK° Polypropylene

- / Ideal for storing non-human sample material and cultivating bacteria / yeast
- / Uniform external dimensions and tolerances
- / Alphanumeric well coding
- / Sealable with adhesive films and heat sealers
- / Sealable with CapMats

Well format: 96, Bottom: solid, Raw material: PP


Item no.	Well profile	Product colour	Total volume (well)	Lid	Sterile	Qty. inner / outer
780201	U-bottom	○ natural	1 ml	CapMat 381070, 381061		1/50
780261	U-bottom	onatural on the same of the sa	1 ml	CapMat 381070, 381061	+	1/50
780215	U-bottom	○ natural	1 ml	CapMat 381070, 381061		5/50
786201	V-bottom	natural	0.5 ml	CapMat 381070, 381061		8 / 80
786261	V-bottom	○ natural	0.5 ml	CapMat 381070, 381061	+	1/80
780270	V-bottom	o natural	2 ml	CapMat 381080, 381081		1/50
780271	V-bottom	○ natural	2 ml	CapMat 381080, 381081	+	1/50
780285	V-bottom	onatural on the state of the st	2 ml	CapMat 381080, 381081		5/50

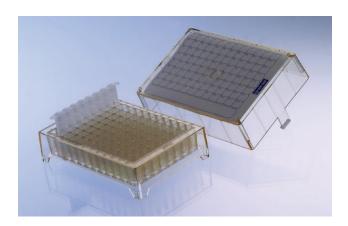

96 Well Storage Box

- / Storage box made of polycarbonate
- / With / without 96 polypropylene vessels
- / Coding card included
- / All components autoclavable

Item no.	Sterile	Qty. inner / outer
975502		1 / 120
975561	+	1/50
975570		1/50

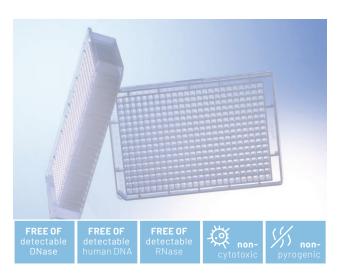


384 Deep Well MASTERBLOCK® Polypropylene

- / Ideal for compound libraries and applications with larger volumes
- / Conical well shape for precise pipetting
- / Alphanumeric well coding
- / Sealable with adhesive films and heat sealers

Well format: 384, Well profile: V-bottom, Bottom: solid, Raw material: PP, Plate geometry: Deep Well, Lid: no

Item no.	Product colour	Sterile	Qty. inner / outer
781270	○ natural		6 / 60
781271	natural	+	6 / 60

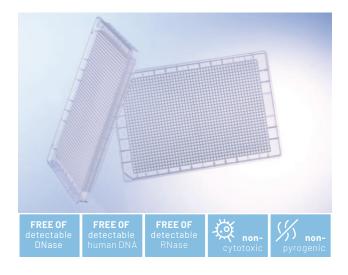


1536 Deep Well Microplates Polypropylene

- / Uniform external dimensions and tolerances
- / Total volume of 18 µl
- / Alphanumeric well coding
- / Sealable with adhesive films and heat sealers

Well format: 1536, Well profile: V-bottom, Bottom: solid, Raw material: PP, Plate geometry: Deep Well, Working volume: 3 μl - 15 μl, Lid: no

Item no.	Product colour	Sterile	Qty. inner / outer
782261	○ natural	+	15 / 60
782270	onatural on the same of the sa		15 / 60

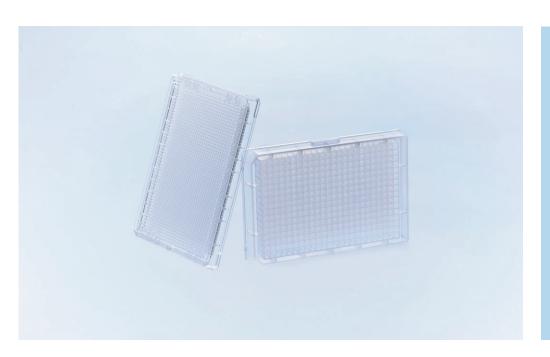



/ Watch our video:



MICROPLATES FOR COMPOUND STORAGE

Polypropylene is still the material of choice for storage plates, but the material class of cycloolefins is becoming more routinely used because of its unsurpassed performance for a wide range of applications.

In compound storage, plates made from cycloolefins offer the

best combination of chemical resistance to polar solvents, like DMSO, and optical clarity. In addition, the dimensional stability and glass-like optical properties make this material ideally suited for plates in fully automated systems.

- / Suited for compound storage
- / Resistant against polar solvents
- / Excellent water and vapour barrier function
- / Almost no leachables
- / Low biomolecule binding
- / Glass-like optical properties



Microplates for Compound Storage 384 and 1536 Well

Microplates for Acoustic Liquid Handling

- / Stringent production specifications for a constant bottom quality
- / Microplates are deionised and packed in antistatic bags

Well profile: F-bottom, Bottom: solid, Lid: no

Item no.	Well format	Raw material	Surface treatment	Product colour	Plate geometry	Plate design	Working volume (well)	Oty. inner / outer
781201-906	384	PP	untreated	○natural			15 μΙ - 145 μΙ	10 / 100
793855	384	cycloolefin		○ clear	HiBase	Small Volume	1 μΙ - 25 μΙ	15 / 60
782855	1536	cycloolefin		○clear	HiBase		1 μΙ - 10 μΙ	15 / 60
792870-906	1536	cycloolefin		○ clear			1 μΙ - 14 μΙ	15 / 60





/ For more detailed information on well profiles and technical details please refer to our data sheets on our website: www.gbo.com

NON-BINDING MICROPLATES

Polystyrene microplates with non-treated surfaces are commonly used for homogeneous biochemical HTS assays and demonstrate low and reproducible biomolecule binding.

However, even low amounts of biomolecular binding (e. g. DNA, RNA, proteins, peptides) can cause an undesirable increase in background, resulting in decreased signal-to-noise ratio.

Greiner Bio-One's non-binding microplate surfaces prevent unwanted non-specific binding, especially advantageous for sensitive biochemical assays. Characterised by low protein, DNA, RNA and peptide binding properties the non-binding surfaces significantly increase assay sensitivity by reducing background and improving signal-to-noise ratio.

- / 96 / 384 / 1536 well format
- / Clear / black / white
- / Solid bottom or µClear° film bottom
- / Ultra low non-specific biomolecular binding properties
- / Long-term surface stability without degradation or leaching



Non-binding Microplates 96 Well

- / Clear / black / white
- / U-bottom, V-bottom or F-bottom/chimney well
- / Solid bottom or $\mu Clear^{\circ}$ film bottom

Well format: 96, Raw material: PS, Binding characterisitc: Non-binding, Lid: no

Item no.	Well profile	Well profile Bottom Product colour		Working volume (well)	Oty. inner / outer	
650901	U-bottom	solid	○clear	40 µl - 280 µl	10 / 40	
651901	V-bottom	solid	○clear	40 μΙ - 200 μΙ	10 / 40	
655901	F-bottom / Chimney Well	solid	○clear	25 μΙ - 340 μΙ	10 / 40	
655904	F-bottom / Chimney Well	solid	○ white	25 μΙ - 340 μΙ	10 / 40	
655900	F-bottom / Chimney Well	solid	● black	25 μΙ - 340 μΙ	10 / 40	
655903	F-bottom / Chimney Well	µClear®	○ white	25 μΙ - 340 μΙ	10 / 40	
655906	F-bottom / Chimney Well	µClear®	black	25 μΙ - 340 μΙ	10 / 40	



Non-binding Microplates 384 and 1536 Well

- / Clear / black / white
- / With F-bottom or as Small Volume option
- / Solid bottom or µClear° film bottom

Well profile: F-bottom, Raw material: PS, Binding characterisitc: Non-binding, Lid: no


Item no.	Well format	Bottom	Product colour	Plate geometry	Plate design	Working volume (well)	Oty. inner / outer
781901	384	solid	○ clear			15 μΙ - 110 μΙ	10 / 40
781904	384	solid	○white			15 μΙ - 110 μΙ	10 / 40
781900	384	solid	● black			15 μΙ - 110 μΙ	10 / 40
781903	384	µClear®	○white			15 μΙ - 110 μΙ	10 / 40

HTS-Microplates Non-binding Microplates

Item no.	Well format	Bottom	Product colour	Plate geometry	Plate design	Working volume (well)	Qty. inner / outer
781906	384	µClear®	● black			15 μΙ - 110 μΙ	10 / 40
784904	384	solid	○ white	HiBase	Small Volume	4 μΙ - 25 μΙ	10 / 40
784900	384	solid	● black	HiBase	Small Volume	4 μΙ - 25 μΙ	10 / 40
782904	1536	solid	○ white	HiBase		3 μΙ - 10 μΙ	15 / 60
782900	1536	solid	● black	HiBase		3 μΙ - 10 μΙ	15 / 60



/ For more detailed information on well profiles and technical details please refer to our data sheets on our website: www.gbo.com

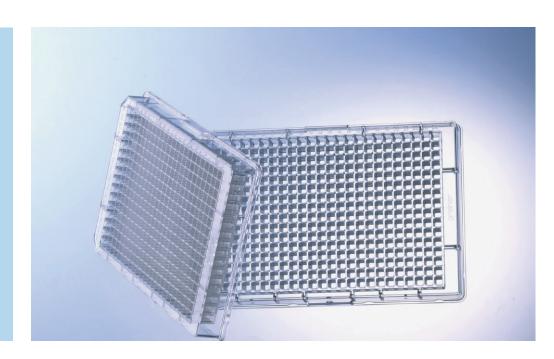

STREPTAVIDIN-COATED MICROPLATES

Streptavidin-coated solid phases serve as reliable binding surfaces for all types of biotinylated molecules.

Numerous ligands can be biotinylated simply and due to the low molecular weight of biotin (244 Da) the functionality of the molecules is normally not impaired. Thus streptavidin-coated solid phases make it possible to rapidly isolate, determine and quantify components from a reaction mixture. By immobilising the bi-

otinylated substance, it is also possible to reproduce complete reaction chains on a streptavidin solid phase, e. g. enzyme immunoassays, enzyme activity assays, DNA hybridisation techniques and quantification of PCR products.

The high-purity streptavidin is bound to the plate surface in a uniform and stable layer. The coefficient of variation from well to well is under 5 % for 96 well microplates and under 8 % for 384 well microplates.

- / Available in 96 well and 384 well format
- / Clear / black / white
- / 3-year shelf life at room temperature
- / All plates pre-blocked and ready-to-use



Streptavidin-coated Microplates 96 and 384 Well

- / Long shelf life at room temperature
- / Lot number on each pack
- / Pre-blocked with BSA

Bottom: solid, Raw material: PS, Surface treatment: Streptavidin, Lid: no

Item no.	Well format	Well profile	Product colour	Qty. inner / outer
655990	96	C-bottom	○ clear	5 / 40
655995	96	C-bottom	○white	5 / 40
655997	96	C-bottom	● black	5 / 40
781990	384	F-bottom	○clear	5 / 40
781995	384	F-bottom	○white	5 / 40
781997	384	F-bottom	● black	5 / 40

 $^{{\}it I} \quad {\sf Further \, strept avidin-coated \, microplates \, are \, available \, on \, request.}$



/ Application Note: UV/ VIS Spectroscopy in Microplates UV-Star®, µClear®, MICROLON and CELLSTAR®



UV-STAR® MICROPLATES

UV / VIS spectroscopy is a classical analytical method for determining the chemical constitution of a substance and its concentration in aqueous solution.

UV / VIS spectroscopy is usually conducted in quartz glass cuvettes. However, cuvettes do not provide sufficient throughput when dealing with large amounts of samples, and microplates can be used to speed up work.

Standard polystyrene microplates are only partially suitable for transmission measurements in the UV. Polystyrene absorbs UV especially in the short-wavelength range (< 320 nm). μ Clear $^{\circ}$ microplates with a thin polystyrene film base al-

ready have much lower background values and can be used up to 340 nm without any problem. The adaptation of the patented μClear^* process technology to a new, innovative UV-transparent material has made it possible to produce microplates that extend the transmission range up to 230 nm.

For the determination of nucleic acid and protein concentrations at 260 nm or 280 nm without background interference UV-Star® microplates are the ideal alternative to expensive and fragile quartz glass plates or cuvettes. UV-Star® plates are also DMSO-resistant and can be stored at -20 °C without any problem.

- / Available in 96 well and 384 well format
- / With cycloolefin film bottom
- / Optical window down to 230 nm ideal for nucleic acid determinations at 260 nm/280 nm
- / For measurements of protein concentration at 280 nm



UV-Star® Microplates


- / Available in 96 well and 384 well format
- / With cycloolefin film bottom
- / Optical window down to 230 nm ideal for nucleic acid determinations at 260 nm/280 nm
- / For measurements of protein concentration at 280 nm

Well profile: F-bottom, Bottom: Cycloolefin film, Raw material: COC, Surface treatment: untreated, Lid: no

Item no.	Well format	Product colour	Plate design	Working volume (well)	Qty. inner / outer
675801	96	○ clear	half area	15 μΙ - 175 μΙ	10 / 40
655801	96	○ clear		25 μΙ - 340 μΙ	10 / 40
781801	384	○clear		15 μΙ - 110 μΙ	10 / 40



/ Forum No 6: Sealers for microplates and their areas of application in molecular biology and cell culture (F073013)

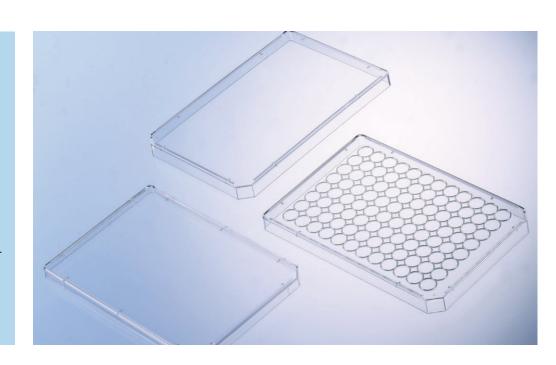

LIDS / SEALERS / CAPMATS

- / Polystyrene lids in three different profile heights with and without condensation rings
- / Classical adhesive films coated with acrylate adhesive as well as advanced adhesive films with pressure-sensitive silicone adhesive
- / EVA CapMats for sealing of 96 well plates


Lids offer protection against contamination and evaporation during sample storage and cell cultivation.

In addition, Greiner Bio-One offers a full set of adhesive **sealers** for a wide range of routine and specialised applications, e.g. high-throughput screen-

ing, immunology, microbiology, molecular biology and PCR.

An alternative method for sealing 96 well plates are CapMats.

They are made of ethyl vinyl acetate (EVA), are resistant to DMSO and can be used in a temperature range between -20 °C and +60 °C. CapMats are not pierceable.

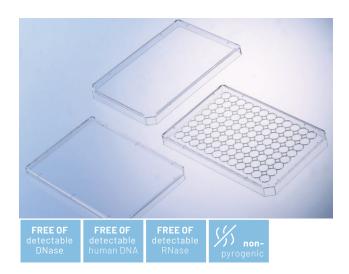

Lids

All sterile lids are non-cytotoxic.

Description: Lid, Raw material: PS

Item no.	Height	Condensation rings	Lid type	Sterile	Qty. inner / outer
656101	9 mm	no	high		1 / 100
656161	9 mm	no	high	+	1 / 100
656170	9 mm	yes	high		1 / 100
656171	9 mm	yes	high	+	1 / 100
656190	6 mm	no	flat		20 / 200
656191	6 mm	no	flat	+	20 / 200
691101	4.8 mm	no	ultra low		25 / 100
691161	4.8 mm	no	ultra low	+	5 / 100



Sealers

The classical sealers such as EASYseal, AMPLIseal, SILVERseal and BREATHseal are coated with an acrylate adhesive. The advanced sealer VIEWseal is coated with a pressure-sensitive silicone adhesive.

Item no.	Description	Feature	Pierceable	Sterile	Qty. inner / outer
676001	1 EASYseal clear				100 / 2,000
676090	SILVERseal	aluminium foil	yes		100 / 1,200
676070	VIEWseal	clear			100 / 1,200
676040 AMPLIseal		clear			100 / 2,000
676050	BREATHseal	gas-permeable			50 / 2.500

Item no.	Description	Feature	Pierceable	Sterile	Qty. inner / outer
676051	BREATHseal	gas-permeable		+	50 / 500

/ Forum No. 6: Sealers for microplates and their areas of application in molecular biology and cell culture (F073013)




CapMats

- / Made of ethyl vinyl acetate (EVA)
- / Available for 96 well microplates and MASTER- BLOCK°
- / Sterile / non-sterile

Description: 96 Well CapMat, Pierceable: no, Raw material: EVA

Item no.	Nap shape	Sterile	Qty. inner / outer
381070	round		10 / 50
381061	round	+	1/50
381080	square		10 / 50
381081	square	+	1/50

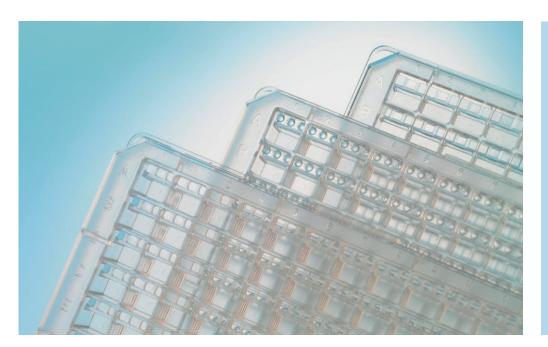



/ Forum No. 7: Advanced high-throughput platforms for protein crystallisation (F073016)

PROTEIN CRYSTALLISATION PLATES


Detailed knowledge of interactions between the structure and function of biomolecules is fundamental for biological, medical and pharmaceutical research. A reliable method is the X-ray structure analysis of protein crystals.

In order to show three-dimensional molecular structures, the proteins must first be crystallised. Since every protein is different and a variety of factors influence protein crystallisation, the optimal conditions are usually tested with high-throughput technologies (pipetting robots).

Greiner Bio-One has specially developed microplates and accessories of the CrystalStar series to meet the requirements of high-throughput crystallisation in a short time and with relatively small amounts of protein. Techniques such as vapour diffusion (sitting drop, hanging drop) and microbatch under oil can be carried out effectively.

Plates with hydrophobic surface are particularly well suited for nanolitre crystallisation of membrane proteins. LBR (low birefringence) plates are specifically designed for the use of polarised light.

- / For vapour diffusion or microbatch under oil
- / Made of polystyrene or polyolefin (LBR plates)
- / Footprint conforming to ANSI 1-2004 (except Terasaki plates)
- / Barcode labelling on request
- / CrystalQuick plates with alphanumeric well coding

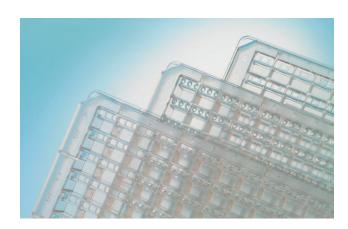

96 Well CrystalQuick / CrystalQuick Plus Plates

- / Crystallisation plates for sitting drop applications with different well profiles and material properties
- / Optimised for sealing with VIEWseal or AMPLIseal
- / Alphanumeric well coding

Item no.	Height	Material properties	Well profile	Plate design	Well / reservoir	Well volume, max	Volume per reservoir	Oty. inner / outer
609171	8 mm	standard	flat	square (LP)	1	3.9 µl	140 μΙ	20/80
609180	8 mm	hydrophobic	flat	square(LP)	1	3.9 µl	140 μΙ	20 / 80
609871	8 mm	LBR	flat	square(LP)	1	3.9 µl	140 μΙ	20 / 80
609101	14.4 mm	standard	flat	square (SW)	3	4.1 µI	320 µI	10 / 40
609120	14.4 mm	standard	concave	round (RW)	3	1.9 μΙ	320 µI	10 / 40
609130	14.4 mm	hydrophobic	flat	square(SW)	3	4.1 µI	320 µI	10 / 40
609801	14.4 mm	LBR	flat	square(SW)	3	4.1 µI	320 µI	10 / 40
609820	14.4 mm	LBR	concave	round (RW)	3	1.9 μΙ	320 µI	10 / 40
609830	14.4 mm	LBR, hydropho- bic	flat	square(SW)	3	4.1 µI	320 µI	10 / 40




24 Well ComboPlate CrystalBridge / Coverslips

- / Universal 24 well crystallisation plate with coverslips and CrystalBridge for sitting drop applications
- / Siliconised coverslips available

Item no.	Description	Feature	Well format	Ø well	Bottom shape	Well volume, max	Lid	Qty. inner / outer
662150	ComboPlate		24	16.3 mm	flat	3,300 µI	yes	6 / 24
662145	CrystalBridge		1	4.6 mm	concave	45 µI	no	-/250
501870	round coverslips 18 mm	siliconised glass 0.19 - 0,22 mm						100 / 1,000

Item no.	Description	Feature	Well format	Ø well	Bottom shape	Well volume, max	Lid	Qty. inner / outer
503870	round coverslips 22 mm	siliconised glass 0.19 - 0,22 mm						100 / 1.000
503850	round coverslips 22 mm	siliconised glass 0.5 - 0,6 mm						100 / 1,000

 $\emph{\emph{I}} \hspace{0.2cm} \textbf{Siliconised coverslips for Linbro plates (round, \& 22\,mm) are available on request.}$

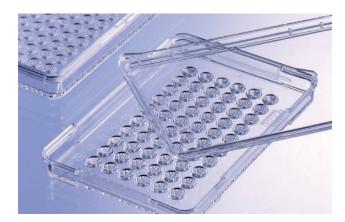

Terasaki Plates 60 and 72 Well


- / For microbatch under oil applications
- Conical well geometry
- Plate rim facilitates filling of all wells at the same time
- / With lid

Height: 83.3 mm, Length: 58 mm, Width: 10 mm, Ø well bottom: 1.3 mm, Well profile: flat, Well volume, max: 11.5 μl, Working volume (well): ≤10 μl

Item no.	Well format	Oty. inner / outer
653102	60	10 / 580
654102	72	10 / 270

 $\emph{\emph{I}} \quad \text{For Terasaki plates with TC treatment, please refer to chapter Immunology}.$

HTS-MicroplatesProtein Crystallisation Plates

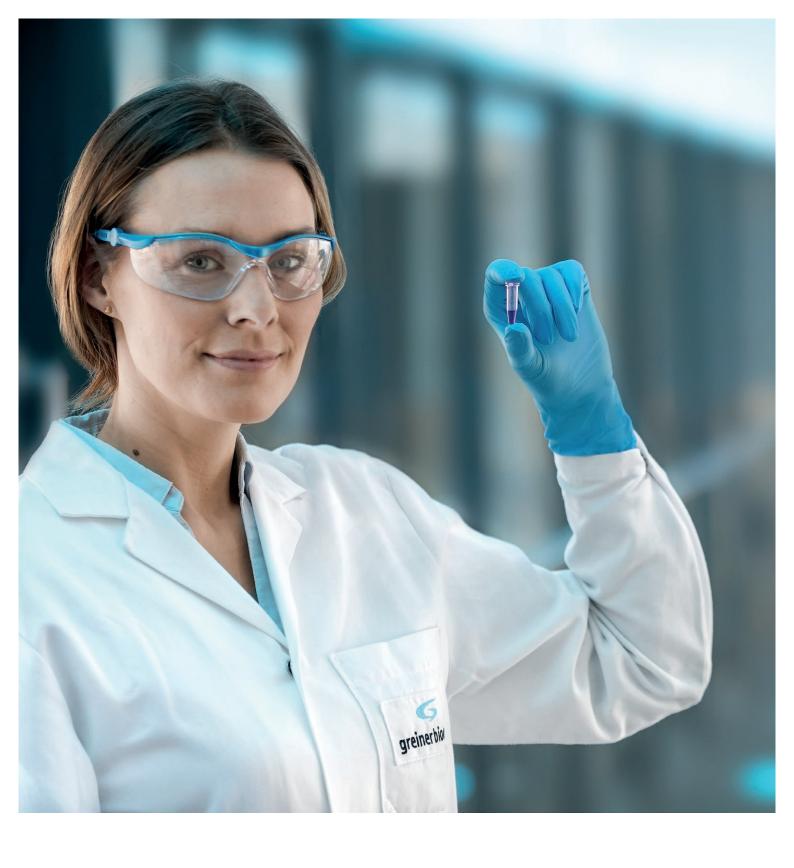

For the handling and analysis of chemical and biological samples with small or medium volumes, Greiner Bio-One offers a wide range of different analyser cups and reaction vessels with corresponding closures and cuvettes.

REACTION TUBES / ANALYSER CUPS

/	Reaction Tubes Analyser Cups .	208
	Reaction Tubes	209
	Reaction Tubes 5 ml	209
	Closures for Reaction Tubes	210
	Analyser Cups	21
/	Semi-micro / Macro Cuvette	212
	Sami-miero / Maero Cuvetto	217





/ For the max. relative centrifugal force (RCF), please refer to the technical appendix.

REACTION TUBES ANALYSER CUPS


Reaction tubes from Greiner Bio-One are available in many different versions. They have a high chemical and temperature resistance and are available from 0.5 ml to 5 ml for small and medium volumes. The reaction tubes are made of high-quality polypropylene and are available with or without lid. The lids of the corresponding vessels are flat and frosted to ensure easy labelling.

Brown reaction tubes have been specially developed for working with light-sensitive sample material. For working with liquid volumes from 1 to 5 ml, Greiner Bio-One offers 5 ml reaction tubes with screw cap

or snap cap. The tubes are also made of highly transparent polypropylene and allow contamination-free access when pipetting medium sample volumes. They also have a frosted, writable surface. This ensures reliable sample identification in everyday laboratory work.

Analyser cups are available with conical or flat bottom design and with or without cap. All analyser cups are manufactured of highly transparent polystyrene. The volumes range from 1.7 ml up to 25 ml. We offer analyser cups that are suitable for the analytical systems of Hitachi or Coulter/Hycel.

- / For volumes from 0.5 ml up to 5 ml
- / With or without cap
- / With or without graduation

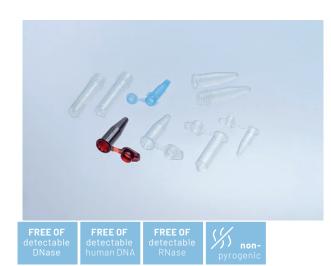

Reaction Tubes

- High chemical resistance and temperature tolerance
- Flat, frosted lids and caps for easy labelling
- Brown tube for light-sensitive materials

Description: reaction tube

Item no.	Feature	Support skirt	Graduation	Product colour	Volume range	Caps attached	Cap design	Qty. inner / outer
667201	Vitatron		-	○ natural	≤0.5 ml	yes	snap cap	1,000 / 10,000
616201	universal		yes	onatural on the same of the sa	≤1.5 mI	yes	snap cap	500 / 4,000
616283	universal		yes	brown	≤1.5 ml	yes	snap cap	500 / 4,000
623201	universal		yes	onatural on the same of the sa	≤2 ml	yes	snap cap	500 / 4,000
618201		no	-	○ natural	≤1.5 mI	no	-	500/3,000
716201		no	yes	onatural on the same of the sa	≤1.5 mI		screw cap	500/5,000
717201		yes	yes	○ natural	≤1.5 mI		screw cap	500/5,000
722201		yes	yes	onatural on the same of the sa	≤2 ml		screw cap	500/5,000

 ${\it I}$ Coloured versions of item no. 616201 and 623201 as well as sterile versions of all tubes are available on request.



Reaction Tubes 5 ml

5 ml reaction tubes with either a snap cap for fast opening and closing or a screw cap for secure sample storage.

- Outer diameter and shape similar to 15 ml tube
- Resealable packaging

Description: reaction tube, Support skirt: no, Graduation: yes, Raw material: PP

Item no.	Product colour	Cap colour	Caps attached	Cap design	Sterile	Qty. inner / outer
725201	○ natural	blue		screw cap		100 / 500
725261	natural	blue		screw cap	+	50 / 500

Item no.	Product colour	Cap colour	Caps attached	Cap design	Sterile	Qty. inner / outer
622201	○ natural		yes	snap cap		100 / 1,000
622261	onatural		yes	snap cap	+	100 / 1,000



Closures for Reaction Tubes

- / Suitable for item no. 716XXX, 717XXX and 722XXX
- / Available in natural, red, blue, green and yellow
- With gasket

Description: screw cap, Feature: with sealing ring, Suitable for tubes Ø: 12 mm, Cap design: screw cap

Item no.	Product colour	Qty. inner / outer
366380	○ natural	500 / 5,000
366383	● red	500 / 5,000
366384	blue	500 / 5,000
366385	● green	500 / 5,000
366386	yellow	500 / 5,000



Analyser Cups

- / For handling and analysis of chemical and biological samples
- / Available for different analytical systems

Description: analyser cup, Raw material: PS, Caps attached: no

Item no.	Feature	Support skirt	Bottom shape	Volume range	Lid type	Qty. inner / outer
668102	Coulter / Hycel	no	flat	≤25 ml	with lid	250 / 1,250
729101	Hitachi	yes	conical	≤1.7 ml	without lid	250 / 5,000





/ For further information, please refer to our website: www.gbo.com



SEMI-MICRO / MACRO CUVETTE

Cuvettes are used for UV/VIS spectroscopy. Greiner Bio-One offers semi-micro cuvettes with a total volume of 1.6 ml and macro cuvettes with a total volume of 4 ml. Both versions are made of crystal-clear polystyrene and are particularly suitable for enzymatic determinations, as the

very thin wall thickness enables fast and even temperature control. Our cuvettes are characterised by their low light scattering combined with a high transmission rate. Semi-micro and macro cuvettes can be used in a wavelength range from 340 nm to 900 nm.

- / Available as semi-micro or macro cuvettes
- / Very thin walls for fast and even temperature control
- / Low light scattering and high transmission
- / Suitable for a wavelength range from 340-900 nm



Semi-micro / Macro Cuvette

- / Ideal for enzymatic determinations
- / Manufactured from crystal clear polystyrene
- / Minimal light scatter but high transmission rate
- / Applicable wavelength from 340 to 900 nm

Height: 45 mm, Length: 12.5 mm, Width: 12.5 mm, Pathlength: 10 mm, Raw material: PS


Item no.	Description	Working volume	Total volume	Qty. inner / outer	
613101	semi-micro cuvette	≥0.95 ml	1.6 ml	100 / 1,000	
614101	macro cuvette	≥2.5 ml	4 ml	100 / 1,000	



The ELISA technique (Enzyme-Linked Immunosorbent Assay) is one of the most widely used biochemical methods in analytical laboratories and in diagnostics. With this method, analytes such as peptides, proteins, antibodies and hormones are detected and quantified selectively and in low concentrations. In addition, ELISAs can be automated and carried out with a high sample throughput, are relatively inexpensive and the test results are available promptly.

IMMUNOLOGY / HLA

/	ELISA Microplates216
	96 Well ELISA Microplates U-Bottom
	/ V-Bottom / F-Bottom / Half Area217
/	ELISA Strip Plates218
	96 Well ELISA Strip Plates219
	Single-break Strip Plates220
/	Immuno Tubes221
	Immuno Tubes222
/	Terasaki Plates
	Torocoki Platos

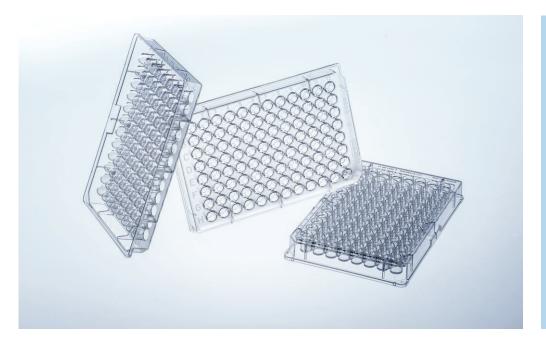



Watch our video:



ELISA MICROPLATES

Greiner Bio-One has been manufacturing microplates for diagnostics and immunological research for over 30 years.

The microplate footprint is compatible with automated systems. A key step in ELISA is the binding of one assay component – antigen or antibody – to the solid surface by passive adsorption. Therefore, the features of this surface are crucial for the performance of the assay. All ELISA microplates from Greiner Bio-One are made out of

high-quality virgin polystyrene. The resin is highly transparent and therefore ideally suited for optical measurements.

Beside products made of clear polystyrene for colorimetric measurements, Greiner Bio-One offers a wide variety of black and white ELISA microplates for luminescence and fluorescence measurements.

The quality of our immunological products is constantly controlled in our quality assurance laboratory by means of ELISA.

- / With U-Bottom / V-Bottom / F-Bottom standard (ST) / F-bottom chimney well /half area well geometry
- / With medium and high-binding surfaces
- / Made of clear / black /
 white polystyrene for
 transmission, fluorescence or luminescence
 measurements




96 Well ELISA Microplates U-Bottom / V-Bottom / F-Bottom / Half

/ Manufactured from crystal clear polystyrene

Well format: 96, Bottom: solid, Raw material: PS, Lid: no

Item no.	Well profile	Binding characterisitc	Binding brand name	Product colour	Plate design	Working volume (well)	Qty. inner / outer
650001	U-bottom	med. binding	MICROLON 200	○ clear		40 µl - 280 µl	10 / 40
650061	U-bottom	high-binding	MICROLON 600	○ clear		40 µl - 280 µl	10 / 40
651001	V-bottom	med. binding	MICROLON 200	○ clear		40 µl - 200 µl	10 / 40
651061	V-bottom	high-binding	MICROLON 600	○ clear		40 µl - 200 µl	10 / 40
655001	F-bottom/ ST	med. binding	MICROLON 200	○ clear		25 μΙ - 340 μΙ	10 / 40
655061	F-bottom/ ST	high-binding	MICROLON 600	○ clear		25 μΙ - 340 μΙ	10 / 40
655080	F-bottom / Chimney Well	med. binding	MICROLON 200	○ clear		25 μΙ - 340 μΙ	10 / 40
655081	F-bottom / Chimney Well	high-binding	MICROLON 600	○ clear		25 μΙ - 340 μΙ	10 / 40
675001	F-bottom	med. binding	MICROLON 200	○ clear	half area	15 μΙ - 175 μΙ	10 / 40
675061	F-bottom	high-binding	MICROLON 600	○clear	half area	15 µl - 175 µl	10 / 40

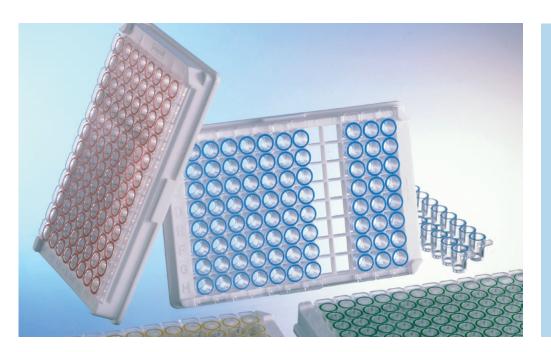



Watch our video:




ELISA STRIP PLATES

Greiner Bio-One is offering a large variety of 96 well strip microplates for diagnostic and immunological research applications.

Microplates in strip format offer the advantage of greater flexibility in diagnostics. Individual strips can be removed from the support frame so that the number of tests to be performed can be adjusted to the number of samples and is not predetermined by the microplate format used. In addition, the individual strips can be subjected to a wide variety of different test conditions. The quality of our immunological products is constantly

controlled in our quality assurance laboratory by means of ELI-SA. Clear MICROLON products are tested in a colorimetric immunoassay, black FLUOTRAC plates in a fluorescence immunoassay, white LUMITRAC plates in a luminescence immuno assay. The portfolio is completed by C8 single-break strip plates for immunological applications. These are supplied as twelve 8-well strips in a support frame with 96 spaces (12 x 8 matrix). The individual wells can be broken off separately ("single-break" option) and the number of tests performed can thus be precisely adjusted to the number of samples.

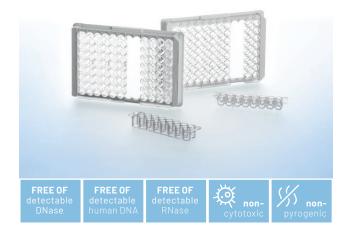


96 Well ELISA Strip Plates

- / Strips mounted in frame
- / Black and white F16 and U16 strip plates on request

Well format: 96, Bottom: solid, Raw material: PS, Lid: no

Item no.	Well profile	Binding characterisitc	Binding brand name	Product colour	Plate design	Working volume (well)	Oty. inner / outer
767070	U-bottom	med. binding	MICROLON 200	○clear	strip plate 12x8	50 µl - 280 µl	10 / 100
767071	U-bottom	high-binding	MICROLON 600	○ clear	strip plate 12x8	50 µl - 280 µl	10 / 100
762070	F-bottom	med. binding	MICROLON 200	○ clear	strip plate 12x8	20 μΙ - 350 μΙ	10 / 100
762071	F-bottom	high-binding	MICROLON 600	○ clear	strip plate 12x8	20 μΙ - 350 μΙ	10 / 100
762075	F-bottom	med. binding	LUMITRAC 200	○white	strip plate 12x8	20 μΙ - 350 μΙ	10 / 100
762074	F-bottom	high-binding	LUMITRAC 600	○white	strip plate 12x8	20 μΙ - 350 μΙ	10 / 100
762076	F-bottom	med. binding	FLUOTRAC 200	● black	strip plate 12x8	20 μΙ - 350 μΙ	10 / 100
762077	F-bottom	high-binding	FLUOTRAC 600	● black	strip plate 12x8	20 μΙ - 350 μΙ	10 / 100
756070	F-bottom	med. binding	MICROLON 200	○clear	strip plate 6x18	20 μΙ - 350 μΙ	10 / 100
756071	F-bottom	high-binding	MICROLON 600	○ clear	strip plate 6x18	20 μΙ - 350 μΙ	10 / 100
754070	U-bottom	med. binding	MICROLON 200	○ clear	strip plate 6x18	50 µl - 280 µl	10 / 100
754061	U-bottom	high-binding	MICROLON 600	○ clear	strip plate 6x18	50 µl - 280 µl	10 / 100

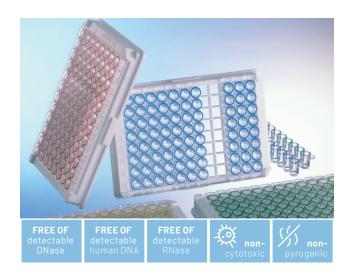

Single-break Strip Plates


- / Clear with / without colour coding
- / Strips mounted in frame
- / C8 Single-break strip plates without colour coding cell culture treated on request

Well format: 96, Well profile: C-bottom, Bottom: solid, Raw material: PS, Plate design: strip plate 12x8, Working volume (well): 20 μl - 300 μl, Lid: no

Item no.	Binding characterisitc	Binding brand name	Colour stripe	Colour coding well rim	Oty. inner / outer
705070	med. binding	MICROLON 200	○ clear		10 / 100
705071	high-binding	MICROLON 600	○ clear		10 / 100
705063	med. binding	MICROLON 200	○ clear	● red	10 / 100
705073	high-binding	MICROLON 600	○ clear	● red	10 / 100
705074	high-binding	MICROLON 600	○ clear	blue	10 / 100
705065	med. binding	MICROLON 200	○ clear	● green	10 / 100
705075	high-binding	MICROLON 600	○ clear	● green	10 / 100
705066	med. binding	MICROLON 200	○ clear	yellow	10 / 100
705076	high-binding	MICROLON 600	○ clear	yellow	10 / 100



/ More tubes can be found in the chapter Tubes / Multipurpose containers

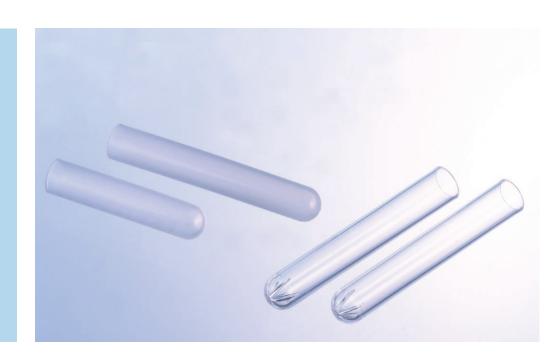

IMMUNO TUBES

Immuno tubes are often used for determining hormone levels, for example TSH (thyroid stimulating hormone). With a length of 75 mm and a diameter of 12 mm with and without a "star", they are available in both MICROLON 200 and MICROLON 600 quality. The so-called "star" at the bottom of

the tubes serves to increase the surface and thus makes it possible to bind larger amounts of antigens or antibodies.

The quality of our immunological products is constantly controlled in our quality assurance laboratory by means of ELISA.



Immuno Tubes


/ More tubes can be found in the chapter Tubes / Multipurpose containers

Height: 75 mm, Ø: 12 mm, Raw material: PS, Nominal volume: 5 ml

Item no.	Binding characterisitc	Binding brand name	Bottom shape	Oty. inner / outer
115001	med. binding	MICROLON 200	round	250 / 2.000
115061	high-binding	MICROLON 600	round	250 / 2.000
115070	med. binding	MICROLON 200	round with star	250 / 2.000
115071	high-binding	MICROLON 600	round with star	250 / 2.000



/ We provide Terasaki plates with one years' stability on the surface treatment.

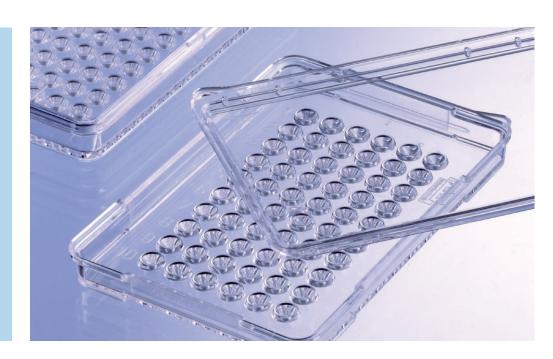

TERASAKI PLATES

Greiner Bio-One Terasaki plates are suitable for all applications for serological determination of HLA antigens. The plates are supplied either with 60 or 72 wells, the plate dimensions stay the same.

The human leucocyte antigen (HLA) system is the major histo-compatibility complex (MHC) of humans and is composed of the two polymorphic classes HLA-I (A, B, and C) and HLA-II (DR, DQ, and DP). Basically, four different areas of indication can be distinguished for HLA typing: transplantation, transfusion, disease association and

forensic.

The serological determination of HLA proteins of the HLA-A, -B, -C and -DR genetic positions is primarily performed with the complement-dependent microlymphocytotoxicity test(LCT) or Terasaki test, which has been standardised since 1964. The basis for this test method is the cytolysis of the lymphocytes to be tested, which is caused by the antibody-antigen mediated activation of the complement system. Permeabilised lymphocytes are generally stained with chromophores or fluorophores and evaluated microscopically.



Terasaki Plates

- / 60 well and 72 well Terasaki plates
- / Manufactured from crystal clear polystyrene
- / Item 659180 contains one lid per bag



Surface treatment: TC, Well volume, max: 11.5 μl, Working volume (well): ≤10 μl, Lid: yes

Item no.	Well format	Stackable	Qty. inner / outer
653180	60	no	10 / 270
653190	60	no	120 / 480
659180	60	yes	10 / 200
659190	60	yes	150 / 1,200
654180	72	no	10 / 270
769190	72	yes	150 / 1.200


Immunology / HLA Terasaki Plates



PCR (Polymerase Chain Reaction) is an important molecular biological method to determine processes in living organisms at the molecular level or genetic changes in the genetic material. It is used for the detection of diseases and parentage relationships as well as in forensics.

MOLECULAR BIOLOGY

/	Sapphire PCR Tubes228
	PCR Tubes229
	PCR 8-Tube Strips230
	Sapphire PCR 8-Cap Strips231
/	Sapphire PCR Microplates232
	96 Well PCR Microplates233
	384 Well PCR Microplates233

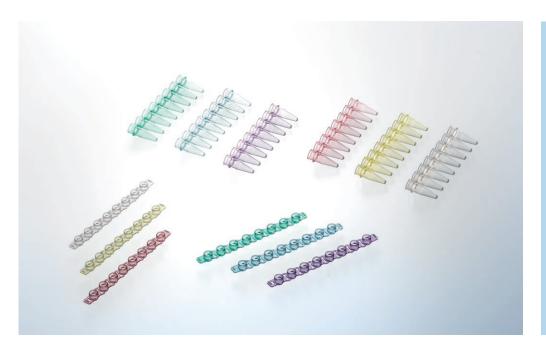



/ For standard reaction tubes please refer to chapter reaction tubes / analyser cups.

SAPPHIRE PCR TUBES

The Sapphire PCR tube portfolio comprises a full range of PCR reaction tubes (0.2 ml and 0.5 ml) as well as PCR 8-tube strips and the matching 8-cap strips.

The thin wall construction optimises the heat transfer from the block to the reaction solution. PCR tubes with flat caps are suitable for real-time / quantitative PCR and have frosted caps for easy labelling.

The classical Sapphire PCR 8-tube strips are also available

with individually attached caps. The angular attached domed or flat caps can be closed and opened individually minimising the risk of cross contamination. Moreover, individual closing and opening helps to avoid pipetting errors and facilitates sample processing.

PCR 8-tube strips are also available as **low profile tubes** with a well volume of 0.1 ml ideal for FAST, standard and real-time PCR supporting the reduction of cycle time without compromises in accuracy and efficiency.

- Manufactured out of medical-grade polypropylene without interfering additives
- / Thin-walled for efficient heat transfer during amplification
- / Comprehensive thermal cycler compatibility



PCR Tubes

Tubes with flat caps are suitable for realtime/ quantitative PCR and have frosted caps for easy labelling.


/ Packaging labelled with product symbols

Description: PCR tubes, Raw material: PP

Item no.	Product colour	Volume	Volume range	Caps attached	Cap design	Lid type	Qty. inner / outer
671201	○ natural	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671273	● red	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671274	blue	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671275	● green	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671276	yellow	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671277	■ violet	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671281	sorted	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
683201	onatural	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683273	●red	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683274	blue	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683275	● green	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683276	yellow	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683277	■ violet	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683271	sorted	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
682201	○ natural	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682273	● red	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682274	blue	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000

Item no.	Product colour	Volume	Volume range	Caps attached	Cap design	Lid type	Qty. inner / outer
682275	● green	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682276	yellow	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682277	■ violet	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682281	sorted	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
684201	natural	0.2 ml	≤0.2 ml	no		without lid	1,000 / 10,000

/ Sorted colour contains the following colours: red, blue, green, yellow, violet

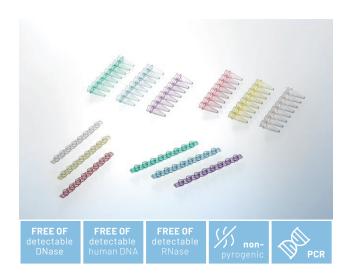

PCR 8-Tube Strips

- / Thin wall for optimal heat transfer
- / Low evaporation rate in PCR
- / Also available with individually attached caps or as low-profile version
- / Packaging labelled with product symbols

Raw material: PP

Item no.	Description	Feature	Product colour	Volume range	Caps attached	Cap design	Lid type	Oty. inner / outer
673210	PCR 8-Tube Strips		○natural	≤0.2 ml	no	separate cap strip		125 / 1,250
673273	PCR 8-Tube Strips		● red	≤0.2 ml	no	separate cap strip		125 / 1,250
673274	PCR 8-Tube Strips		blue	≤0.2 ml	no	separate cap strip		125 / 1,250
673275	PCR 8-Tube Strips		green	≤0.2 ml	no	separate cap strip		125 / 1,250
673276	PCR 8-Tube Strips		yellow	≤0.2 ml	no	separate cap strip		125 / 1,250
673277	PCR 8-Tube Strips		● violet	≤0.2 ml	no	separate cap strip		125 / 1,250
673271	PCR 8-Tube Strips		sorted	≤0.2 ml	no	separate cap strip		125 / 1,250
671221	Low profile 8-tube strips	low profile	onatural	≤0.1 ml	yes	indi- vidually attached	flat	120 / 1,200

Item no.	Description	Feature	Product colour	Volume range	Caps attached	Cap design	Lid type	Oty. inner / outer
673283	8-tube strips with individually attached caps		○natural	≤0.2 ml	yes	indi- vidually attached	domed	120 / 1,200
608281	8-tube strips with individually attached caps		○ natural	≤0.2 ml	yes	indi- vidually attached	flat	120 / 1,200
673281	Low profile 8-tube strips	low profile	○ natural	≤0.1 ml	no	separate cap strip	flat	125 / 1,250

[/] Sorted colour contains the following colours: red, blue, green, yellow, violet




Sapphire PCR 8-Cap Strips

- / Real-time PCR cap strips (Item No. 373250) have a flat lid and are made of highly transparent polypropylene
- / These are ideal for real-time PCR applications and are compatible with most real-time PCR devices

Description: 8-cap strip, Raw material: PP

Item no.	Feature	Product colour	Qty. inner / outer
373270		○ natural	125 / 1,250
373273		● red	125 / 1,250
373274		blue	125 / 1,250
373275		● green	125 / 1,250
373276		yellow	125 / 1,250
373277		◆ violet	125 / 1,250
373281		sorted	125 / 1,250
373250	for RT PCR	natural	125 / 1,250

 $^{{\}it I} \quad {\tt Sorted \, colour \, contains \, the \, following \, colours: \, red, \, blue, \, green, \, yellow, \, violet}$

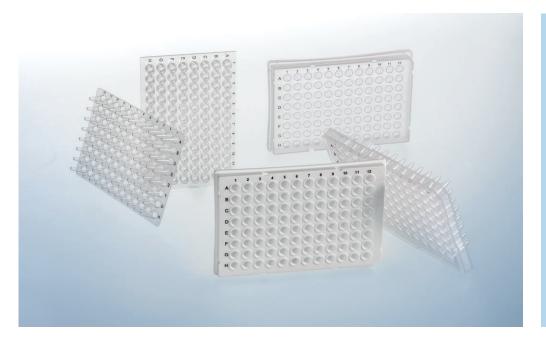



/ For details regarding plate geometry as well as compability with thermal cyclers, please refer to the technical appendix.


SAPPHIRE PCR MICROPLATES

The use of the 96 well format allows the scale up of basic PCR work, while the 384 well format is ideal for high-throughput screening projects.

All microplates are made of thinwalled polypropylene. This op-

timises the heat transfer from the thermoblock to the reaction solution. Our heat-resistant sealers AMPLIseal, VIEW-seal and SILVERseal are ideal for sealing the microplates during PCR, and the 96 well microplate may also be sealed with 8-cap strips.

- / Available in 96 well and 384 well format
- / Ultra-thin polypropylene for optimal heat transfer
- / Sealable with sealers SILVERseal, VIEWseal and AMPLIseal

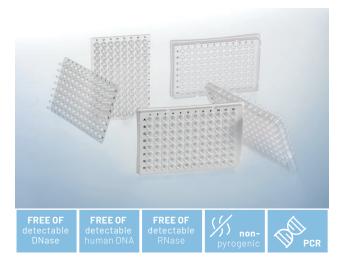


96 Well PCR Microplates

- / Without skirt, half-skirt and full-skirt
- Alphanumeric printing for rapid sample identification
- / Sealable with sealers SILVERseal , VIEWseal and AMPLIseal or with compatible 8-cap strips
- / Low profile microplate for reduced sample volume and cycle time

Well format: 96, Raw material: PP

Item no.	Feature	Skirt	Product colour	Colour alphanumeric code	Well volume, max	Qty. inner / outer
652201		without skirt	○natural	● black	0.2 ml	10 / 100
652250	flat, universal	without skirt	natural	● black	0.2 ml	10 / 100
652210	low profile	without skirt	○ natural	● black	0.1 ml	10 / 100
669285	suitable for LightCycler®	half-skirt	○white	● black	0.2 ml	10 / 100
652290	suitable for ABI	half-skirt	○ natural	● black	0.2 ml	10 / 100
652260	ABI design	half-skirt	natural	● black	0.2 ml	10 / 100
652270		yes	○ natural	● black	0.2 ml	10 / 100

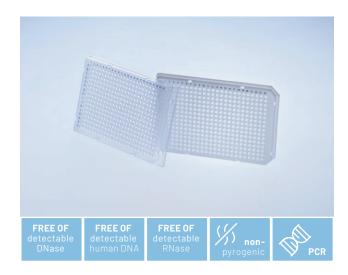


384 Well PCR Microplates

- / Thin wall for optimal heat transfer
- / Alphanumeric well coding
- / Sealable with sealers SILVERseal , VIEWseal and AMPLIseal
- / White and black microplate versions are available on request

Well format: 384, Skirt: yes, Raw material: PP, Total volume (well): 25 μl

Item no.	Feature	Product colour	Colour alphanumeric code	Qty. inner / outer		
785201		○ natural	blue	15 / 60		
785290	suitable for ABI	natural	blue	15 / 60		
785285	suitable for LightCycler®	○white	● black	50 / 100		

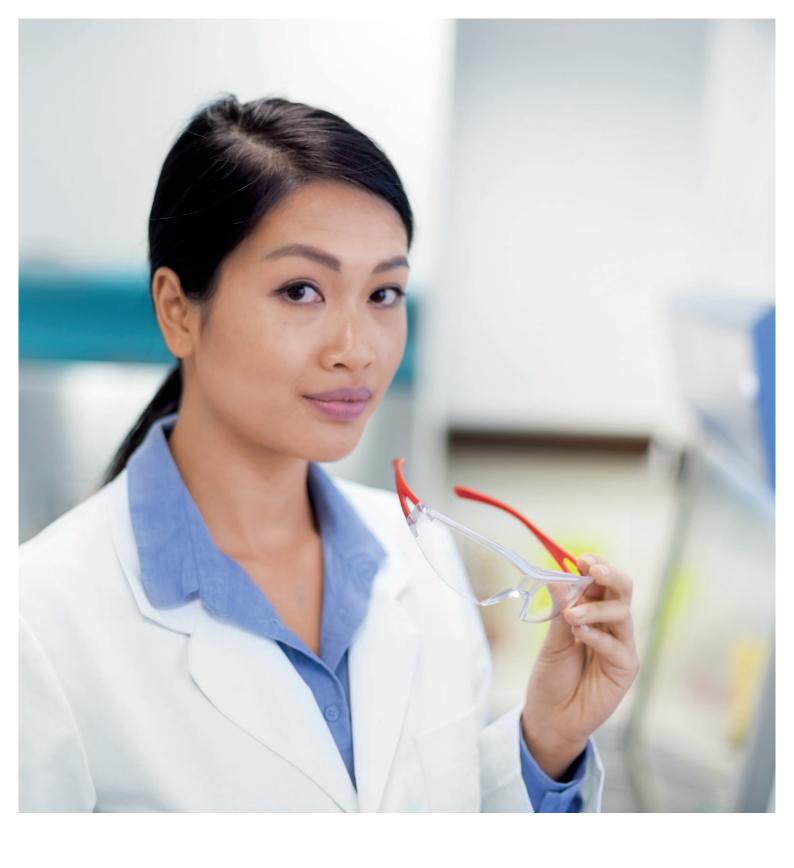

The Greiner Bio-One portfolio includes small laboratory equipment to match the consumables such as PCR tubes, plates and tubes, in order to provide our customers with comprehensive service, compatibility and highest quality: Mini-centrifuges, Vortex Mixers, Block Heaters, Microplate Centrifuges and ergonomic pipettes are used for modern handling in everyday laboratory work. All from one source - for highest quality requirements.

LAB EQUIPMENT

/	Lab Equipment
	Mini Centrifuge237
	Microplate Centrifuge237
	Mini Vortex Mixer238
	Vortex Mixer238
	Mini Block Heater239
	Inserts for Mini Block Heater239





- / For Sapphire MaxiPette pipette controller please refer to chapter Liquid Handling
- / For further information, please refer to our website: www.gbo.com

LAB EQUIPMENT

Greiner Bio-One, a leading provider of labware, completes its product portfolio by compatible lab equipment.

Mini centrifuges are ideal for quick spin-downs of microtubes and PCR tubes. Braking system is activated when the lid is opened, smoothly decelerating to a complete stop in just one second.

The microplate centrifuge allows centrifugation of two plates, even without a lid or seal.

Our mini vortex mixer can vortex even the largest samples instantly with its 4 mm orbit and speed of 2800 rpm. Despite its powerful motor, it has a size smaller than 10×10 cm, allowing it to fit on even the most crowded bench.

The Mini Bock Heater is the perfect tool for labortory incubations, with its compact construction, broad temperature range and multiple block options.

- / Centrifuges
- / Vortex Mixer
- / Mini Block Heater



Mini Centrifuge

- / Ideal for quick spin-downs of microtubes and PCR tubes
- / Rotor instantly reaches 6,000 rpm
- / Storage compartment for PCR tube rotor
- / Decelerates in just one second
- / Compatible with all Greiner Bio-One PCR and reaction tubes

Feature: capacity 8 x 1.5 / 2.0 ml tubes, 4 x PCR strips (8 x 0.2 ml), 32 x 0.2 ml PCR tubes, Height: 11.4 cm, Length: 11.4 cm, Width: 15 cm, Speed: 6,000 RPM, Operating temperature: 4 - 45 °C, Max. radius: 4.9 cm

Item no.	Plug	Qty. inner / outer
843070	EU	1/1



Microplate Centrifuge

- / Rapid spin-down of microplates with 2,550 rpm
- / Unique rotor design holds two plates without spillage
- / 50 % smaller than traditional centrifuges
- / Compatible with most Greiner Bio-One microplates

Description: Microplate Centrifuge, Feature: capacity x 96 well microplates, 2 x 96 well PCR microplates, Height: 19.7 cm, Length: 23 cm, Width: 26 cm, Speed: 2,550 RPM, Operating temperature: 4 - 45 °C, Max. radius: 8 cm

Item no.	Plug	Qty. inner / outer
846070	EU	1/1

/ Device with US or UK plug available on request



Mini Vortex Mixer

- / Powerful, quiet vortexing for tubes up to 50 ml
- / Instant pressure-activated operation
- / Compact, fits in the palm of your hand
- / Compatible with all Greiner Bio-One tubes

Description: Mini Vortex Mixer, Feature: head: Standard rubber lid for mixing single tubes, Height: 6.6 cm, Length: 9.4 cm, Width: 9.9 cm, Mode: TOUCH Mix, Orbit: 4 mm, Operating temperature: 4 - 45 °C

Item no.	Plug	Oty. inner / outer
845070	EU	1/1

/ Device with US or UK plug available on request



Vortex Mixer

- / Powerful motor for instant vortexing of up to 50 ml tubes
- / Dynamic balance system to minimise noise and vibration
- / Touch or continuous operation
- / Variable speed control from 200 to 3,200 rpm
- / Compatible with all Greiner Bio-One tubes

Description: Vortex Mixer, Feature: head: standard rubber cap single tube, base with four suction cups, Height: 17 cm, Length: 13 cm, Width: 16 cm, Mode: ON (continuous), OFF and TOUCH mix, Orbit: 3, Operating temperature: 4 - 45 °C

Item no.	Plug	Oty. inner / outer
844070	EU	1/1

/ Device with US or UK plug available on request



Mini Block Heater

- / Simple touchpad operation
- / Easy-to-read digital display
- / Compact, fits in the palm of your hand
- / Exchangeable blocks for tubes from 0.2 to 50 ml

Description: Mini block heater (without insert), block options, high-grade aluminium, Feature: ambient temperature from 5 - 100 °C, Height: 10 cm, Length: 11.2 cm, Width: 15 cm, Temperature increments: +/- 0.5 °C, Temperature uniformity: 0.2 °C, Temperature incements: 0.1 °C, Operating temperature: 4 - 45 °C

Item no.	Plug	Oty. inner / outer
848070	EU	1/1

/ Device with US or UK plug available on request



Inserts for Mini Block Heater

/ Exchangeable blocks for tubes from 0.2 to 50 ml

Item no.	Feature	Qty. inner / outer
848916	suitable for 15 pcs. 1.5 ml reaction tube, 0.5 ml PCR tube	1/1
848923	suitable for 15 pcs. 1.5 / 2 ml reaction tube, 0.5 ml PCR tube	1/1
848902	suitable for 40 pcs. 0.2 ml PCR tube, PCR 8-tube strip	1/1
848913	suitable for 15 pcs. 4 / 5 ml Cryo.s, 4.5 / 5 / 7 ml tube	1/1
848921	suitable for 15 pcs. 1 / 2 ml Cryo.s, 4 ml tube	1/1
848915	suitable for 4 pcs. 15 ml conical tube, 12 / 14 / 20 ml tube	1/1
848950	suitable for 2 pcs. 50 ml conical tube	1/1



DIN EN ISO 50001 Certification

QUALITY STANDARDS AT GREINER BIO ONE

Greiner Bio-One is certified according to the international standards DIN EN ISO 9001 and EN ISO 13485 for Medical Devices. Since 2013,

Greiner Bio-One Frickenhausen (Germany) is also certified according to DIN EN ISO 50001 (systematic energy management).

TECHNICAL APPENDIX

TECHNICAL APPENDIX

Quality Standards at Greiner Bio One240

/	General Information for the Lab	. 242
	Chemical Resistance of Various Materials	242
	Chemical Resistance of Various Materials	243
	Chemical Resistance of	
	Cycloolefins (COC/COP)	244
	Chemical Resistance of Polyethylene	
	Terephthalate(PET)	245
	Capillary Pore Membranes	
	(ThinCert® Cell Culture Inserts)	245
	Chemical Resistance of Sealers	246
	Physical Properties of Various Materials	247
/	Laboratory Information for	
	Microplates	.250
	Well Profiles of 96 Well Microplates	
	·	
/	Laboratory Information for Liquid	
	Handling	.252
	Compatibility for SAPPHIRE Pipette Tips	
	and pipettes	252
	Overview PCR Microplates	
	Compatibility for Mini Block Heater Inserts	
	, , , , , , , , , , , , , , , , , , , ,	
/	Laboratory Information for	
	Centrifugation	.260
	Centrifugation - Principle and Calculation	
	of the RCF	260
	(Relative Centrifugal Force)	
	Alignment Chart	
	Maximal Centrifuge Capacity of Tubes,	
	maximar continuage duputity of Tubes,	U _

REACTION TUBES AND MICROPLATES	262
Polystyrene Tubes	26
Polyethylene Tubes	264
Microplates	26
Laboratory Information for Cryo.s	
Sample Storage	266
Laboratory Information for	
Immunology	268
Volume-Dependent Wetting of	
Immunological Products	268
	Polystyrene Tubes Polyethylene Tubes Microplates Laboratory Information for Cryo.s Sample Storage Laboratory Information for Immunology Volume-Dependent Wetting of

GENERAL INFORMATION FOR THE LAB

CHEMICAL RESISTANCE OF VARIOUS MATERIALS

	PS 20°C	PS 50°C	PP 20 °C	PP50°C	HDPE 20°C	HDPE 50°C	LDPE 20°C	LDPE 50°C	PC 20°C	PC 50°C
Acetic acid 10 %	1	1	1	1	1	1	1	1	1	2
Acetic acid 50 %	2	2	1	1	1	1	1	1	1	2
Acetic acid 90 %	4	4	1	2	1	1	1	2	4	4
Acetone	4	4	1	3	1	1	3	3	4	4
Acetonitrile	4	4	3	4	1	1	1	1	4	4
Ammonia 25 %	2	2	1	1	1	1	1	1	4	4
Ammonium acetate	1	1	1	1	1	1	1	1	1	1
Amyl alcohol	1	1	1	1	1	1	1	2	1	
Ascorbic acid			1	1	1		1		2	2
Benzene	4	4	4	4	4	4	4	4	4	4
Benzyl alcohol	4	4	4	4	3	4	4	4	4	4
Boric acid 10 %	1	1	1	1	1	1	1	1	1	1
Carbon tetrachloride	4	4	4	4	3	4	4	4	4	4
Carbonic acid	1	1	1	1	1	3	1	1	1	
Chloroform 100 %	4	4	3	4	3		3		4	4
Citric acid 10 %	1	1	1	1	1	1	1	1	1	2
Cyclohexanol	3	3	1	3	1	1	1	1	3	
Detergents			1	1						
Dichloroacetic acid			1	1	1	1			4	4
Diethyl ether	4	4	4	4	3	4	4	4	4	4
Dimethyl acetamide	4	4	1	1	1	1	3	4		
Dimethylsulfox. (DMSO)	1	2	1	1	1	1	1	1	4	4
Emulsifier	·	_	1	1	·	-	-	-	•	·
Ethanol 50 %	1	1	1		1	1	1	2	1	1
Ethanol 96 %	1	1	1	1	1		1		1	3
Ether	4	4	4	4	3	4	4	4	4	4
Formaldehyde 10 %	3	4	1	1	1	1	1	1	1	2
Formaldehyde 40 %	4	4	1	2	1	2	2	3	1	2
Formamide	1	1	1	1	1	1	1	1	3	3
Formic acid 50 %	3	3	1	2	1	1	1	2	3	3
Glucose	1	1	1	1	1	1	1	1	1	1
Glycerine	1	1	1	1	1	1	1	1	3	3
Heptane	4	4	3	3	2	3	3	4	1	2
Hexanol			1		1		1		2	2
Hydrochloric acid 20 %	1	1	1	1	1	1	1	1	2	3
Hydrochloric acid conc.	3	3	1	1	1	1	1	1	4	4
Hydrogen peroxide 3 %	1	1	1	1	1	1	1	1	3	3
Hydroquinone	4	4	1	•	·	·	1	3	3	3
Isoamyl alcohol	1	1	•				·	Ü	3	0
Isobutanol	2	2	1	1	1	1	1	1	1	2
Isopropanol	2	2	1	1	1	1	1	1	1	2
Isopropyl acetate	4	4	2	3	1	2	2	3	4	4
Isopropyl benzene	4	4	3	4	2	3	3	4	4	4
Isopropyl ether	4	4	4	4	4	4	4	4	4	4
Lactic acid 3 %	2	2	1	2	1	1	1	2	1	2
Lactic acid 85 %	2	2	1	2	1	1	1	1	1	2
Liquid paraffin	1	1	1	3	1	1	1	3	1	_
ridaia hai ai iiii	'	1	'	J	'	ı .		J	1	

CHEMICAL RESISTANCE OF VARIOUS MATERIALS

	PS 20°C	PS 50°C	PP 20°C	PP50°C	HDPE 20°C	HDPE 50°C	LDPE 20°C	LDPE50°C	PC 20°C	PC 50°C
Methanol	3	4	1	1	1	1	1	1	4	4
Methyl acetate	4	4	2	3	3	3	3	4	4	4
Methyl phenyl ether 100%	4	4	3				3		4	4
Methyl propyl ketone	4	4	2	3	1	2	2	3	4	4
Methylamine 32 %			1		1		1		4	4
Methylene chloride	4	4	3	4	4	4	4	4	4	4
Naphthalene			1		1	3			3	3
Nitrobenzene	4	4	4	4	3	4	4	4	4	4
Oxalic acid	1	1	1	1	1	1	1	1	3	4
Ozone	3	3	1	2	1	1	1	2	1	2
Palmitic acid	1	1	3	4	3		2		2	2
Phenol 10 %	4	4	1	1	1	1	1	1	4	4
Phenol 100 %	4	4	1	1	2	3	3	3	4	4
Phosphoric acid 1 - 5 %	2	2	1	1	1	1	1	1	1	1
Phosphoric acid 85 %	1	1	1	2	1	1	1	1	1	2
Phthalic acid	1	1	1	1	1	1	1	1	3	3
Potassium carbonate	1	1	1	1	1	1	1	1	3	3
Potassium chromate	1	1	1	1	1	1	1		2	2
Potass. permanganate	2	3	1	1	1	1	1	1	1	
Propanol	3	3	1	1	1	1	1	1	1	
Sodium acetate	2	2	1	1	1	1	1	1	1	2
Sodium chloride	1	1	1	1	1	1	1	1	1	1
Sodium hydroxide 30 %	1	1	1	1	1	1	1	1	4	4
Sodium hydroxide 45 %	1	1	1	1	1	1	1	1	4	4
Sodium hydroxide 60 %	1	1	1	1					4	4
Sodium hypochloride	1	1	2	3	2	3	2	3	2	3
Sodium permanganate	2	3	1	1	1	1	1	1		
Sodium thiosulfate	1	1	1	1	1	1	1	1	2	2
Stearic acid	1	2	1	1	1	1	1	1	1	2
Sulphuric acid 1 - 6 %	1	2	1	1	1	1	1	1	1	1
Sulphuric acid 60 %	2	4	1	3	1	3	1	3	3	3
Sulphuric acid conc.	4	4	4	4	4	4	4	4	4	4
Tannin acid	1	1	1	1					3	3
Terpentine oil					3	4	3	4	4	4
Tetrahydrofuran	4	4	3	4	3	4	4	4	4	4
Toluene	4	4	3	4	3	4	3	4	4	4
Trichloroacetic acid	4	4	3	4	3	3	3	4	4	4
Urea	1	2	1	1	1	1	1	1	1	1
Uric acid			1		1		1		1	
Urine	3	3	1	1	1	1	1	1	1	
Xylene	4	4	4	4	2	3	2	4	4	4

¹⁼resistant 2=limited resistant 3=moderate resistant 4=no resistance

 $These \ tables \ are \ a \ general \ guide \ only. \ As \ many factors \ can \ affect \ the \ chemical \ resistance \ of \ a \ given \ product, its \ suitability \ for \ a \ specific \ application$ should be tested.

CHEMICAL RESISTANCE OF CYCLOOLEFINS (COC / COP)

	Cycloolefin		Cycloolefin		Cycloolefin
Acetic acid 99 %	1	Dibutyl ether	4	Isopropanol	1
Acetone	1	Dichloroethane	4	Methanol	1
Acrylonitrile	1	Dichloromethane	4	Methylene chloride	4
Ammonia 33 %	1	Diethyl ether	4	Nitric acid (HNO ₃)	1
Benzaldehyde	3	Dimethyl sulfoxide	1	Octane	4
Benzene	4	DMS0	1	Pentane	4
Benzine	4	Ethanol 50 %	1	Sodium hydrox. (NaOH) 50 %	1
Butanon	1	Ethanol 96 %	1	Sulphuric acid (H ₂ SO ₄) 40 %	1
Carbon tetrachloride	4	Fatty acid	4	Sulphuric acid (H ₂ SO ₄) 95 %	1
Chloroform	4	Heptane (n-Heptane)	4	Toluene	4
Cyclohexane	4	Hexane	4	Xylene	4
Cyclohexanone	4	Hydrochloric acid (HCI) 36 %	1		
Detergents	1	Hydrogen peroxide water 30 %	1		

^{1 =} resistant 2 = limited resistant 3 = moderate resistant 4 = no resistance

This table is a general guide only. As many factors can affect the chemical resistance of a given product, its suitability for a specific application should be tested.

CHEMICAL RESISTANCE OF POLYETHYLENE TEREPHTHALATE (PE CAPILLARY PORE MEMBRANES (THINCERT® CELL CULTURE INSER

	PET		PET		PET
Acetaldehyde	1	Ethanol	1	Monochlorbenzene	1
Acetic acid (10 %)	1	Ethyl acetate	1	Nitric acid (30 %)	1
Acetic acid (100 %)	3	Ethyl ether	1	Nitrobenzene	1
Acetone	1	Ethylendichloride	1	Nitropropane	1
Ammonium hydroxide (5 %)	1	Ethylene glycol	1	n-Propanol	1
Amyl acetate	1	Fluoric acid (35 %)	1	Pentane	1
Amyl alcohol	1	Formaldehyde	1	Perchlorethylene	1
Aniline	1	Formic acid (50 %)	1	Petroleum ether	1
Benzene	3	Freon	1	Phosphoric acid (85 %)	3
Benzyl alcohol	1	Glutaraldehyde	1	Potassium hydroxide	4
Benzyl benzoate	1	Glycerol	1	Propyl acetate	1
Boric acid (5 %)	1	H ₂ O ₂ (30 %)	1	Pyridine	1
Butanol	1	Halogenated phenoles	4	Silicon oil	1
Butyl acetate	1	Hexane	1	Sodium hydroxide	4
Butyl cellusolve	1	Hydrochloric acid (20 %)	1	Sulphuric acid (25 %)	1
Carbon tetrachloride	1	i-Propanol	1	Terpentine oil	1
Chloroform	1	Isopropyl myristate	1	Tetrahydrofurane	1
Concentrated strong acids	4	Methanol	1	Tetraline	1
Cyclohexane	1	Methyl acetate	1	Toluene	3
Cyclohexanone	3	Methyl cellusolve	1	Trichlorbenzene	1
Dekaline	1	Methylenchloride	3	Trichlorethylene	1
Dimethylacetamide	1	Methylethylketone	1	Triethanolamin	1
Dimethylformamide	1	Methylglycol acetate	1	Trikresyl phosphate	1
Dimethylsulfoxide	1	Methylisobutylketone	1	Xylene	3
Dioxane	1	Mineral oils	1		

For the solvents effecting slight changes the user should test the compatibility under the specific application conditions. All tests have been performed at RT. Please be aware that ThinCert® cell culture inserts are made of PET membranes sealed on polystyrene housings. Therefore, solvents shown compatible with PET membranes in the above table might be incompatible with the polystyrene housing. Please check solvent compatibility with polystyrene.

Resistance scale from 1 to 4

1=resistant $i.e.\ the\ plastics\ may\ be\ treated\ with\ the\ chemical\ compound\ at\ mentioned\ temperature\ over\ several\ years$ without

any significant alterations in its physical, optical and chemical properties

2 = limited resistant i.e. the plastics may be treated with the chemical compound at mentioned temperature over several weeks without any significant alterations in its physical, optical and chemical properties

 $3 = moderate \, resistant \, i.e. \, the \, plastics \, may \, be \, treated \, with \, the \, chemical \, compound \, at \, mentioned \, temperature \, for \, short \, time \, only \, (several \, minutes \, to \, plastics \, may \, be \, treated \, with \, the \, chemical \, compound \, at \, mentioned \, temperature \, for \, short \, time \, only \, (several \, minutes \, to \, plastics \, may \, be \, treated \, with \, the \, chemical \, compound \, at \, mentioned \, temperature \, for \, short \, time \, only \, (several \, minutes \, to \, plastics \, may \, be \, treated \, with \, the \, chemical \, compound \, at \, mentioned \, temperature \, for \, short \, time \, only \, (several \, minutes \, to \, plastics \, may \, be \, treated \, with \, the \, chemical \, compound \, at \, mentioned \, temperature \, for \, short \, time \, only \, (several \, minutes \, to \, plastic \, may \, be \, treated \, with \, the \, chemical \, compound \, at \, mentioned \, temperature \, for \, short \, time \, only \, (several \, minutes \, to \, plastic \, may \, be \, treated \, with \, the \, chemical \, compound \, at \, mentioned \, temperature \, for \, short \, chemical \, compound \, at \, mentioned \, temperature \, for \, short \, chemical \, ch$ one hour)

without any alterations in physical, optical and chemical properties (mixing and measuring is possible)

4 = no resistance i.e. treating the plastics with the substance named may cause alterations in physical, optical and chemical properties within seconds

CHEMICAL RESISTANCE OF SEALERS

	EASYseal (Art. No. 676001)	VIEWseal (Art. No. 676070)	AMPLIseal (Art. No. 676040)	SILVERseal (Art. No. 676090)
Acetone	4	4	4	3
Acetonitrile	3	3	4	1
Acetic acid 1 %	3	1	4	3
Glacial acetic acid	1	3	4	3
Chloroform	4	4	4	4
DMS0	3	3	3	1
Ethanol	3	1	1	1
Hydrochloric acid 32 %	3	1	3	4
Isopropanol	3	1	1	1
Methanol	3	1	4	1
Phenol	3	3	4	3
Sulphuric acid 0.5 M	1	1	1	1

¹⁼Stable no visible change in the sealer after one week's incubation

 $This table \ can only \ be \ used \ as \ an \ orientation \ aid \ for \ the \ suitability \ of \ the \ respective \ sealers, since \ their \ behaviour \ against \ chemicals \ depends \ on \ the$ respective application. Tests under practical conditions are absolutely essential in many cases.

TEMPERATURE STABILITY OF SEALERS

	Temperature Stability
EASYseal	-40 °C to + 120 °C
VIEWseal	-70 °C to +100 °C
AMPLIseal	-80 °C to + 110 °C
SILVERseal	-70 °C to +100 °C
BREATHseal	n.a. / Evaporation rate 4200 g H ₂ 0/m² in 24 h

This table can basically be used as an orientation aid for the temperature stability of the respective sealers, since the behaviour of the product depends on the respective application. Tests under practical conditions are absolutely essential in many cases.

^{3 =} Moderately stable after one week, optical and physical changes in the sealer (clouding tears on removal)

^{4 =} Unstable adhesive and foil are dissolved, wells not leak-tight

PHYSICAL PROPERTIES OF VARIOUS MATERIALS

Material	Sterilisation by			Auto- clavability	Thermal Stability[°C]	Transparency	Gas Permeability²		WVTR ³		
	gamma irradiation	chemicals (formalin, ethanol)	dry heat	gas¹				02	N ₂	CO ₂	
Polystyrene		•		•		-20 to +60	clear	4.7	853	17.8	108 – 155
Polypropylene	•	•		•	•	-196 to +121	translucent	3.7	744	12.4	3.9
HDPE	•	•		•		-50 to +100	translucent	2.9	651	9	4.6 - 6.2
LDPE	•	•		•		-50 to +80	translucent	7.8	2.8	41.9	15.5 -23.3
UV-Star®	•			•		-20 to +40	clear				
PETG	•	•		•		-40 to +60	clear	388	155	1.2	62
PET	•	some		•		-40 to +60	clear	46.5	10.9	236	15 – 20
Cycloolefin	•			•		-80 to +100	clear				

Exemptions are mentioned in the respective product data sheets.

Material	Refractive Index
Polystyrene	1.59
UV-Star®	1.53
Cycloolefin	1.53
Glass	1.53

 $These \ tables \ are \ a \ general \ guide \ only. \ As \ many factors \ can \ affect \ the \ resistance \ of \ a \ given \ product, its \ suitability \ for \ a \ specific \ application \ should \ be \ an \ application \ should \ be \ application \ should \$ tested.

¹ Ethylene oxide, formaldehyde

 $^{^{2}}$ [cc x mm/m 2 x 24 h x Bar]

 $^{^3}$ at 37°C, 90 % humidity[g x mm/m² x 24 h $\,$ x Bar]

MANUAL CALCULATION


COEFFICIENT OF VARIATION (CV)

The coefficient of variation compares the variability of several random samples with different means, taking into account the different dimensions of means:

/ where S is the standard deviation and IXI is the absolute value of the arithmetic mean.

VOLUME OF DIVERSE BODIES

VOLUME OF A CYLINDER:

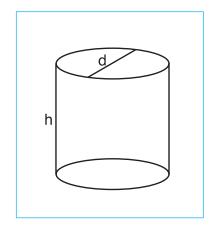


$$V = \frac{\pi \cdot d^2 \cdot h}{4}$$

$$h = \frac{4 \cdot V}{\pi}$$

/ This formula can be used for calculating the filling level in relation to the filling volume in a 96 well microplate with cylindrical wells.

VOLUME OF A CUBOID:

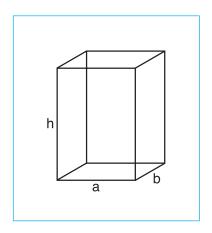

$$V = a \cdot b \cdot h$$

$$h = \frac{V}{a \cdot b}$$

/ This formula can be used for calculating the filling level in relation to the filling volume in 384 and 1536 well microplates with rectangular wells.


OVERVIEW

METRIC PREFIXES

```
G = giga
             = 10 9
M = mega = 10^6
k = kilo
             = 10^{3}
c = centi
             = 10<sup>-2</sup>
             = 10^{-3}
m = milli
\mu = micro = 10^{-6}
             = 10^{-9}
n = nano
             = 10^{-12}
p = pico
f = femto = 10^{-15}
             = 10^{-18}
a = atto
z = zepto = 10^{-21}
```

LABORATORY INFORMATION FOR MICROPLATES

WELL PROFILES OF 96 WELL MICROPLATES

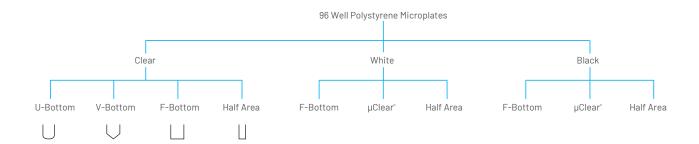

WELL PROFILE

The well profile is a critical aspect in a 96 well microplate. Different well shapes are available for each application (Fig. 1 – Fig. 4):

U-BOTTOM

The "U" describes the round bottom shape (Fig. 1). U-bottom microplates are ideally suited for agglutination tests.

- / No sharp corners to facilitate easy and residue-free pipetting
- / Suitable for +/- analyses

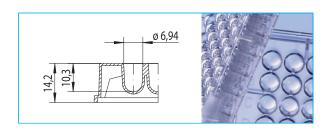


Figure 1:

Well profile: 96 well U-bottom, polystyrene

Total volume: 323 µl Working volume: 40 – 280 µl

V-BOTTOM

The "V" stands for the conically tapered well bottom (Fig. 2). These microplates are ideally suited for applications in which the entire sample volume must be pipetted off.

- / For precise pipetting
- / Ideally suited for the storage of samples
- / Suitable for +/- analyses

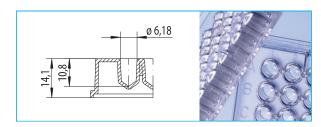


Figure 2:

Well profile: 96 well V-bottom, polystyrene

Total volume: 234 µl Working volume: 40 – 200 µl

F-BOTTOM/STANDARD (ST)

The "F" refers to the flat well bottom (Fig. 3). This well type is ideal for precise optical measurements. The measuring light source is not deflected by the well profile.

- / Excellent optical properties
- / For precise optical measurements
- For microscopic applications (bottom reading)

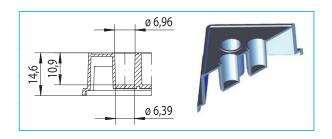


Figure 3: Well profile: 96 well F-bottom/ST, polystyrene

Total volume: 382 µl Working volume: 25 – 340 µl Growth area: 32 mm²

F-BOTTOM/CHIMNEY WELL

The standard flat bottom microplate (Fig. 3) has the same well profile as the chimney well microplate (Fig. 4). The difference from the standard plate is the chimney-like arrangement of the wells. Each well stands on its own. Therefore the risk of sample carryover and cross contamination is minimised.

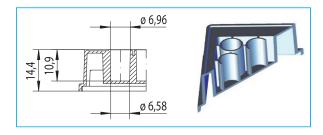


Figure 4:
Well profile: 96 well F-bottom/ chimney well,
polystyrene

Total volume: 392 µl Working volume: 25 – 340 µl Growth area: 34 mm²

LABORATORY INFORMATION FOR LIQUID HANDLING COMPATIBILITY FOR SAPPHIRE PIPETTE TIPS AND PIPETTES

		PIPETTE TIPS (STANDARD / LOW RETENTION)							
Volume [µI]		10	10 extended	20	100	200	300	1000	1250
Volume range [µI]		0.2-10	0.5-10	2-20	2-100	5-200	10-300	50-1000	50-1250
Tip version		Tip/ Filter tip	Tip/ Filter tip	Filtertip	Filtertip	Tip/ Filter tip	Tip/ Filter tip	Tip	Tip/ Filter tip
Article no.		7713XX	7723XX	7733XX	7743XX	7753XX	7763XX	7773XX	7783XX
PIPETTE MODELS									
GREINER BIO-ONE									
GREINER BIO-ONE			Single-	channel pipet	tos				
Sapphire	0.2-2 μΙ	•	• Siligle	chamile piper	163				
очррии с	1-10 µI	•	•						
	2-20 µl		, and the second	•		•			
	10-100 μΙ			•	•	T● I F-			
	20-200 µI				•	•	•		
	20-200 μl			-	-	·	•	•	
	500-5000 μI							•	
	1000-10000 µl		M. det						
) b i	8/12 CH 0.5-10 µl	•	Multi-C	channel pipett	es				
Sapphire		•	•	•		T			
	8/12 CH 2-20 µI			_	_	T● I F-	•		
	8/12 CH 20-200 μI			-	_	•	•		
	8/12 CH 20-300 μl			•		•	•		
BRAND									
			Single-	channel pipet	tes				
Transferpette® S	0.5-10 μΙ	•	•						
	2-20 μΙ	•	•	•		T● I F-			
	10-100 μΙ			•		T● I F-			
	20-200 μΙ				•	•	•		
	100-1000 μΙ							•	T● I F-
EPPENDORF									
			Single-	channel pipet	tes				
Reference®	0.5-10 μl	•	•	onamici piper					
	2-20 µl			•	•	T● I F-			
	10-100 µl				•	T●IF-			
	50-200 μI				•	•	•		
	100-1000 μl				_	•	•	•	•
Research®	0.5-10 μl	•	•					- J	,
nescarell	2-20 µl			•	•	T● I F-			
	10-100 μΙ			•	•	T • I F-			
	20-200 µI				•		•		
				-	-	•	•	•	
Doggarch® =!··-	100-1000 µl	•							•
Research® plus	0.5-10 µl	•	•	_		Tele			
	10-100 µI			•	•	T●IF-			
	20-200 µl				•	•	•	_	_
	100-1000 μΙ							•	•

		PIPETTE TIPS (STANDARD / LOW RETENTION)								
Volume [µl]		10	10 extended	20	100	200	300	1000	1250	
Volume range [µI]		0.2-10	0.5-10	2-20	2-100	5-200	10-300	50-1000	50-1250	
Tip version		Tip/ Filter tip	Tip/ Filter tip	Filtertip	Filtertip	Tip/ Filter tip	Tip/ Filter tip	Tip	Tip/ Filter tip	
Article no.		7713XX	7723XX	7733XX	7743XX	7753XX	7763XX	7773XX	7783XX	
PIPETTE MODELS										
Research® pro	0.5-10 μΙ	•	•							
	5-100 µl				•	T● IF-				
	20-300 μΙ						•			
	·		Fixed-	volume pipett	es					
Reference®	50 μl				•	T● I F-				
	100 µl				•	T● I F-				
	500 µl							•	•	
	1000 μΙ							•	•	
Research®	500 μl							•	•	
EPPENDORF (cont	inued)									
			Multi-	channel pipett	es	1	ı	ı	ı	
Eppendorf	8CH 10-200 μI			•	•	•	•			
Research®	12CH 0.5-10 µI	•	•							
Research® plus	8CH 0.5-10 μI	•	•							
	8CH 10-100 μI			•	•	•	•			
	8CH 20-300 μI			•	•	•	•			
			Elec	tronic pipettes	3					
Xplorer [®]	50-1000 μΙ							•	•	
	8CH 15-300 μI			•	•	•	•			
GILSON										
OLEOGIA .			Single-	channel pipet	tes					
Pipetman [®]	P2 (0.5-2 µI)	•	• Onigic	onamici piper						
ripetinan	P10 (1-10 μl)	•	•							
	P20 (2-20 µl)	•	•	•		T● I F-				
	P100 (20-100 µI)			•	•	•				
	P200 (50-200 µI)				•	•	•			
					-	•	•	•	•	
Dia -+	P1000 (200-1000 μl)							•	•	
Pipetman® L	P2L (0.2-2 μl)	•								
	P10L (1-10 µl)	•	•			T = 1 =				
	P20L (2-20 µl)			•	_	T● I F-				
	P200L (20-200 µl)				•	•	•	_		
	P1000L (100-1000 μl)							•	•	
			Fixed-	volume pipett	es					
Pipetman®	F5 (5 µI)			•		T●IF-				
	F10 (10 µI)			•		T● I F-				
	F20 (20 µI)			•		T●IF-				
	F25 (25 µI)				•	T● I F-				

T = Pipette tip

• compatible - not compatible

F= Filter tip

 \blacksquare Pipetting volume limited to the maximum volume of the pipette tips

				PIPETT	TIPS/STAND	ARD / LOW RET	ENTION)		
Volume [µI]		10	10 extended	20	100	200	300	1000	1250
Volume range [µl]		0.2-10	0.5-10	2-20	2-100	5-200	10-300	50-1000	50-1250
		Tip/	Tip/			Tip/	Tip/	Tip	Tip/
Tip version		Filtertip	Filtertip	Filtertip	Filtertip	Filtertip	Filtertip		Filterti
Article no.		7713XX	7723XX	7733XX	7743XX	7753XX	7763XX	7773XX	7783X)
PIPETTE MODELS									
GILSON (continued)									
			Fixed	-volume pipett	es				
Pipetman®	F100 (100 µI)						T●IF-		
	F200 (200 µI)					•	•		
	F250 (250 µI)							•	
	F300 (300 µI)							•	•
	F500 (500 µI)							•	
	F1000 (1000 µl)							•	•
				channel pipett	es				
Pipetman [®] L	P8X 10L (0.5-10 μl)	•	•						
	P8X 200L (20-200 μI)			•	•	•	•		
	P8X 300L (20-300 μI)					•	•		
IKA									
			Single	-channel pipet	tes		ı	I	
IKA Pette vario	100-1000 μΙ							•	
METTLER TOLEDO									
			Single	-channel pipet	tes				
Rainin Pipette-Lite™	SL200 (20-200 µI)			-	•	•	•		
Rainin Pipette-Lite™	SL1000 (100-1000 µI)							•	•
NICHIRYO									
Nichikito			Single	-channel pipet	tes				
Nichipet	0.5-10 μl	•	• Omgic	channer piper					
Monipet	2-20 μΙ	•				T● I F-			
	10-100 µl			•	•	•	•		
	20-200 μl			-	•	•	•		
				-	-			•	
	100-1000 μΙ		Multi	channel pipett	•				
Nichipet	8K 30-300 μI		riuiti-	channer pipett	es ■	•	•		
	ок 30-300 µ				_	_			
SARTORIUS									
				-channel pipet	tes				
Biohit Proline®	0.5-10 μΙ	•	•						
Biohit Proline® plus	0.5-10 μΙ	•	•						
	2-20 μΙ			•		T●IF-			
	10-100 μΙ			•	•	•	•		
	20-200 μΙ			•	•	•	•		
	100-1000 μΙ							•	•
				channel pipett	es				
Biohit Proline®	8CH 0.5-10 μI	•	•						
	8CH 25-250 μI					•	•		
	8CH 50-300 μI						•		
SOCOREX									
			Single	-channel pipet	tes				
Acura® 825	0.5-10 μΙ	•	•						

				PIPETTI	E TIPS (STAND	ARD / LOW RET	ENTION)		
Volume [µI]		10	10 extended	20	100	200	300	1000	1250
Volume range [µI]		0.2-10	0.5-10	2-20	2-100	5-200	10-300	50-1000	50-1250
Tip version		Tip/	Tip/			Tip/	Tip/	Tip	Tip/
		Filter tip	Filtertip	Filtertip	Filtertip	Filter tip	Filtertip		Filtertip
Article no. PIPETTE MODELS		7713XX	7723XX	7733XX	7743XX	7753XX	7763XX	7773XX	7783XX
PIPET TE MODELS	2-20 µI			•		T● I F-			
	5-50 μl				•				
				•	•	T● I F-			
	10-100 μl 20-200 μl			-	•	T●IF-	•		
	100-1000 µI			-	-		<u> </u>	•	
Calibra® 822			•						•
Calibra® 822	1-10 µl		•	_					
	2-20 μΙ			•		T●IF-			
	10-100 μΙ			•	_	T●IF-			
	20-200 μΙ			•	•	•	•		
	100-1000 μl							•	•
			Multi-	channel pipett					
Acura® 855	8CH 5-50μI			•	•	T●IF-			
Calibra® 852	8CH 1-10 μI	•	•						
	8CH 20-200 µI			-	•	T● I F-			
	12CH 10-100 μI					T● I F-			
STARLAB									
			Single	-channel pipet	tes				
ErgoOne®	0.1-2.5 µl		•						
	0.5-10 μΙ	•	•						
	10-100 µl			•	•	T● I F-			
	100-1000 µl							•	•
			Multi-	channel pipett	es				
ErgoOne®	8 CH 30-300 µI				•		•		
		l							
THERMO FISHER SCIE	INTIFIC								
			Single	-channel pipet	tes				
Finnpipette [™] F1	1-10 μΙ	•	•						
	2-20 μΙ			•		T● IF-	•		
	20-200 μΙ			-	•	•	•		
	30-300 μl						•		
	100-1000 μΙ							•	•
Finnpipette [™] 4500	200-1000 μΙ							•	•
			Multi-	channel pipett	es				
Finnpipette [™] F1 MCP8	8 CH 1-10 μI	•	•						
	8 CH 10-100 μI			•	•	•	•		
	8CH 30-300 μI				•	-	•		
Finnpipette [™] 4510	8CH 50-300 μI				•	•	•		
Finnpipette™ 4500	8CH 200-1000 μI							•	•
VWR									
			Single	-channel pipet	tes				
Erg. High Perform.	0.5-10 µl	•	•	- I a i i i i i i i i i i i i i i i i i i					
Erg. High Perform.	2-20 µI	-	-	•		T● I F-			
Erg. High Perform.	10-100 μΙ			•	•	•			
Erg. High Perform.	20-200 μI			-	•	•	•		
Erg. High Perform.	100-1000 µI			-	_	_	_	•	•
Erg. nigir erroriii.	100-1000 μΙ							•	•

LABORATORY INFORMATION FOR PCR

OVERVIEW PCR MICROPLATES

The use of the 96 well format allows the scale up of basic PCR work, while the 384 well format is ideal for high-throughput screening projects. All microplates are made of thin-walled polypropylene. This optimises the heat transfer from the thermoblock to

the reaction solution. Our heat-resistant sealers AMPLIseal, VIEWseal and SILVERseal are ideal for sealing the microplates during PCR, and the 96 well microplate may also be sealed with 8-cap strips.

96 WELL POLYPROPYLENE MICROPLATES FOR PCR

1/ NON-SKIRTED MICROPLATES

Non-skirted microplates may be used in all commonly available thermocyclers with a 96 well block.

1a) Non-skirted microplate with raised well rims (Art. No. 652201)



-CHARADADADA A A A A A A

Figure 1a: View of a non-skirted microplate with raised well rims

1b) Non-skirted microplate with flat surface (Art. No. 652250)

Black alphanumeric coding enables a quick identification of samples



#HAAAAAAAAAAA

Figure 1b: View of a non-skirted microplate with flat surface

1c) Non-skirted microplate, low profile (Art. No. 652210)



AAAAAAAAAAAA

Figure 1c: View of a non-skirted low profile microplate

2/ HALF-SKIRTED MICROPLATES

2a) Half-skirted microplate with one notch suitable for real-time PCR systems such as LightCycler® 480 (Art. No. 669285)

- / Maximal pigmented white polypropylene and therefore most suitable for sensitive real-time PCR reactions
- / Black alphanumeric coding enables a quick identification of samples
- Notches in the rim facilitate automation due to better gripping in robotic systems



-77777777

Figure 2a: View of a half-skirted microplate with one notch suitable for real-time PCR systems such as LightCycler® 480 from Roche

 $The \ registered \ trademarks \ of \ the \ mentioned \ manufacturers \ belong \ to \ the \ above \ mentioned \ companies.$

2b) Half-skirted microplate with one notch suitable for ABI (Art. No. 652290)





Figure 2b: View of a half-skirted microplate with one notch suitable for ABI

2c) Half-skirted microplate, recessed rim, ABI design with one notch (Art. No. 652260)





Figure 2c: View of a half-skirted microplate, recessed rim, ABI design with one notch

3/ FULL-SKIRTED MICROPLATE WITH ONE NOTCH (ART. NO. 652270)





Figure 3: View of full-skirted microplate with one notch

384 WELL POLYPROPYLENE MICROPLATES FOR PCR

The 384 well PCR microplates from Greiner Bio-One are manufactured in an advanced injection moulding process following stringent quality criteria. Minimal distortion and sagging curvature, homogeneous heat transfer and sealing of the individual wells are essential quality criteria here. The footprint of all 384 well PCR microplates is compatible with automated systems.

1/ FULL-SKIRTED 384 WELL MICROPLATE WITH ONE NOTCH SUITABLE FOR ABI (ART. NO. 785290)





Figure 4: Full-skirted 384 well microplate with one notch and alphanumeric coding suitable for ABI

2/ FULL-SKIRTED 384 WELL MICROPLATE WITH TWO NOTCHES (ART. NO. 785201)





Figure 5: Full-skirted 384 well microplate with two notches and alphanumeric coding

3/ FULL-SKIRTED 384 WELL MICRO-PLATE WITH TWO NOTCHES FOR REAL-TIME PCR SYSTEMS (ART. NO. 785285)

- / White pigmentation boosts real-time PCR signal
- Black alphanumeric coding enables a quick identification of samples





Figure 6: Full-skirted 384 well microplate with two notches for real-time PCR systems such as LightCycler® 480 from Roche

 $The \ registered \ trademarks \ of \ the \ mentioned \ manufacturers \ belong \ to \ the \ above \ mentioned \ companies.$

COMPATIBILITY FOR PCR MICROPLATES

	▲ Optimal Fit	Model	652201 unskirted, chimneytop	652250 unskirted, universal	652210 unskirted, low-profile	652260 semi-skirted, ABI-Design	652270 skirted	652290 semi-skirted, suitable for ABI	669285 semi-skirted, RT-PCR	785201 skirted	785285 skirted, RT-PCR	785290 skirted, suitable for ABI
						96 well					384 well	
		2700	•	•		•		•				•
		2720				A						•
	Thermal	6100				A						
	Cyclers	9600	•	•		•		•				
		9700	•	•		A		•				A
		Veriti 0.2 ml				A		A				•
		5700		•		•		•				
JIS		PRISM 7000		•		A		A				
ster	Real Time	7300		•		A		A				
iosy	Thermal	7500		•		A		A				
d B	Cyclers	7700		•		•		A				
Applied Biosystems	0,0.0.0	7900HT		•		A						•
A		ViiA 7				A						•
		Quant Studio 12K Flex				•						•
		PRISM 310	•	•		•		A				
		PRISM 3100	•	•		A		•				A
	Sequencers	3130(XL)	•	•		A		A				A
	Sequencers	3700	•	•		A		A				A
		PRISM 3730 (XL)	•	•		A		A				A
		3500(XL)				•						•
		Gene Cycler	•	•		•						
		PTC-100	•	•	•		A	•				•
		PTC-200	•	•	•		A	•				•
		PTC-225 Tetrad	•	•	•		A	•				•
		Dyad/Dyad Disciple	•	•	•		A	•				•
	Thermal	iCycler	•	•			•	A				
	Cyclers	MyCycler	•	•								
c		Mini Gradient		•	•							
Bio-Rad/MJResearch		Personal		•				•				
Res		T100		A			A					
Σ		DNA Engine family		•	•		•	•		•		•
ad/		C1000/S1000		•	A		•	•		•		•
-0-F		Opticon/Opticon2		•	•		A					
Δ.		Chromo-4			•		•					
	Real Time	iCycler	•	•			•	A				
	Thermal	MyiQ	٠	•			•	A				
	Cyclers	iQ5	•	•			•	A				
		CFX Connect			•		•	•				
		CFX96					A					
		CFX384								٠		•
	Sequencers	BaseStation					•					
		Mastercycler	•	•	•		•	•				
		Mastercycler ep Gradient	•	•		•	A	•				
orf	Thermal	Mastercycler M384								٠		٠
Eppendorf	Cyclers	Mastercycler Nexus		•	•		•	•				
pdd:		Mastercycler Nexus Gradient		•	•		A	•				
ш		Mastercycler Nexus Eco		•	•		•	•				
		Mastercycler Pro		•			•	•		٠		٠
	Real Time T.C.	Mastercycler ep Realplex	•	•			A	•				

The registered trademarks of the mentioned manufacturers belong to the above mentioned companies. This compatibility chart is a general guide only and subject to error and modifications. We cannot accept any liability or responsibility for the above information.

	▲ Optimal Fit	Model	652201 unskirted, chimneytop	652250 unskirted, universal	652210 unskirted, low-profile	652260 semi-skirted, ABI-Design	652270 skirted	652290 semi-skirted, suitable for ABI	669285 semi-skirted, RT-PCR	785201 skirted	785285 skirted, RT-PCR	785290 skirted, suitable for ABI
						96 well					384 well	
Amers- ham	Sequencers	MegaBACE 500 MegaBACE 1000 MegaBACE 4000					•					•
Beck- mann	Sequencers	CEQ		•								
Biometra	Thermal Cyclers	Uno Uno II T1 Thermal Cycler Tgradient Trobot T3000	•	•	•		•	•				•
		T Professional		•	•		•					•
Cor- bett	Real Time T.C. Thermal Cyclers	Toptical PalmCycler 96		•	•		•	•				•
Ericom	Thermal Cyclers	PalmCycler 384 Power Block I Deltacycler I Deltacycler II Single Block Twin Block	•	•	•	•		•				•
Esco	Thermal Cyclers	Swift Gene Genius	•	•			•	•				•
G-Storm	Thermal Cyclers	GS1 GS2 GS4 GSX	•	•	•	•		•				
MWG	Thermal Cyclers	Primus 96 Primus 384	٠	•			•					•
Strate- gene	Thermal Cyclers Real Time T.C.	Robocycler 96 Robocycler Gradient Mx4000 and Mx3005P	•	•	•	•	•	•				
ТаКа- Ra	Thermal Cyclers	TP240 TP3000					•					
Techne	Thermal Cyclers	Touchgene Cyclogene Genius Genius Quad Genius(TC412)	•	•	•	•	•	•				
Te	Paul Tima T C	Flexigene Touchgene X Touchgene Gradient (TC512)	•	•	•	•	•	•				•
Thermo Hybaid	Thermal Cyclers	Quantica PCR Sprint MBS Satelite (Multiblock) Px2 and PxE PCR Express and Omni-E Touchdown Omnigene	•	•	•	•	•	•				•
Trans- gen- omic	Sequencers	WAVE System					•					
Ro- che	Real Time T.C.	LightCycler 96 Light Cycler 480							•	•		•

 $The \ registered\ trademarks\ of\ the\ mentioned\ manufacturers\ belong\ to\ the\ above\ mentioned\ companies.\ This\ compatibility\ chart\ is\ a\ general\ guide$ only and subject to error and modifications. We cannot accept any liability or responsibility for the above information.

COMPATIBILITY FOR MINI BLOCK HEATER INSERTS

Art. no.	848916	848923	848902	848913	848921	848915	848950
No. of Tubes per Insert	15x	15x	40x	15x	15x	4x	2x
1ml Cryo.s					•		
2 ml Cryo.s					•		
4 ml Cryo.s				•			
5 ml Cryo.s				•			
1.5 ml Reaction Tube	•	•					
2 ml Reaction Tube		•					
0.5 ml PCR Tube	•	•					
0.2 ml PCR Tube			•				
1x8 PCR Tube Strip			•				
15 ml Conical Tube						•	
50 ml Conical Tube							•
4 ml, 12 x 55 mm Tube					•		
5 ml, 12 x 75 mm Tube				•			
4.5 ml, 12 x 75 mm Tube				•			
7 ml, 13 x 100 mm Tube				•			
12 ml, 16 x 100 mm Tube						•	
12 ml, 17 x 100 mm Tube						•	
20 ml, 16 x 152 mm Tube						•	
14 ml, 17 x 95 mm Tube						•	
14 ml, 18 x 95 mm Tube						•	

This compatibility chart is a general guide only and subject to error and modifications. We cannot accept any liability or responsibility for the above information.

LABORATORY INFORMATION FOR CENTRIFUGATION

CENTRIFUGATION - PRINCIPLE AND CALCULATION OF THE RCF (RELATIVE CENTRIFUGAL FORCE)

Sedimentation of particles in a gravitational field

If a mixture of sand and water is shaken thoroughly and then left to stand, the sedimentation of the solid particles takes place according to their size. As a result of gravitational acceleration $(g = 9.81 \text{m/s}^2)$, all of the particles are located in a gravitational field under the

influence of which the coarse grains of sand collect at the bottom first and the smaller grains of sand are deposited later. After around 10-20 minutes, the following layering is produced (from bottom to top): coarse grains of sand - fine grains of sand - water. However, other particles (proteins, nucleic acids, viruses, pro- or eucaryotic cells) do not necessarily

precipitate or only sediment out after they have been exposed to higher forces than the force of gravity resulting from the gravitational acceleration. If these forces exceed the counter-forces resulting from convection (heat circulation) and Brownian molecular motion, both of which cause constant mixing of solutions and suspensions, sedimentation takes place.

The sedimentation rate can be calculated on the basis of Stoke's law as follows:

$$V = \frac{d^2(p_p - p_L)g}{18\mu}$$

However, a particle will only sediment out if $p_P > p_L$. If $p_P < p_L$, V becomes negative, consequently the particle floats rather than sedimenting out.

/ V = sedimentation rate

 p_{P} = density of the particle

 p_i = density of the liquid

 $g = 9.81 \,\mathrm{m/s^2}$

 μ = viscosity of the liquid

Influence of the Centrifugation and Calculation of the RCF respectively RPM

A centrifuge can be used to create a transient gravitational field under the influence of which the sedimentation of cells, cellular components and macro-molecules takes place. In a centrifuge, a suspension located in a centrifuge tube

rotates around a rotational axis. Each particle of the suspension is subject to centrifugal force, which moves it radially away from the rotational axis.

The centrifugal force F_c is calculated as follows:

$$F_{c} = m_{p} w^{2} r$$

/ where m_p = mass of the particle, w= angular velocity (s⁻¹) and r = distance of the particle from the rotational axis

The force acting on a particle in a centrifugal field is stated relative to gravitational acceleration, usually as so-called relative centrifugal force (RCF) or g-force (x g). It is calculated as follows:


$$RCF = 11.18r \left(\frac{R}{1000} \right)^2$$

/ where R = rotor revolutions per min and r = distance of the particle from the rotational axis (cm)

For easier conversion of the *RCF* or *g*-force into revolutions per min, the equation can be transformed as follows:

$$R = 299 \sqrt{\frac{RCF}{r}}$$

ALIGNMENT CHART

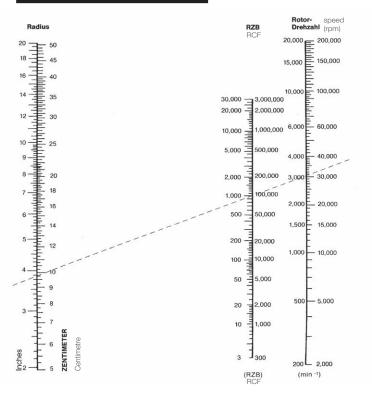

/ By use of a ruler, the third value relating to two known scale values can be read from the alignment chart.

MAXIMAL CENTRIFUGE CAPACITY OF TUBES, REACTION TUBES AND MICROPLATES

The maximum centrifuge capacity for Greiner Bio-One tubes, reaction tubes and microplates is listed in the form of the RCF in the tables below.

For centrifugation, all products were filled with water up to their maximum filling volume. Determination of the maximum RCF in a swinging-bucket rotor was conducted in a Thermo Scientific Centrifuge (Heraeus Multifuge BSR

Plus). Determination of the maximum RCF in a fixed angle rotor was conducted in a Sorvall Centrifuge (Evolution RC). Therefore special rotor inserts for different vessel shapes and sizes were used for a stable fit.

REACTION TUBES

Art. no.	Volume[ml]	max. RCF [g] fixed-angle rotor
6162XX	1.5	18000
6232XX	2.0	16000
6672XX	0.5	15000
6932XX	0.5	18000
7162XX	1.5	20000
7172XX	1.5	20000
7222XX	2.0	22000

The stated maximum RCF values are guidelines only, depending on a variety of factors such as rotor, temperature, density, pH and type of liquid. The suitability of tubes for a specific application using high centrifugation forces has to be tested.

POLYSTYRENE TUBES

Art. no.	Dimensions ø [mm] x height [mm]	max. RCF[g]swinging-bucket rotor	max. RCF[g]fixed-angle rotor
1031XX	10.5×40	5800	_1
1121XX	12×55	5800	6200
1151XX	12×75	5800	5800
1161XX	12×75	5800	_1
1201XX	12.4×75	4800	5000
1251XX	13×100	4000	7500
1361XX	14 x 100	4000	5200
1601XX	16 x 100	3500	5000
1631XX	17×100	3000	3000
1641XX	16.8×100	5000	5000
1661XX	16×110	2500	3200
1861XX	17 x 120	2500	2800
1871XX	17×100	5200	6600
1881XX	17 x 120	2500	4500
1911XX	18×95	4000	5500
2011XX	24×90	1000	3500

¹ No fitting rotor inserts available.

POLYPROPYLENE TUBES

Art. no.	Dimensions ø [mm] x height [mm]	max. RCF[g]swinging-bucket rotor	max. RCF[g]fixed-angle rotor
1022XX	8.5 x 44	5800	20000
1122XX	12×55	5800	30000
1152XX	12×75	5800	34000
1212XX	12.5×48	5800	34000
1222XX	12.5×48	5800	34000
1232XX	12.5 x 42	5800	26000
1242XX	12.5×86	5800	34000
1262XX	12.4×47	5800	26000
1272XX	12.4×83	5800	34000
1602XX	16×100	5800	33000
160297	16×100	3500	33000
1632XX	16×100	5000	26000
184261	17×77	4800	34000
187201	17×100	4800	34000
187261	18×95	4800	34000
1882XX	17×120	4000	15000
1912XX	18×95	4800	34000
2102XX	30×115	2800	9500
2272XX	30×115	3200	9500
227261	30×115	3200	17000
227270	30×115	3200	17000
227281/227285	30×115	3200	9500
227280/227283	30 x 115	3200	9000

 $The stated \, maximum \, RCF \, values \, are \, guidelines \, only, \, depending \, on \, a \, variety \, of \, factors \, such \, as \, rotor, \, temperature, \, density, \, pH \, and \, type \, of \, liquid.$ $The \ suitability \ of \ tubes \ for \ a \ specific \ application \ using \ high \ centrifugation \ forces \ has \ to \ be \ tested.$

POLYETHYLENE TUBES

Art. no.	Dimensions ø [mm] x height [mm]	max. RCF[g]swinging-bucket rotor	max. RCF[g]fixed-angle rotor
1123XX	12×55	4200	22000
1153XX	12×75	4200	20000
1603XX	16×100	3500	30000
1873XX	17×100	5800	20000

MULTIWELL PLATES

Art. no.	Multiwell Plate	max. RCF[g] swinging-bucket rotor
657160	6 well, PS, clear	4800
665102	12 well, PS, clear	4800
662160	24 well, PS, clear	4800
677180	48 well, PS, clear	4800

MICROPLATES

Art. no.	Microplate	max. RCF[g]swinging-bucket rotor
650101	96 well, PS, U-bottom, clear	1000
651101	96 well, PS, V-bottom, clear	1000
655101	96 well, PS, F-bottom, clear	4800
650201	96 well, PP, U-bottom, natural	4800
651201	96 well, PP, V-bottom, natural	4800
655201	96 well, PP, F-bottom, natural	4800
655209	96 well, PP, U-bottom, black	4800
655074	96 well, PS, F-bottom, white	4800
655076	96 well, PS, F-bottom, black	4800
655094	96 well, PS, µClear®, white	4800
655096	96 well, PS, µClear®, black	4800
655801	96 well, PS, UV-Star®	4800
780201	96 well, PP, MASTERBLOCK® 1 ml	4800
780270	96 well, PP, MASTERBLOCK® 2 ml	4800
786201	96 well, PP, MASTERBLOCK® 0.5 ml	4800
781101	384 well, PS, clear	4800
781073	384 well, PS, white	4800
781077	384 well, PS, black	4800
781094	384 well, PS, μClear®, white	4000
781096	384 well, PS, µClear®, black	3000
781201	384 well, PP, F-bottom, natural	4800
781280	384 well, PP, V-bottom, natural	4800
781270	384 well, PP, V-bottom, Deep Well, natural	4800
781801	384 well, PS, UV-Star®	4800

 $The stated \ maximum \ RCF \ values \ in \ these \ tables \ are \ guidelines \ only, \ depending \ on \ a \ variety \ of factors \ such \ as \ rotor, \ temperature, \ density, \ pH \ and \ type$ $of \ liquid. \ The \ suitability \ of \ tubes \ and \ plates \ for \ a \ specific \ application \ using \ high \ centrifugation forces \ has \ to \ be \ tested.$

MICROPLATES

Art. no.	Microplate	max. RCF [g] swinging-bucket rotor
784101	384 well, PS, Small Volume, clear	800
784075	384 well, PS, Small Volume, white	800
784076	384 well, PS, Small Volume, black	800
784201	384 well, PP, Small Volume, natural	4800
782101	1536 well, PS, HiBase, clear	1800
782074	1536 well, PS, HiBase, white	1500
782077	1536 well, PS, HiBase, black	1500
782094	1536 well, PS, µClear®, HiBase, white	1000
782096	1536 well, PS, µClear®, HiBase, black	1500
782270	1536 well, PP, V-bottom, Deep Well, natural	4800
783101	1536 well, PS, LoBase, clear	4800
783075	1536 well, PS, LoBase, white	4800
783076	1536 well, PS, LoBase, black	4800
783094	1536 well, PS, µClear®, LoBase, white	4800
783096	1536 well, PS, µClear®, LoBase, black	4800

PCR PLATES

Art. no.	PCR Plate	max. RCF[g] swinging-bucket rotor
652270	96 well, PP, natural, full-skirt	4800
652290	96 well, PP, natural, half-skirt, suitable for ABI	4800
785201	384 well, PP, natural, full-skirt	4800
785290	384 well, PP, natural, full-skirt, suitable for ABI	4800

For centrifugation the plates were filled with water as follows: $96 \text{ well} = 300 \, \mu\text{I} + 384 \, \text{well} = 50 \, \mu\text{I} + 1536 \, \text{well} = 5 \, \mu\text{I} + 160 \, \text{m}$

The stated maximum RCF values in these tables are guidelines only, depending on a variety of factors such as rotor, temperature, density, pH and type of liquid. The suitability of tubes and plates for a specific application using high centrifugation forces has to be tested.

LABORATORY INFORMATION FOR CRYO.S SAMPLE STORAGE

FREEZING PROTOCOL

0

Wash the cells with warm PBS solution, aspirate the solution and cover the cells with a solution containing trypsin and EDTA (a thin liquid film is enough; the concentration should be evaluated for each cell line).

2

Incubate the cells for max. 3-5 min at 37° C.

3

Once the cells detach from the bottom, stop incubation by adding cell culture medium supplemented with serum and slightly suspend cells using a pipette.

4

Spin down the suspension (500 x g, 5 min) and resuspend the pellet with medium containing serum.

5

Determine the cell number (using a Neubauer chamber).

6

Spin down the cells for 5 min at 500 x g and discard the supernatant. Resuspend the pellet with an adequate volume of cell culture medium containing serum.

7

Mix the cell suspension 1:1 with freezing medium (60 % medium, 20 % FCS, 20 % DMSO) and transfer it in Cryo.s. For freezing in Cryo.s, the concentration of cells should be $1-5 \times 10^6$ cells / ml.

Cryo.s containing cells should

8

be frozen at a cooling rate of -1 K/min. This can be achieved by placing them into an isopropanol-filled chamber at -70 °C. If other types of samples are contained, Cryo.s may be frozen directly at -20 °C, -70 °C or in the gas phase of liquid nitrogen. In order to assure even freezing of the sample, 4 and 5 ml Cryo.s should be frozen at -20 °C overnight before transferring them to -70 °C or to the gas phase of liquid nitrogen.

9

Then transfer the Cryo.s into the nitrogen tank. To avoid contamination (e.g. mycoplasma) and due to safety precautions Cryo.s must only be stored in the gas phase above and not in the liquid nitrogen.

THAWING PROTOCOL



Immediately after removing them out of the nitrogen tank the frozen cells are thawed in about 1-2 min brandishing the Cryo.s in a water bath at 37°C. The thawing process should be performed as fast as possible.



Transfer the thawed cell suspension into a 15 ml tube and mix it immediately with copious amounts of cell culture medium containing serum.



After spinning down the cells (500 x g, 5 min) discard the supernatant and resuspend the pellet in an appropriate cell culture medium supplemented with serum and transfer it into one or more cell culture flasks.



Follow the recommended cell concentration for seeding.



During the next 12 hours cells should rest.



A change of medium is recommended after 24 resp. 48 hours.

SAFETY ADVISORY FOR **WORKING WITH CRYO.S**

Cryo.s, tubes are intended for sample storage exclusively in the gas phase over liquid nitrogen or in freezers! If Cryo.s are stored in the liquid phase, nitrogen can seep into the tubes. Then upon thawing the vaporising nitrogen can generate high pressure, ultimately resulting in an explosion, as well as the release of any infectious material.

Always take appropriate personal safety measures when working with Cryo.s, including wearing safety clothing, using goggles and working at a safety laboratory bench. When undertaking cryogenic preservation, Cryo.s must be evenly exposed to freezing temperatures. Uneven temperature exposures can cause formation of ice plugs (i.e. at tube top) that inhibit the expansion of freezing liquid (i.e. at tube bottom), resulting in dangerous high pressure and subsequent harm or damage of tubes.

Never exceed maximum working volumes as specified.

LABORATORY INFORMATION FOR IMMUNOLOGY

VOLUME-DEPENDENT WETTING OF IMMUNOLOGICAL PRODUCTS

Liquid volume [µl]	Covered area	Liquid height	Area / volume	Liquid volume [µl]	Covered area	Liquid height	Area / volume
		roplate, U-Bot			LISA Micropla		
25		•			•		
	34	1.7	13.6	25	38	1.65	15.2
50	52	2.6	10.4	50	60	3.2	12.0
75	68	3.4	9.1	75	81.5	4.7	10.9
100	84.6	4.2	8.5	100	103.6	6.2	10.4
125	99	4.9	7.9	125	124.5	7.6	10.0
150	115.5	5.7	7.7	150	144	8.9	9.6
175	130	6.4	7.4	175	165.8	10.3	9.5
200	145	7.1	7.3	200	181.7	11.5	9.1
225	160	7.8	7.1		C8 Stri	p Plate	
250	174.7	8.5	7.0	25	39	1.0	15.6
275	190	9.2	6.9	50	56	1.9	11.2
300	205	9.9	6.8	75	73	2.8	9.7
96 V	Vell ELISA Mic	roplate, V-Bot	tom	100	88.6	3.6	8.9
25	35	2.3	14.0	125	104.3	4.4	8.3
50	54.6	3.4	10.9	150	120	5.2	8.0
75	72.4	4.4	9.7	175	136.5	6.0	7.8
100	88.6	5.3	8.9	200	150.8	6.7	7.5
125	105	6.2	8.4	225	165.4	7.4	7.4
150	123.8	7.2	8.3	250	181	8.15	7.2
175	140.8	8.1	8.0	275	196	8.85	7.1
200	156	8.9	7.8	300	211	9.55	7.0
		te, F-Bottom/			F8 Stri		7.0
25	47	0.8	18.8	25	50.4	0.8	20.2
50	62		12.4	50	64		12.8
75	77.5	1.55 2.3	12.4	75	79.7	1.45 2.2	10.6
100	92	3.0	9.2	100	93.5	2.2	9.4
		3.8			108.3		8.7
125	108		8.6	125		3.55	
150	123	4.5	8.2	150	123	4.25	8.2
175	137.6	5.2	7.9	175	138	4.95	7.9
200	152.3	5.9	7.6	200	153	5.65	7.7
225	168	6.65	7.5	225	167	6.3	7.4
250	183	7.35	7.3	250	182	7.0	7.3
275	197	8.0	7.2	275	196	7.65	7.1
300	212	8.7	7.1	300	211.5	8.35	7.1
96 Well ELI	SA Microplate,	F-Bottom/Ch	imney Well		U8 Stri	p Plate	
25	47	0.7	18.8	25	34	1.7	13.6
50	64	1.5	12.8	50	52	2.6	10.4
75	78.5	2.2	10.5	75	68	3.4	9.1
100	93	2.9	9.3	100	84	4.2	8.4
125	108	3.6	8.6	125	99.6	4.95	8.0
150	122.6	4.3	8.2	150	115	5.75	7.7
175	137.5	5.0	7.9	175	129.6	6.45	7.4
200	152	5.7	7.6	200	144	7.15	7.2
225	167	6.4	7.4	225	159	7.85	7.1
250	182	7.1	7.3	250	174	8.55	7.0
275	197	7.8	7.2	275	189	9.25	6.9
300	212	8.4	7.1	300	204	9.95	6.8
000	- 12	0.1	7.1	000	201	0.00	0.0

Liquid volume [µl]	Covered area [mm²]	Liquid height [mm]	Area / volume ratio [cm² / cm³]				
	U16 Strip Plate						
25	35	1.75	14.0				
50	52	2.6	10.4				
75	68	3.4	9.1				
100	84	4.2	8.4				
125	98.6	4.9	7.9				
150	115	5.7	7.7				
175	129.6	6.4	7.4				
200	144	7.1	7.2				
225	159	7.8	7.1				
250	174	8.5	7.0				
275	189	9.2	6.9				
300	204	9.9	6.8				

Liquid volume [µl]	Covered area [mm²]	Liquid height [mm]	Area / volume ratio [cm²/cm³]
	F16 Stri	p Plate	
25	49	0.8	19.6
50	63	1.5	12.6
75	79.8	2.3	10.6
100	94.3	3.0	9.4
125	108	3.7	8.6
150	123.5	4.4	8.2
175	138	5.1	7.9
200	153	5.8	7.7
225	168	6.5	7.5
250	183	7.2	7.3
275	198	7.9	7.2
300	213	8.6	7.1
3	84 Well Microp	late, F-Botton	n
25	39.07	2.50	15.6
50	66.60	4.8	13.3
75	94.03	7.00	12.5
100	119.63	9.05	12.0
125	145.6	11.05	11.6
132	152.6	11.50	11.6

Abbreviations

ANSI	American National Standards Institute	PLL	Poly-L-Lysine
COC	Cycloolefin co-polymer	PP	Polypropylene
COP	Cycloolefin polymer	PS	Polystyrene
CV	Coefficient of Variation	PTFE	Polytetrafluoroethylene
DMS0	Dimethyl Sulphoxide	RNA	Ribonucleic Acid
DNA	Deoxyribonucleic Acid	RNase	Ribonuclease
DNase	Deoxyribonuclease	rRNA	Ribosomal RNA
ECM	Extracellular Matrix	RT	Room Temperature
EL-Rack	EasyLoad® Rack	SBS	Society for Biomolecular Sciences
ELISA	Enzyme Linked Immuno Sorbent Assay	SPA	Scintillation Proximity Assays
EVA	Ethyl Vinyl Acetate	ST-Rack	Standard Rack
FDA	Food and Drug Administration	TC	Tissue Culture
FIA	Fluorescence Immuno Assay	USP	United States Pharmacopoeia
F-Rack	Filter Tip Rack	UV Spectrum	Ultraviolet Spectrum
HDPE	High Density Polyethylene	VIS Spectrum	Visible Spectrum
HLA	Human Leucocyte Antigen		
HTS	High-Throughput Screening	Units	
IgG	Immunoglobulin G	°C	Degree Centigrade
ID-Card	Identity Card	Da	Dalton, the unit of molecular mass
LAL	Limulus Amoebocyte Lysate	g	Gram or Gravitational
LIA	Luminescence Immuno Assay		Acceleration (about 9.81 m/s²)
med.	Medium	Gy	Gray, Radiation Unit
NMWCO	Nominal Molecular Weight Cut-Offs	h	Hour
PC	Polycarbonate	1	Liter
PCR	Polymerase Chain Reaction	M	Molarity, moles of solute
PDL	Poly-D-Lysine		per litre of solution
PET	Polyethylene Terephthalate	m	Meter
PETG	Polyethylene Terephthalate Copolymer	min	Minute
рН	pH Value	Mol	Absolute Amount of Substance
PLA	Polylactate	S	Second

GLOSSARY

Advanced TC is a polymer modification increasing the cellular primary and long-term adhesion of Greiner Bio-One cell culture vessels. Based on the innovative technique the surface of the cell culture vessels is modified to positively influence cellular features and functions. Enhanced cellular adhesion and higher proliferation rates improve cell expansion and cultivation of sensitive cells or cells under restricted growth conditions.

Biobanking Tubes are $300\,\mu$ l, $600\,\mu$ l and $1000\,\mu$ l Cryo.s tubes for the efficient storage of biological samples in large-scale biorepositories. The design of tubes and racks allows for a very space-efficient storage with up to 30~% better utilisation of storage space in freezers or liquid nitrogen tanks. In addition, they are optimised for sample storage at extremely low temperatures over long periods of time.

Bioburden is used to describe the colonisation of viable microorganisms on a material or product and is the basis for determining the necessary radiation dose for sterilisation.

C-bottom stands for the well profile of a flat well bottom with rounded corners.

CELLCOAT® is the Greiner Bio-One brand name for all protein-coated cell culture vessels for adherent cell culture.

CELLMASTER is a quality term that refers to all roller hottles.

CELLreactor is a 15 ml/50 ml polypropylene tube with filter screw cap for the cultivation of suspension and spheroid cells, expansion of aerobic bacteria, yeast or other microorganisms as well as storage of components and liquids requiring gas exchange.

Cell-repellent surface reliably prevents cell attachment in suspension cultures of semi-adherent and adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.

CELLSTAR® is a Greiner Bio-One brand name and includes culture vessels with physically modified surfaces for adherent or suspension cell cultures.

CELLview is a quality term for cell culture products with glass bottom for high-resolution microscopic applications.

Datamatrix Code is a 2D barcode which can also be used for tracking biological and medical reagents and samples. Its small footprint provides nearly infinite scalability and large data capacity. Datamatrix codes can be scanned independent of their orientation and are very accurate due to the Reed-Solomon error correction method.

Deep Well microplates have conical bottom wells and are ideally suited for the storage of non-human samples.

EASYstrainer are cell strainers for the fast and safe filtration of cell suspensions such as those from tissue dissociation or for flow cytometry.

F-bottom stands for a flat bottom well profile.

F-bottom/chimney well stands for the well profile of a flat well bottom in a chimney-like arrangement. In other words, each well stands on its own. The risk of contamination from sample material being carried over is minimised.

FLUOTRAC is a quality term for immunological products, referring to black microplates (fluorescence measurement).

FourWell Plate is a subdivided plate for microscopic applications facilitating the cultivation of cells and the storage of microscopic slides in an HTS-compatible plate complying with ANSI standards.

Hanging Drop is a technique for protein crystallisation based on > vapour diffusion, where droplets literally hang from the top of an upper substrate.

HiBase is a special plate profile of 384 well > Small Volume and 1536 well microplates. In contrast to the > LoBase profile, the HiBase profile is particularly well suited for top-reading systems, since the measuring optic has a minimal separation from the upper edge of the well in this plate profile.

High binding microplates (= MICROLON 600, FLUOTRAC 600 and LUMITRAC 600) are immunological microplates with a high-binding polystyrene surface. Hydrophilic groups are introduced to the polystyrene surface by physical treatment. The high binding surface contains more hydrophilic groups than the less hydrophilic > medium binding surface.

LoBase is a special plate profile in 1536 well microplates. In contrast to the > HiBase profile, the LoBase profile is particularly well suited for bottom-reading sytems, since the measuring optic has a minimal separation from the well bottom in this plate profile.

LUMITRAC is a quality term for immunological products, referring to white microplates (luminescence measurement).

MASTERBLOCK® is a brand name that stands for polypropylene microplates that are suitable for the storage of non-human sample material. They are also ideally suited for cultivating bacteria or yeast.

Med. binding (medium binding) microplates (= MICROLON 200, FLUOTRAC 200 and LUMITRAC 200) are immunological microplates with a less hydrophobic surface than > high binding microplates.

Microbatch under oil is a method for protein crystallisation where the droplet is covered with oil. The oil generally used is paraffin wax and/or silicone oil. Paraffin wax allows little to no diffusion of water out of the droplet. Hence, all the reagents involved in the crystallisation process, as well as the protein, are present at defined concentrations, and no significant increase of concentration occurs within the crystallisation droplet. When paraffin wax is mixed with silicone oil, it is possible for water to diffuse out of the droplet through the oil and both protein and reagent concentrations increase within the droplet.

 $\label{eq:policy} \begin{array}{l} \mu Clear^{@}(\mbox{Micro-Clear}) \mbox{microplates, in contrast to standard} \\ \mbox{microplates with a solid bottom, have a very thin foil bottom.} \\ \mu Clear^{@}\mbox{microplates are ideal for cell-based test} \\ \mbox{systems, microscopic analyses, as well as for bottom-reading systems.} \end{array}$

MICROLON is a quality term for immunological products, referring to clear microplates (transmission measurement).

Non-binding microplates are characterised by low protein, DNA, RNA and peptide binding properties.

OneWell Plate is a non-divided HTS plate for tissue culture applications complying with the ANSI standards. The plate is also available in a non-TC-treated version for bacteriology.

Sapphire is a quality term for pipette and filter tips. The product family comprises standard pipette tips, standard filter tips as well as a low-retention version of both. All tips are transparent, graduated and allow precise pipetting with maximal recovery. They can be used with all common pipettes.

SCREENSTAR is a quality term for microplates manufactured out of high-quality cycloolefin with an ultra-clear film bottom for high-content and high-throughput screening.

Sitting Drop is a technique for protein crystallisation based on > vapour diffusion, where droplets sit on the bottom of a substrate.

Small Volume is a well profile that was developed in 384 well format for reducing the sample volume. In contrast to the 384 well standard microplate, the sample volume can be considerably reduced, while the detection limit remains the same or is even improved.

TC surface treatment stands for a special physical procedure with which the surfaces of CELLSTAR® products for

adherent cell culture are treated. This treatment leads to the incorporation of polar groups, such as carboxy and hydroxy groups, into the plastic surface making it hydrophilic. This enables the adhesion of cells to the plastic surface.

 $\mbox{\sc U-bottom}$ stands for well profile of wells with round bottom.

UV-Star[®] microplates are made of polyolefin and have a film bottom. In contrast to standard microplates with a solid bottom, they are characterised by an extended transparency range to as low as 200 nm.

Vapour diffusion is the most commonly used method for protein crystallisation. In this method a crystallisation droplet, formed by combining a protein solution with a reagent solution, is incubated together with a larger volume of the same reagent solution within a closed system. The reagent solution can contain a wide range of chemicals, e.g. buffers, salts or precipitating agents. Due to mixing the reagent and protein solutions, the concentration of reagents within the crystallisation droplet becomes lower than the concentration of the reagent solution itself. This causes water to evaporate out of the droplet until equilibrium is reached. During this process, the concentration of protein and chemicals in the crystallisation droplet is continuously rising, and, if optimal conditions have been chosen, protein crystals will begin to form. Vapour diffusion experiments are most often set up as > hanging or > sitting drop.

 $\mbox{\sc V-bottom}$ stands for the well profile of wells with a conically tapered well bottom.

GLOSSARY OF SYMBOLS



NUMERIAL INDEX

400004				
102201118	127279135	227289125	606190164	639160-TRI48
102261118	127280135	227290125	607107163	63996049
102270118	131202142	303321118	607160163	64320146
112101117	131263142	307321118	607160-TRI164	64340146
112201117	132202143	310321118	607180163	64420146
115001222	132263143	318321118	607190164	64440146
115061222	133202144	323070129	608281231	64620146
115070222	133263144	330070129	609101203	64640146
115071222	136101117	332070129	609120203	64920146
115101117	160101117	354070129	609130203	64940146
115201117	160201117	366380210	609171203	650001217
115261123	16316069	366383210	609180203	650061217
115262123	163177122	366384210	609801203	650101179
12016069	163288125	366385210	609820203	65016061
120161123	163289125	366386210	609830203	650161179
120180123	163290125	373250231	609871203	65018061
12019069	16416069	373270231	612301165	65018564
121261133	164161122	373273231	612361165	650201180
		373274231		
121263134	164180122		612362165	650209181
121263-128139	169101117	373275231	613101213	650261181
121277134	186171123	373276231	614101213	650901193
121278134	187101117	373277231	616201209	65097088
121279134	187201117	373281231	616283209	65097988
121280134	187261123	381061201	618201209	651001217
122261134	187262123	381070201	622201210	651061217
122263134	188161123	381080201	622261210	651101179
122263-2D3139	188171123	381081201	623201209	65116061
122263-2DG138	18824169	385270145	627102169	651161179
122263-TRI134	188241122	385273145	62716048	65118061
122277134	188261120	385274145	62717048	651201181
122278134	188261-N120	385275145	62786050	651209181
122279134		385276145		65152493
	188271120		627860105	
122280134	188271-N120	385277145	62786150	65159693
123261133	188271-TRI 120	385278145	627861105	651901193
123263133	188280120	385279145	62787051	65197088
123263-2D3 138	188281120	420161170	627870105	652201233
123263-2DG138	188283120	420180170	62787150	652210233
123263-TRI133	188285120	421084170	627871105	652250233
123277133	19116069	421161170	62796048	652260233
123278133	191161123	421180170	62796550	652270233
123279133	191180123	501870203	627965105	652290233
123280133	201150127	503850204	62797551	653102204
124261135	201152127	503870204	627975105	653180224
124263135	201170127	54107045	62797950	653190224
124263-128139	201172127	54108045	62797987	654102204
124273136	205101129	54200071	628102169	654180224
124274136	210261121	54204071	62816048	655001217
124275136	210270121	54207071	628160-TRI48	655061217
124276136	217101129	54210071	62891049	65507361
126261134	224170127	54212071	62892049	655074179
126263134	225170127	54214071	62893049	655075179
126263-2D1139	225180127	54217071	62894049	655076179
126263-2DG138	22724569	543078107	62895049	655077179
126263-TRI134	227245122	543079107	62896048	65507961
126277134	227250125	543978107	62897950	655080217
126278134	227255125	543979107	62897987	655081217
126279134	227261121	604107163	629102169	65508361
126280134	227261-TRI121	604160163	629161169	65508661
127261135	227270121	604160-TRI164	629180169	65508762
127263135	227280121	604181163	632180169	65508862
127263-2D1139	227281121	606107163	633180169	65509062
127263-2DG138	227283121	606160163	635102169	655094179
127277135	227285121	606160-TRI164	639102169	655095179
127278135	227288125	606180163	63916048	655096179

655097179	65764095	66261058	673277230	68167275
65509862	65764157	66261096	673281231	68167575
655101179	65764195	66263058	673283231	682201229
65516061	65782493	66263096	675001217	682273229
655161179	65782591	66263158	675061217	682274229
65516261	65783091	66263196	675074180	682275230
65518061	65784091	66263858	675075180	682276230
65518261	65784193	66263896	675076180	682277230
65518564	65784393	66264058	675077180	682281230
655201181	65784693	66264096	67508362	68261275
				68261575
655209181	65784793	66264158	67508662	
655801198	65785093	66264196	67509062	68262475
65583090	65785293	66282591	675096180	68262575
65584090	65786092	66283091	675101180	68266075
65584190	65789693	66284091	675161180	68267075
65584692	65793054	662892113	67518062	68267275
65584992	65794054	66293054	675801198	68267875
65585090	65795054	66294054	676001200	683201229
655866111	65796054	66295054	676040200	683271229
655891109			676050200	
	65797055	66296054		683273229
655892113	65797088	66297055	676051201	683274229
655900193	65817041	66297088	676070200	683275229
655901193	658170-TRI41	664102169	676090200	683276229
655903193	65817541	66416048	67710253	683277229
655904193	658175-TRI41	664160-TRI48	67718053	684201230
655906193	65819043	66491049	67797055	688102170
65593063	658190-TRI43	66492049	67797088	69016041
65593663	65819543	66494049	67798054	690160-TRI41
65594063	658195-TRI43	66495049	67810178	69017041
65594463				69017541
	65891043	66496049	678101-CF180	
65594663	65892043	66497050	678101-EXF79	690175-TRI41
65594863	65894043	66497087	67810478	69019043
65595063	65895042	66510253	678104-CF180	690190-TRI43
65595663	65897042	66511059	678104-CF280	69019543
65597088	65897542	66511097	678104-EXF79	690195-TRI43
65597688	65898044	66518053	67810878	69091043
655976-SIN88	65898087	66561057	678108-CF180	69092043
65598063	65898544	66561095	678108-CF280	69094043
655981109	65898587	66563058	678108-EXF79	69095042
65598363	659180224	66563096	67811278	69096042
65598663				69097542
	659190224	66563158	678112-CF180	
655990196	66016041	66563196	678112-CF280	69098044
655995196	660160-TRI41	66563858	678112-EXF79	69098087
655997196	66017541	66563896	67811678	69098544
65610166	660175-TRI41	66564057	678116-CF180	69098587
656101200	66019043	66564095	678116-CF280	69110167
65616166	660190-TRI43	66564157	678116-EXF79	691101200
656161200	66096042	66564195	67812478	69116167
65617066	66097542	66597088	678124-CF180	691161200
656170200	66098044	66598054	678124-CF280	700361165
65617166	66098087	667201209	678124-EXF79	700370165
656171200	66098544	668102211	67814078	705063220
65619067	66098587	669285233	678140-CF180	705065220
656190200	66116041	670102170	678140-CF280	705066220
65619167	66117541	67018051	678140-EXF79	705070220
656191200	66119043	67019051	67890478	705071220
657102170	66119543	671201229	67890878	705073220
65711059	661195-TRI43	671221230	67891278	705074220
65711097	66191043	671273229	67891678	705075220
65716053	66192043	671274229	67892478	705076220
65718553	66194043	671275229	67894078	710107163
65761057	66195042	671276229	68019076	710160163
65761095	66198044	671277229	68019576	710160-TRI164
65763057	66198087	671281229	68064575	710180163
65763095	66198544	673210230	68064875	710183164
65763157	66198587	673271230	68065875	716201209
65763195	66210253	673273230	68066075	717201209
65763857	662145203	673274230	68066575	722201209
65763895	662150203	673275230	68066875	724170 118
65764057	66216053	673276230	68167075	724411 128
00/070	552 100	0,02,0200	551070	/ ∠¬¬ ∠0

724412128	775352157	781186183	784201184	976587142
724413 128	775353157	781201184	784900194	976588142
724414128	775354157	781201-906 184	784904194	976589142
725201209	775355157	781201-906 191	78494665	977501145
725261209	775361157	781209184	785201233	977561143
729101211	775362157	781270189	785285233	977563143
731101171	775363157	781271189	785290233	977564143
731161171		781280184		977565143
	775364157		786201188	
731165171	775365157	781801198	786261188	977566143
731170171	776350158	78183091	78797988	977567143
731171171	776351158	78184090	789866111	977568143
731175171	776352158	78184190	792870-906191	977569143
731180171	776353158	78184692	793855191	977570143
731181171	776354158	78184992	802202136	977580143
731185171	776355158	78185090	802203136	977583143
745290161	776361158	781866111	802204136	977584143
754061219	776362158	781892113	802206136	977585143
754070219	776363158	781900193	802225136	977586143
756070219	776364158	781901193	802501136	977587143
756071219	776365158	781903193	802506140	977588143
760107163	777350159	781904193	802576140	977589143
760160163	777351159	781906194	803202139	978501145
	777352159	78193065		978561144
760160-TRI164			803270139	
760180163	777354159	78193665	803277139	978563144
762070219	777355159	78194065	843070237	978564144
762071219	777361159	78194465	844070238	978565144
762074219	777362159	78194565	845070238	978566144
762075219	777364159	78194665	846070237	978567144
762076219	777365159	78194865	847070165	978568144
762077219	778350160	78195065	848070239	978569144
767070219	778351160	78195665	848902239	978570144
767071219	778352160	78197088	848913239	978580144
768160163	778353160	78197488	848915239	978583144
768160-TRI164	778354160	78197688	848916239	978584144
768180163	778355160	781976-SIN88	848921239	978585144
769190224	778361160	78198365	848923239	978586144
770310161	778362160	78198665	848950239	978587144
770320161	778363160	781990196	849050146	978588144
770330161	778364160	781995196	849070146	978589144
770340161	778365160	781997196	852070146	89000002 151
770370161	77916044	782061186	87807182	89000010151
770380161	77919044	78207366	87807281	89000020151
771350154	780201188	782074186	87807381	89000099152
771351154	780215188	782075186	87807481	89000100151
771352154	780261188	782076186	87807582	89000200151
771353154	780270188	782077186	95070045	89000500151
771354154	780271188	78207866	95170045	89000810151
771355154	780285188	78208066	95270045	89000820151
771361154	781061183	78208666	960161128	89001000151
771362154	78107364	78209266	960177129	89001210152
771363154	781074183	78209366	968162128	89001220152
771364154				
	781075183	782095186	968177128	89008200151
771365154	781076183	782096186	975502188	89008300151
772350155	781077183	782097186	975561188	89010000151
772351155	78107964	782101186	975570188	89012200152
772352155	78108064	78218066	976501145	89012300152
772353155	78108664	782261189	976561142	9607730751
				30077007
772354155	78109065	782270189	976563142	
772355155	78109164	782855191	976564142	
772361155	78109264	782900194	976565142	
772362155	78109364	782904194	976566142	
772363155	781095183	78294666	976567142	
772364155	781096183	783892113	976568142	
	781097183		976569142	
772365155		784075184		
773353156	78109864	784075-25 184	976570142	
773363156	781101183	784076184	976580142	
774353156	781162183	784076-25 184	976583142	
774363156	78116564	78408065	976584142	
775350157	78118264	78408665	976585142	
	781185183		976586142	
775351157	701100100	784101184	070300142	

GLOSSARY

1536 Well Microplates	186
384 Well Microplates	
3D Cell Culture	
48-way Datamatrix Cryo Rack	
81-way Datamatrix Cryo Rack	
96 Well ELISA Microplates	
·	
96 Well ELISA Strip Plates	
96 Well Microplates	
96-way Datamatrix Cryo Rack	145
Advanced TC	54
Analyser Cups	211
AutoFlask - Cell Culture Flask	44
Biobanking Tubes	142,143,144,145
CELLCOAT®	54
CELLdisc	
CELLevator	
CELLhandle	
CELLreactor	
CELLring	
CELLstage	
CELLview Dish	
CELLview Distr	
CELLview Slide	
Ceaprene stopper	
Cell Culture	
Cell Culture Dishes	
Cell Culture Dishes - Advanced TC	
Cell Culture Dishes - CELLCOAT®	
Cell Culture Dishes - Cell-Repellent Surface	
Cell Culture Flasks	
Cell Culture Flasks - Advanced TC	
Cell Culture Flasks - CELLCOAT®	
Cell Culture Inserts	
Cell Culture Microplates	64,65,66,88
Cell Culture Roller Bottles	74,75,76
Cell Culture Tubes	
Cell Scraper	45
Cell Strainer	71
Cell-Repellent Surface	44,55,87,88
Centrifuge Tube	120,121
Containers for Plant Cultures	128
Cryo Tubes	
Cryo storage box	
Cryo.s	
Cryo.s Biobanking Tubes	
Cryo.s Decapper	
Cryo.s Rack Scanner	
Cryo.s with Barcode	
Drosophila Containers	129
FACVatrainer Call Strain	74
EASYstrainer Cell Strainers	
ELISA Strip Plates	219,220

FourWell Plate	51
Grip stopper	118
HTS-Microplates	179,180,183,184,186
Immuno Tubes	222
Inserts for Mini Block Heater	239
Leucosep	125
Macro Pipette Tip	
Magnetic 3D Cell Culture	90,91,92,93
Mass Cell Culture	
MaxiPette	
Media Bottles	
Microplate Centrifuge	
Microplates	
Mini Vortex Mixer	
Multi-Channel Pipettes	
Multipurpose Containers / Beakers	
Multiwell Plates	
OncoQuick®	125
OneWell Plate	
Pasteur / Serum Pipettes	165
Pipette Carrousel	
Polypropylene Tubes	
Polystyrene Tubes	117,122,123
Polystyrene drosophila containers	
Products for Microscopy	
Reaction Tubes	209
Roller Bottles	
SCREENSTAR Microplates	111
Screw Caps	
Semi-micro / Macro Cuvette	
SensoPlate	113
Serological Pipettes	
Single-Channel Pipettes	151
Support Rack	136
Suspension Culture Flasks	43
Terasaki Plates	224
ThinCert® Cell Culture Inserts	57,58
ThinCert® Plate	59
Transfer Pipettes	165
Tube	69,122,123
Tubes / Multipurpose Beakers	
Tubes with closures	122,123
VtMin	0.70



making a difference



GREINER BIO-ONE GMBH FRICKENHAUSEN, GERMANY

PHONE +49 7022 948-0 FAX +49 7022 948-514 E-MAIL info@de.gbo.com



GREINER BIO-ONE IS A GLOBAL PLAYER.

FIND THE CONTACT DETAILS OF YOUR LOCAL PARTNER ON OUR WEBSITE.



Devices of Greiner Bio-One are to be used by properly qualified persons only in accordance with the relevant Instructions for Use (IFU), where applicable.

For more information contact your local Greiner Bio-One sales representative or visit our website (www.gbo.com).

All information is provided without guarantee despite careful processing. Any liability, warranty or guarantee of Greiner Bio-One GmbH is excluded. All rights, errors and changes are reserved. If not stated otherwise, Greiner Bio-One GmbH has all copyrights and/or other (user-) rights in this documents, in particular to signs such as the mentioned (word-picture-) brands and logos. Any use, duplication or any other use of the rights of Greiner Bio-One GmbH is expressly prohibited. Media owner: Greiner Bio-One GmbH / Represented by Managing Directors Jakob Breuer and Heinz Schmid. The company is registered in the Commercial Register at the first instance court in Stuttgart, HRB 224604 / VAT Number: DE812585719.

