


Plastic Body Nebulizer Parts for AAnalyst AA Instruments

AAnalyst 100/300/200/400/700/800* High-Sensitivity Nebulizer Parts

For Nebulizers with Part Nos. **B0505480** and **B0505590**

Description	Part No.
1 Impact Bead	B0505086
2 Solvent Resistant O-Ring for Nebulizer Body	B0174378
3 Nebulizer Body	B0506666
4 Spring	B0165332
5 Solvent Resistant O-Ring for Capillary Assembly	B0190672
6 Pt/Ir Capillary Assembly	B0501022
7 Locking Ring	B0501025
8 Fine Adjustment Nut	B0501026

*This nebulizer requires the use of end cap assembly

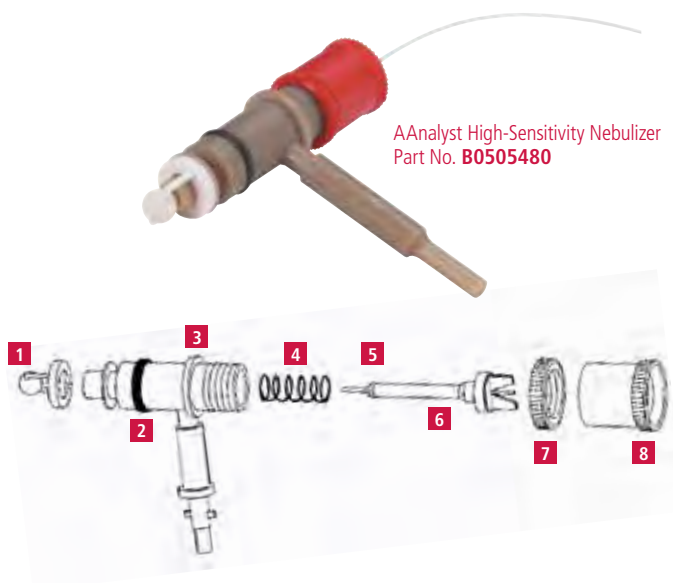

AAnalyst 200/400/700*/800* High-Sensitivity Nebulizer Parts

For Nebulizers with Part Nos. **N3150188** and **N2010250**

Description	Part No.
1 Impact Bead	B0505086
Nebulizer Spacer (not pictured)	B3150531
2 O-Rings for Nebulizer Body (Each nebulizer requires a Qty. of 2)	09902015
Solvent Resistant O-Rings for Nebulizer Body (Each nebulizer requires a Qty. of 2)	09200276
3 Nebulized Body	—
4 Spring	09920483
5 O-Ring for Capillary Assembly	09926127
Solvent Resistant O-Ring for Capillary Assembly	09200272
6 Pt/Ir Capillary Assembly w/O-Ring	B3150525
Tantalum Capillary Assembly w/Solvent Resistant O-Ring and Spring	N3160146
7 Locking Ring	N3151205
8 Fine Adjustment Nut	B3150462
N/A O-Ring for Sample Tube (Each sample tube requires a Qty. 2)	B0507708
N/A Solvent Resistant O-Ring for Sample Tube (Each sample tube a Qty. 2)	09200369
N/A Solvent Resistant O-Ring Kit (Includes 1 x 09200272 , 2 x 09200276 and 2 x 09200369)	N2010246

*Requires End Cap Assembly **N0370392**

