


Nebulizer Assemblies and Spares

Metal Body Nebulizer Parts for PinAAcle AA Instruments

For Nebulizers with Part Nos. **N3160143** and **N3160158**

Description	Part No.
1 O-Ring sealing Nebulizer to End Cap	09902239
Solvent Resistant O-Ring Sealing Nebulizer to End Cap	09200274
2 Cap	N3161057
3 Venturi	03031810
4 O-Ring for Venturi	09902005
Solvent Resistant O-Ring for Venturi	09200273
5 O-Rings for Nebulizer Body (Each nebulizer requires a quantity of 2)	09902015
Solvent Resistant O-Rings for Nebulizer Body (Each nebulizer requires a quantity of 2)	09200276
6 Nebulizer Body	N3160133
7 Nebulizer Insert	03031812
8 O-Ring for Nebulizer Insert	09902047
Solvent Resistant O-Ring for Nebulizer Insert	09200273
9 Locking Ring	N3161006
10 Spring	03031806
11 O-Ring for Capillary Needle Assembly	09902102
Solvent Resistant O-Ring for Capillary Needle Assy.	09200271
12 Capillary Needle Assembly W/O-Ring	03030354
Solvent Resistant Capillary Needle Assembly W/O-Ring	N3160153
13 PTFE Washer	03031802
14 Fine Adjustment Nut	03031460
15 Sample Capillary	09908265

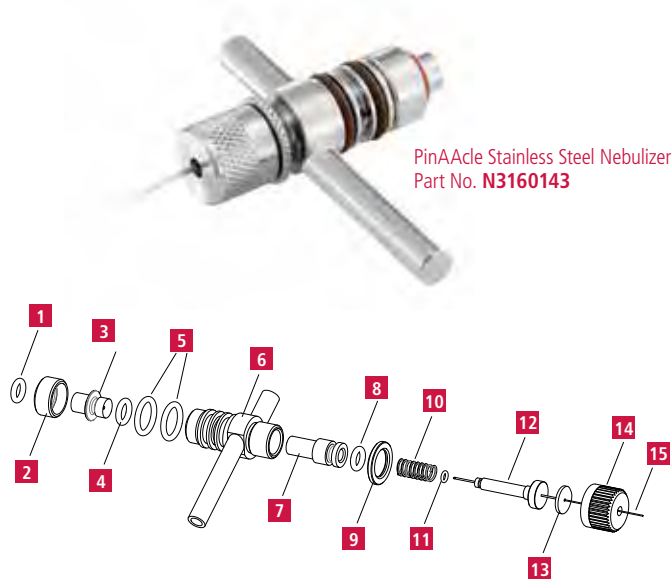


PinAAcle Stainless Steel Nebulizer
Part No. **N3160143**

Plastic Body Nebulizer Parts for PinAAcle AA Instrument

For Nebulizers with Part Nos. **N3160144**, **N3160152** and **N3160159**

Description	Part No.
1 Impact Bead	B0505086
2 Spacer	B3150531
3 O-Rings for Nebulizer Body (Each nebulizer requires a quantity of 2)	09902015
Solvent Resistant O-Rings for Nebulizer Body (Each nebulizer requires a quantity of 2)	09200276
4 Plastic Nebulizer Body	B3161049
5 Spring	09920483
6 Locking Ring	N3151205
7 O-Ring for Capillary Needle Assembly	09926127
Solvent Resistant O-Ring for Capillary Needle Assembly	09200272
8 Pt/Ir Capillary Needle Assembly W/Spring & O-Ring	B3150525
Solvent Resistant Pt/Ir Capillary Needle Assembly W/ Spring & O-Ring	N3160154
9 Fine Adjustment Nut	B3150462
10 Sample Tubing	B3150458

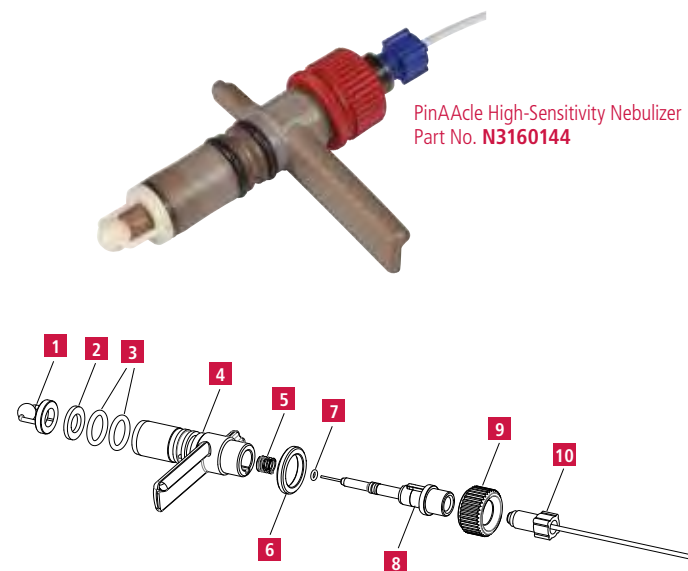

PinAAcle High-Sensitivity Nebulizer
Part No. **N3160144**